QUICK REFERENCE



EVOS

Limited Product Warranty

Life Technologies Corporation and/or its affiliate(s) warrant their products as set forth in the Life Technologies' General Terms and Conditions of Sale found on Life Technologies' website at www.lifetechnologies.com/termsandconditions. If you have any questions, please contact Life Technologies at www.lifetechnologies.com/support.

DISCLAIMER

LIFE TECHNOLOGIES CORPORATION AND/OR ITS AFFILIATE(S) DISCLAIM ALL WARRANTIES WITH RESPECT TO THIS DOCUMENT, EXPRESSED OR IMPLIED, INCLUDING BUT NOT LIMITED TO THOSE OF MERCHANTABILITY, FITNESS FOR A PARTICULAR PURPOSE, OR NON-INFRINGEMENT. TO THE EXTENT ALLOWED BY LAW, IN NO EVENT SHALL LIFE TECHNOLOGIES AND/OR ITS AFFILIATE(S) BE LIABLE, WHETHER IN CONTRACT, TORT, WARRANTY, OR UNDER ANY STATUTE OR ON ANY OTHER BASIS FOR SPECIAL, INCIDENTAL, INDIRECT, PUNITIVE, MULTIPLE OR CONSEQUENTIAL DAMAGES IN CONNECTION WITH OR ARISING FROM THIS DOCUMENT, INCLUDING BUT NOT LIMITED TO THE USE THEREOF.

Limited Use Label License: Research Use Only

The purchase of this product conveys to the purchaser the limited, non-transferable right to use the purchased amount of the product only to perform internal research for the sole benefit of the purchaser. No right to resell this product or any of its components is conveyed expressly, by implication, or by estoppel. This product is for internal research purposes only and is not for use in commercial applications of any kind, including, without limitation, quality control and commercial services such as reporting the results of purchaser's activities for a fee or other form of consideration. For information on obtaining additional righs, please contact outlicensing@lifetech.com or Out Licensing, Life Technologies Corporation, 5791 Van Allen Way, Carlsbad, California 92008.

Trademarks

The trademarks mentioned herein are the property of Life Technologies Corporation and/or its affiliate(s) or their respective owners. © 2013 Life Technologies Corporation. All rights reserved.

Headquarters

5791 Van Allen Way | Carlsbad, CA 92008 USA | Phone +1 760 603 7200 | Toll Free in USA 800 955 6288

For support visit lifetechnologies.com/support or email techsupport@lifetech.com


lifetechnologies.com 18 March 2012

life nlonies





For Research Use Only. Not for use in diagnostic procedures.

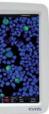

EVOS[®] FL EVOS® FL Color

Imaging Systems

for Fluorescence and Transmitted Light Applications



Catalog Numbers AME4300, AMEFC4300

Doc. Part no. ZP-PKGA-0494 Pub. no. MAN0007717 **Rev.** 2.0



Overview

The EVOS® FL and EVOS® FL Color Imaging **Systems** have two types of controls: manual and onscreen. Manual controls include the the stage X-Y axis knobs, focusing knobs, objective selection wheel, phase annuli selector, and the light cube selection lever. Onscreen controls are located in the **control bar** at the bottom of the display screen. The **channel bar** at the top of the display screen shows the selected light cube or transmitted light position.

~恭


CAUTION! UV HAZARD. Avoid exposure to beam and use protective shields. NEVER look directly at UV light!

Basic Operation

Note: For detailed instructions, refer to the EVOS® FL Imaging System User Guide, which is provided on the USB flash drive. You can also dowload it from the EVOS[®] FL Imaging System product page at www.lifetechnologies.com.

- 1. Turn on the instrument using the **power switch ()** on the right side of the base.
- 2. Plug a USB flash drive into one of the USB ports 2 on the right side of the instrument.
- 3. Place the **sample (3)** on the stage, using a vessel holder if needed.
- 4. Set magnification with the **objective selection** wheel 4 on the front of the instrument.
- 5. Pull the light cube selection lever (5) (left side of base) all the way toward the front of the instrument *(the Channel Bar will* highlight the "Transmitted" position).

2

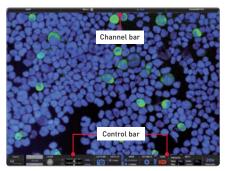





- **6.** Turn on illumination with the **LIGHT ON button** located on the left side of the control bar.
- 7. Focus the sample with focusing knobs 6.
- **8.** *Optional:* To take a picture of the transmitted light image, click the **Capture button** on the control bar.
- 9. Place the light shield box 7 on the stage, over the sample.
- 10. Move the light cube selection lever 5 to the desired fluorescence channel (the Channel Bar will highlight the selected light cube).
- 11. With the Find & Focus tab active, turn on fluorescence illumination using the **LIGHT** ON button.
- **12.** Adjust the focus as necessary.
- 13. Adjust the Illumination Intensity slider on the control bar as needed.
- 14. Click the Capture button.
- **15.** Repeat steps 10–14 to acquire each fluorescence channel.
- 16. Click the **Overlay tab** to show all channels in color overlay mode.
- 17. Adjust Brightness and Contrast for each channel to bring them into desired balance.
- 18. Click the Save button to save the color image (refer to the EVOS® FL Imaging System User Guide).







3



8. Color Option Scalebar/Toolbar Options 9. 10. Setting Control Button

5.

6

2. Login Button

3. Control Bar Tabs

Actual

• Overlay

4. Light ON/OFF Button

Illumination Slider

7. Image Capture Button

Exposure Time Slider

Find & Focus

11. Save Image Button

Helpful Tips

- 12. Info Display Bar
- 13. Selected Objective



In Find & Focus or Actual Mode, the Color option (for EVOS® FL, Cat. no. AME4300) can be turned off to display a grayscale image. This often shows more details than a color image.

For EVOS® FL Color (Cat. no. AMEFC4300), use the Color Adjustment button to fine tune your live image Brightness, Contrast, Saturation, and Hue prior to capture.

▶ In Find & Focus Mode, the exposure time is set to 100 ms to assist real-time focusing, moving the stage, etc. The illumination level is approximately 60% of the amount used for image capture, in order to minimize photobleaching and phototoxicity. Clicking Capture results in brighter illumination and a longer exposure time during image capture to provide a high-quality image.

In Actual Mode, turning on the illumination results in full-powered illumination and actual exposure times for live viewing of the sample. With longer exposure times (more than 200 ms) there will be a lag between moving the focus knob and seeing the focus change onscreen.

Onscreen Controls

1. Active Channel (highlighted)

- Exposure Time Slider

- Overlay Color Dialogue Box

4

