

NanoDrop Micro-UV/Vis Spectrophotometer

# NanoDrop Eight

# **User Guide**

M020 NanoDrop Eight UG Revision Date March 2022



© 2022 Thermo Fisher Scientific Inc. All rights reserved.

	For U.S. Technical Support, please contact:	For International Support, please contact:
3411 S Tatnall	Thermo Fisher Scientific 3411 Silverside Road	http://www.thermofisher.com/ NanoDropSupport
	Tatnall Building, Suite 100 Wilmington, DE 19810 U.S.A.	Contact your local distributor. For contact information go to:
	Telephone: 302 479 7707 Toll Free: 1 877 724 7690 (U.S. & Canada only) E-mail: nanodrop@thermofisher.com	http://www.thermofisher.com/ NanoDropDistributors
	Thermo Fisher Scientific Inc. provides this docum	ent to its customers with a product purchase t

Thermo Fisher Scientific Inc. provides this document to its customers with a product purchase to use in the product operation. This document is copyright protected and any reproduction of the whole or any part of this document is strictly prohibited, except with the written authorization of Thermo Fisher Scientific Inc.

The contents of this document are subject to change without notice. All technical information in this document is for reference purposes only. System configurations and specifications in this document supersede all previous information received by the purchaser.

Thermo Fisher Scientific Inc. makes no representations that this document is complete, accurate or error-free and assumes no responsibility and will not be liable for any errors, omissions, damage or loss that might result from any use of this document, even if the information in the document is followed properly.

This document is not part of any sales contract between Thermo Fisher Scientific Inc. and a purchaser. This document shall in no way govern or modify any Terms and Conditions of Sale, which Terms and Conditions of Sale shall govern all conflicting information between the two documents.

For Research Use Only. This instrument or accessory is not a medical device and is not intended to be used for the prevention, diagnosis, treatment or cure of disease.



**WARNING** Avoid an explosion or fire hazard. This instrument or accessory is not designed for use in an explosive atmosphere.

**C** -

# Contents

Chapter 1	About the Spectrophotometer	7
-	Features	
	USB-B port	
	Accessories	
	PR-1 Pedestal Reconditioning Kit	
	PV-8 Performance Verification Kit	
	Instrument Detection Limits	
Chapter 2	Instrument Set up	11
	Register Your Instrument	11
	Computer Requirements	11
	Update Software	12
	Technical Support	12
	For U.S./Canada Support, please contact:	12
	For International Support, please contact:	12
Chapter 3	Application Measurement Ranges	
	Detection Limits for All Applications	
	Detection limits for pre-defined dyes	14
Chapter 4	Nucleic Acid Applications	15
-	Measure dsDNA, ssDNA or RNA	15
	Measure dsDNA, ssDNA or RNA	15
	Best practices for nucleic acid measurements	17
	Nucleic Acid Reported Results	19
	Settings for Nucleic Acid Measurements	
	Calculations for Nucleic Acid Measurements	
	Measure Microarray	25
	Measure Microarray Samples.	
	Microarray Reported Results	
	Settings for Microarray Measurements	
	Calculations for Microarray Measurements.	

### Contents


	Measure using a Custom Factor	37
	Measure Nucleic Acid using a Custom Factor	37
	Custom Factor Reported Results	39
	Settings for Nucleic Acid Measurements using a Custom Factor	40
	Detection Limits for Nucleic Acid Measurements using a Custom	
	Factor	41
	Measure Oligo DNA or Oligo RNA	43
	Measure Oligo DNA or Oligo RNA	
	Oligo Reported Results	
	Settings for Oligo DNA and Oligo RNA Measurements	
	Detection Limits for Oligo DNA and Oligo RNA Measurements	
	Calculations for Oligo DNA and Oligo RNA Measurements	
	5 5	
Chapter 5	Protein Applications	53
-	Measure Protein A280	53
	Measure Protein Concentration at A280	53
	Best practices for protein measurements	55
	Protein A280 Reported Results	
	Settings for Protein A280 Measurements	
	Protein editor	
	Detection Limits for Protein A280 Measurements	63
	Calculations for Protein A280 Measurements	63
	Measure Protein A205	67
	Measure Protein Concentration at A205	67
	Protein A205 Reported Results	69
	Settings for Protein A205 Measurements	70
	Calculations for Protein A205 Measurements	71
	Measure Proteins and Labels	73
	Measure Labeled Protein Samples	73
	Proteins & Labels Reported Results	75
	Settings for Proteins and Labels Measurements	76
	Detection Limits for Proteins and Labels Measurements	78
	Calculations for Proteins and Labels Measurements	79
Chapter 6	Custom Applications	
	Measure UV-Vis	
	Measure UV-Vis	
	Best practices for UV-Vis measurements	83
	UV-Vis Reported Results	
	Settings for UV-Vis Measurements	
	Measure Custom	87
	Measure using a Custom Method	87
	Delete Custom Method	88
	Custom Method Reported Results	89
	Manage Custom Methods	90

**Thermo Scientific** 

### Contents

Chapter 7	Learning Center	101
-	Micro-Volume Sampling—How it Works	102
	Set Up the Instrument	104
	Connect Power	104
	Connect to a Computer	104
	Operating Specifications	105
	Measure a Micro-Volume Sample	106
	Best practices for micro-volume measurements	106
	Recommended sample volumes	107
	Sample Modes	110
	Module startup	110
	Sample Plate Map	
	Prepare Samples and Blanks	110
	Preparing Samples	
	Run a Blanking Cycle	
	Basic Instrument Operations	
	NanoDrop Eight Home Screen	
	NanoDrop Eight Measurement Screens	
	Measurement Screen Display Options	
	NanoDrop Eight General Operations	
	View Acclaro Sample Intelligence Information.	
	On-Demand Technical Support	
	System Settings	
	Protein Editor	
	Measurement Screen Display Options	148
Chapter 8	Maintenance	151
enapter e	Maintenance Schedule	
	Daily Maintenance	
	Periodic Maintenance	
	Every 6 Months	
	Maintaining the Pedestals	
	Cleaning the Pedestals	
	Reconditioning the Pedestals.	
	Decontaminating the Instrument	
	Instrument Diagnostics	
	Intensity Check	
	Performance Verification	

# **About the Spectrophotometer**

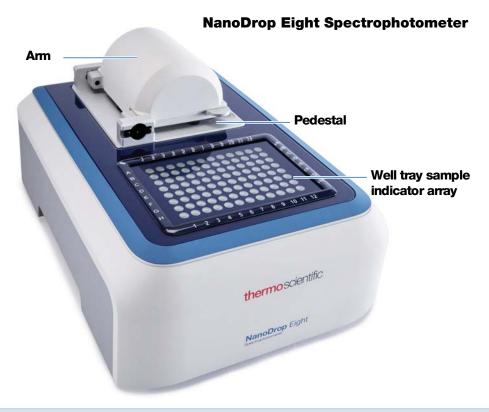

**Note** Locate the instrument away from air vents and exhaust fans to minimize evaporation

The Thermo Scientific<sup>™</sup> NanoDrop<sup>™</sup> Eight is a UV-Visible spectrophotometer developed for micro-volume analysis of a wide variety of analytes. The patented sample retention system enables the measurement of highly concentrated samples without the need for dilutions. The instrument accurately measures up to 8 individual 1 µl samples in one measurement cycle. The software allows the user to measure samples using either the full 8-channel mode or a convenient single sample mode.

The NanoDrop Eight software is installed on a local PC used to control the instrument and view data.

**1** About the Spectrophotometer Features

**Note** Before operating a NanoDrop Eight instrument, please read the safety and operating precautions and then follow their recommendations when using the instrument.

## **Features**

TheNanoDrop Eight spectrophotometer features the patented micro-volume sample retention system.

### **USB-B** port

Connect the NanoDrop Eight to your PC via USB to operate the instrument using the PC software.

### **Accessories**

This section lists the accessories included for use with the NanoDrop Eight.

### **PR-1 Pedestal Reconditioning Kit**



Specially formulated conditioning compound that can be applied to the pedestals to restore them to a hydrophobic state (required to achieve adequate surface tension for accurate sample measurements). The kit includes conditioning compound and applicators. For more information, see Reconditioning the Pedestals.

### **PV-8 Performance Verification Kit**

Kit containing plastic consumables and liquid photometric standard used to check instrument performance. For more information, see Performance Verification.

# **Instrument Detection Limits**



Upper Detection Limit (10 mm Equivalent Absorbance)	
12.5	
90	
200	

2

# **Instrument Set up**

## **Register Your Instrument**

Register your instrument to receive e-mail updates on software and accessories for the NanoDrop Eight instrument. An Internet connection is required for registration.

### To register your instrument

From any PC that is connected to the Internet, use any web browser to navigate to www.thermofisher.com/nanodropsw.

On the website, locate the Register Your Instrument button, and follow the instructions to register the instrument.

### **Computer Requirements**

### **Required Windows Operating System**

• Windows 10 Enterprise or Professional, build 1607 or greater

### **Minimum hardware configuration**

- 2.0 GHz dual-core processor enabled
- 4 GB RAM with system managed memory enabled
- 5 GB available on drive C
- Display resolution 1366 × 768

### **Recommended hardware configuration**

- 2.33 GHz 4-core processor enabled (or greater)
- 8 GB RAM with system managed memory enabled (or greater)
- 200 GB available on drive C (or greater)

# **Update Software**

Quickly and easily download and install the latest NanoDrop Eight software and release notes from our website. Follow the steps to update or upgrade the NanoDrop Eight software on a personal computer (PC). An Internet connection is required to download software. Navigate to www.thermofisher.com/nanodropsw and click on the NanoDrop Eight software tab. Follow the Instrument Software Download instructions.

### To install or update NanoDrop Eight software on a PC

- 1. Insert the USB flash drive containing the installer software into an available USB port on your PC, or open the installation folder downloaded from the internet.
- 2. Launch **Start.exe** and click the **Software** button. The software installer will run.

## **Technical Support**

### For U.S./Canada Support, please contact:

Thermo Fisher Scientific 3411 Silverside Road Tatnall Building, Suite 100 Wilmington, DE 19810 U.S.A.

Telephone: 302 479 7707 Toll Free: 1 877 724 7690 (U.S. & Canada only) Fax: 302 792 7155 E-mail: nanodrop@thermofisher.com Website: www.thermofisher.com/nanodrop

### For International Support, please contact:

Contact your local distributor. For contact information go to:

http://www.thermofisher.com/NanoDropDistributors

If you are experiencing an issue with your system, refer to the troubleshooting information. If the issue persists, contact us. If you are outside the U.S.A. and Canada, please contact your local distributor.

If your instrument requires maintenance or repair, contact us or your local distributor.

# **Application Measurement Ranges**

# **Detection Limits for All Applications**



**Note** Detection limits provided in the following tables are approximate and apply to micro-volume measurements only; they are based on the instrument's photometric absorbance range (10 mm equivalent) of 0.04–200 A.

Sample Type	Lower Detection Limit	Upper Detection Limit	Typical Reproducibility <sup>a</sup>
dsDNA	2.0 ng/µL	10,000 ng/µL	±2.0 ng/μL for sample concentrations between 2.0 and 100 ng/μL samples; ±2% for samples >100 ng/μL
RNA	1.6 ng/μL	8,000 ng/µL	±2.0 ng/μL for sample concentrations between 2.0 and 100 ng/μL samples; ±2% for samples >100 ng/μL
DNA Microarray (ssDNA)	1.3 ng/µL	495 ng/µL	±2.0 ng/μL for sample concentrations between 2.0 and 100 ng/μL samples; ±2% for samples >100 ng/μL

### **3** Application Measurement Ranges Detection Limits for All Applications

Sample Type	Lower Detection Limit	Upper Detection Limit	Typical Reproducibility <sup>a</sup>
Purified BSA by Protein A280	0.06 mg/mL	300 mg/mL	±0.10 mg/mL (for 0.10–10 mg/mL samples); ±2% for samples >10 mg/mL
IgG by Protein A280	0.03 mg/mL	145 mg/mL	
Purified BSA by Proteins & Labels	0.06 mg/mL (pedestal)	19 mg/mL (pedestal)	±0.10 mg/mL for 0.10–10 mg/mL samples
	0.006 mg/mL (cuvette)		

<sup>a</sup> Based on five replicates (SD=ng/µL; CV=%)

## **Detection limits for pre-defined dyes**

Sample Type	Lower Detection Limit	Upper Detection Limit <sup>a</sup>	Typical Reproducibility <sup>b</sup>
Cy3, Cy3.5, Alexa Fluor 555, Alexa Fluor 660	0.2 pmol/µL	100 pmol/µL	±0.20 pmol/µL for sample concentrations between 0.20 and 4.0 pmol/µL; ±2% for samples >4.0 pmol/µL
Cy5, Cy5.5, Alexa Fluor 647	0.12 pmol/µL	60 pmol/µL	±0.12 pmol/µL for sample concentrations between 0.12 and 2.4 pmol/µL; ±2% for samples >2.4 pmol/µL
Alexa Fluor 488, Alexa Fluor 594	0.4 pmol/µL	215 pmol/µL	±0.40 pmol/µL for sample concentrations between 0.40 and 8.0 pmol/µL; ±2% for samples >8.0 pmol/µL
Alexa Fluor 546	0.3 pmol/µL	145 pmol/µL	±0.30 pmol/µL for sample concentrations between 0.30 and 6.0 pmol/µL; ±2% for samples >6.0 pmol/µL

<sup>a</sup> Values are approximate

<sup>b</sup> Based on five replicates (SD=ng/ $\mu$ L; CV=%)

4

# **Nucleic Acid Applications**

# Measure dsDNA, ssDNA or RNA

Measures the concentration of purified dsDNA, ssDNA or RNA samples that absorb at 260 nm.

Measure dsDNA, ssDNA or RNA

**Reported Results** 

Settings

**Detection Limits** 

Calculations



### Measure dsDNA, ssDNA or RNA

Use the dsDNA, ssDNA and RNA applications to quantify purified double-stranded (ds) or RNA samples. These applications report nucleic acid concentration and two absorbance ratios (A260/A280 and A260/A230). A single-point baseline correction can also be used.

### To measure dsDNA, ssDNA or RNA samples

### NOTICE

- Do not use a squirt or spray bottle on or near the instrument as liquids will flow into the instrument and may cause permanent damage.
- Do not use hydrofluoric acid (HF) on the pedestals. Fluoride ions will permanently damage the quartz fiber optic cables.

### Before you begin...

Before taking pedestal measurements with the NanoDrop Eight instrument, lift the instrument arm and clean the upper and lower pedestals. At a minimum, wipe the pedestals with a new laboratory wipe. For more information, see Cleaning the Pedestals.

### To measure nucleic acid


- 1. From the Home screen, select the **Nucleic Acids** tab, select your sample mode, and select **dsDNA**, **ssDNA** or **RNA**, depending on the samples to be measured.
- 2. Specify a baseline correction if desired.
- 3. Pipette 1–2 µL blanking solution onto the lower pedestal and lower the arm.
- 4. Select **Blank** and wait for the measurement to complete.

**Tip**: If Auto-Blank is On, the blank measurement starts automatically after you lower the arm.


- 5. Lift the arm and clean all pedestals with a new laboratory wipe.
- 6. Pipette 1-2 µL sample solution onto the pedestal and lower the arm.
- 7. Start the sample measurement: If Auto-Measure is On, lower arm; if Auto-Measure is off, lower arm and select **Measure**.

When the sample measurement is completed, the spectrum and reported values are displayed (see the next section).

- 8. When you are finished measuring samples, select End Experiment
- 9. Lift the arm and clean all pedestals with a new wipe.

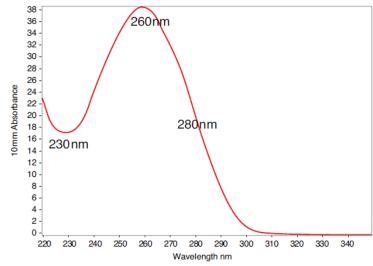

**Typical nucleic acid spectrum** 

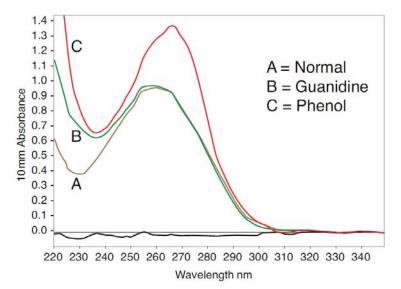



### Best practices for nucleic acid measurements

 Isolate and purify nucleic acid samples before measurement to remove impurities. Depending on the sample, impurities could include DNA, RNA, free nucleotides, proteins, some buffer components and dyes. See Preparing Samples for more information. **Note** Extraction reagents such as guanidine, phenol, and EDTA contribute absorbance between 230 nm and 280 nm and will affect measurement results if present in samples (even residual amounts).

- Ensure the sample absorbance is within the instrument's absorbance detection limits.
- Blank with the same buffer solution used to resuspend the analyte of interest. The blanking solution should be a similar pH and ionic strength as the analyte solution.
- Run a blanking cycle to assess the absorbance contribution of your buffer solution. If the buffer exhibits strong absorbance at or near the analysis wavelength (typically 260 nm), you may need to choose a different buffer or application. See Choosing and Measuring a Blank for more information.
- For micro-volume measurements:
  - Ensure pedestal surfaces are properly cleaned and conditioned.
  - If possible, heat highly concentrated or large molecule samples, such as genomic or lambda DNA, to 63 °C (145 °F) and gently (but thoroughly) vortex before taking a measurement. Avoid introducing bubbles when mixing and pipetting.
  - Follow best practices for micro-volume measurements.
  - Use a 1-2 µL sample volume. See Recommended Sample Volumes for more information.


### **Related Topics**

- Measure a Micro-Volume Sample
- Best Practices for Micro-Volume Measurements
- Prepare Samples and Blanks
- Basic Instrument Operations

### **Nucleic Acid Reported Results**

### dsDNA measurement screen

For each measured sample, the dsDNA, ssDNA and RNA applications show the UV absorbance spectrum and a summary of the results. Below is an example of the measurement screen of the instrument software:

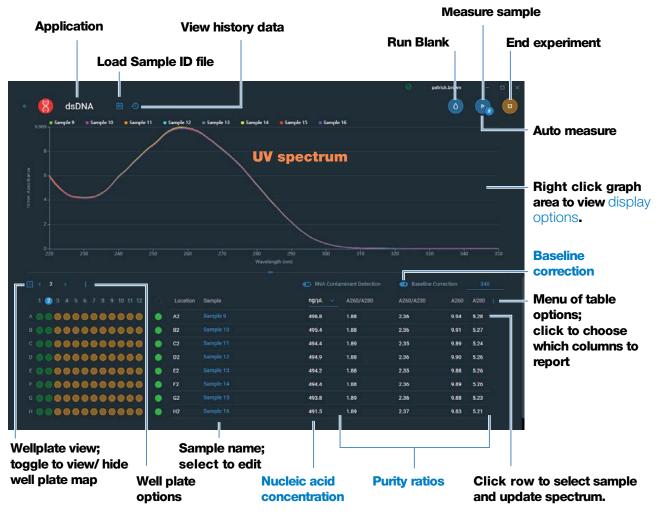

**Note** Micro-volume absorbance measurements are normalized to a 10.0 mm pathlength equivalent.

### dsDNA, ssDNA and RNA reported values

- sample name
- created on (date sample measurement was taken)
- nucleic acid concentration
- Species
- A260/A280
- A260/A230
- A260
- A280
- factor
- baseline correction
- monitor wavelength
- pathlength used
- contaminant
- corrected

### **Settings for Nucleic Acid Measurements**

### **Baseline Correction**

For the dsDNA, ssDNA or RNA, enter baseline correction wavelength in nm or use default value (340 nm)

**Optional user-defined baseline correction**. Can be used to correct for any offset caused by light scattering particulates by subtracting measured absorbance at specified baseline correction wavelength from absorbance values at all wavelengths in sample spectrum. As a result, absorbance of sample spectrum is zero at specified baseline correction wavelength.

# **Calculations for Nucleic Acid Measurements**

The Nucleic Acid applications use a modification of the Beer-Lambert equation (shown at right) to calculate sample concentration where the extinction coefficient and pathlength are combined and referred to as a "factor."

### **Extinction Coefficients vs Factors**

Using the terms in the Beer-Lambert equation, factor (f) is defined as:

factor (f) = 
$$1/(\varepsilon * b)$$

where:

 $\epsilon$  = wavelength-dependent molar extinction coefficient in ng-cm/µL

### b = sample pathlength in cm

As a result, analyte concentration (c) is calculated as:

$$c = A * [1/(\varepsilon * b)]$$

or

where:

c = analyte concentration in ng/ $\mu$ L

A = absorbance in absorbance units (A)

 $\mathbf{f} = \text{factor in ng-cm/}\mu\text{L}$  (see below)

### **Factors Used**

- **dsDNA** (factor = 50 ng-cm/µL)
- ssDNA (factor = 33 ng-cm/µL)
- **RNA** (factor = 40 ng-cm/ $\mu$ L)
- **Custom Factor** (user entered factor between 15 ng-cm/µL and 150 ng-cm/µL

For the dsDNA, ssDNA and RNA applications, the generally accepted factors for nucleic acids are used in conjunction with Beer's Law to calculate sample concentration. For the Custom Factor application, the user-specified factor is used.

#### **4** Nucleic Acid Applications Measure dsDNA, ssDNA or RNA

Calculated nucleic acid concentrations are based on the absorbance value at 260 nm, the factor used and the sample pathlength. A single-point baseline correction (or analysis correction) may also be applied.

Concentration is reported in mass units. Calculators are available on the Internet to convert concentration from mass to molar units based on sample sequence.

Absorbance values at 260 nm, 280 nm and sometimes 230 nm are used to calculate purity ratios for the measured nucleic acid samples. Purity ratios are sensitive to the presence of contaminants in the sample, such as residual solvents and reagents typically used during sample purification.

### **Measured Values**

**Note**: For micro-volume absorbance measurements the spectra are normalized to a 10 mm pathlength equivalent.

### A260 absorbance

- Nucleic acid absorbance values are measured at 260 nm using the normalized spectrum. This is the reported A260 value if Baseline Correction is not selected.
- If Baseline Correction is selected, the absorbance value at the correction wavelength is subtracted from the absorbance at 260 nm. The corrected absorbance at 260 nm is reported and used to calculate nucleic acid concentration.

### A230 and A280 absorbance

• Normalized and baseline-corrected (if selected) absorbance values at 230 nm and 280 nm are used to calculate A260/A230 and A260/A280 ratios.

### **Sample Pathlength**

- For micro-volume measurements, the software selects the optimal pathlength (between 1.0 mm and 0.1 mm) based on sample absorbance at the analysis wavelength.
- Displayed spectra and absorbance values are normalized to a 10 mm pathlength equivalent.

### **Reported Values**

- Nucleic acid concentration. Reported in selected unit (i.e., ng/µL, µg/µl or µg/mL). Calculations are based on modified Beer's Law equation using corrected nucleic acid absorbance value.
- A260/A280 purity ratio. Ratio of corrected absorbance at 260 nm to corrected absorbance at 280 nm. An A260/A280 purity ratio of ~1.8 is generally accepted as "pure" for DNA (~2.0 for RNA). Acidic solutions may under represent the reported value by 0.2-0.3; the opposite is true for basic solutions.
- **A260/A230 purity ratio**. Ratio of corrected absorbance at 260 nm to corrected absorbance at 230 nm. An A260/A230 purity ratio between 1.8 and 2.2 is generally accepted as "pure" for DNA and RNA.

**Note**: Although purity ratios are important indicators of sample quality, the best quality indicator quality is functionality in the downstream application of interest (e.g., real-time PCR).

- **Factor**. Used in conjunction with Beer's Law to calculate sample concentration
- **Contaminant** If a contaminant was identified by the Acclaro software, the contaminant will be displayed in this column.
- **Species** The species of nucleic acid contaminant detected by the Acclaro sample intelligence technology
- A260 absorbance.
- A280 absorbance.
- Baseline correction
- **Monitor wavelength** Enter an additional wavelength whose absorbance value you want included in the report.
- **Corrected** Displays the corrected analyte concentration determined using the Acclaro software, if one is available.

## **Measure Microarray**

Measures the concentration of purified nucleic acids that have been labeled with up to two fluorescent dyes for use in downstream microarray applications.

Measure Microarray Samples

**Reported Results** 

Settings

**Detection Limits** 

Calculations



### **Measure Microarray Samples**

Use the Microarray application to quantify nucleic acids that have been labeled with up to two fluorescent dyes. The application reports nucleic acid concentration, an A260/A280 ratio and the concentrations and measured absorbance values of the dye(s), allowing detection of dye concentrations as low as 0.2 picomole per microliter.

### To measure microarray samples

### NOTICE

- Do not use a squirt or spray bottle on or near the instrument as liquids will flow into the instrument and may cause permanent damage.
- Do not use hydrofluoric acid (HF) on the pedestals. Fluoride ions will permanently damage the quartz fiber optic cables.

### Before you begin...

Before taking pedestal measurements with the NanoDrop Eight instrument, lift the instrument arm and clean the upper and lower pedestals. At a minimum, wipe the pedestals with a new laboratory wipe. For more information, see Cleaning the Pedestals.

### To measure a microarray sample

- 1. From the Home screen, select the Nucleic Acids tab and select Microarray.
- 2. Specify the sample type and factor and the type of dye(s) used.

**Tip**: Select a dye from the pre-defined list or add a custom dye using the Dye/Chromophore Editor.


- 3. Pipette  $1-2 \mu$ L blanking solution onto the lower pedestal and lower the arm.
- 4. Select **Blank** and wait for the measurement to complete.

**Tip**: If Auto-Blank is On, the blank measurement starts automatically after you lower the arm.

- 5. Lift the arm and clean all pedestals with a new laboratory wipe.
- 6. Pipette 1-2 µL sample solution onto the pedestal and lower the arm.
- 7. Start the sample measurement: If Auto-Measure is On, lower arm; if Auto-Measure is off, lower arm and select **Measure**.

When the sample measurement is completed, the spectrum and reported values are displayed (see the next section).

- 8. When you are finished measuring samples, select **End Experiment**.
- 9. Lift the arm and clean all pedestals with a new wipe.

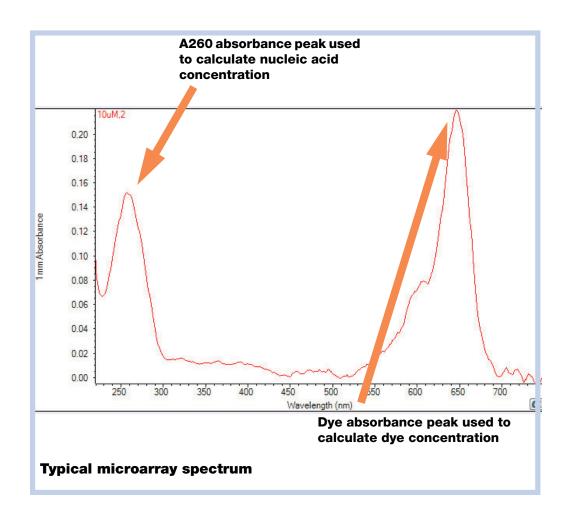


### **Related Topics**

- Best Practices for Nucleic Acid Measurements
- Measure a Micro-Volume Sample
- Best Practices for Micro-Volume Measurements
- Prepare Samples and Blanks
- Basic Instrument Operations

### **Microarray Reported Results**

### **Microarray measurement screen**

For each measured sample, this application shows the absorbance spectrum and a summary of the results. Here is an example:

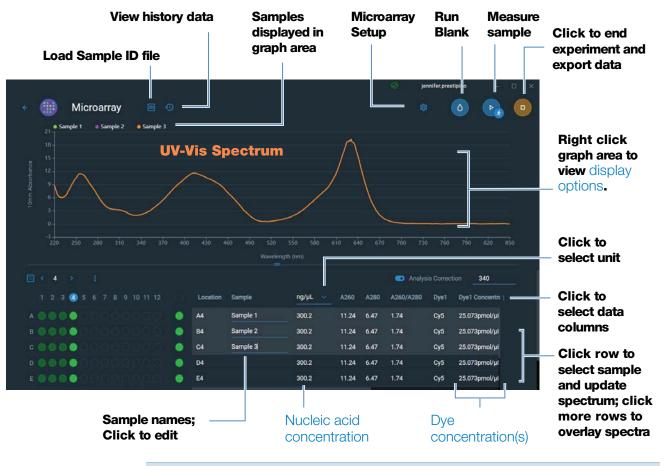

### Note

- Analysis correction is performed at 340 nm (absorbance value at 340 nm is subtracted from absorbance values at all wavelengths in sample spectrum).
- Micro-volume absorbance measurements are normalized to a 10.0 mm pathlength equivalent.

### **Microarray reported values**

The initial screen that appears after each measurement (see previous image) shows a summary of the reported values. To view all reported values, press and hold the sample row. Here is an example:

- sample name
- created on (date sample measurement was taken)
- nucleic acid concentration
- A260
- A260/A280
- dye 1/dye 2 concentration
- sample type
- analysis correction
- factor

### **Settings for Microarray Measurements**

#### **Microarray settings**

The Microarray Setup screen appears after you select the Microarray application from the Nucleic Acids tab on the Home screen. To show the Microarry settings from the Microarray measurement screen, select Microarray Setup

Setting	Available Options	Description
Sample type and Factor	dsDNA (with non-editable factor of 50 ng-cm/µL)	Widely accepted value for double-stranded DNA
	ssDNA (with non-editable factor of 33 ng-cm/µL)	Widely accepted value for single-stranded DNA
	RNA (with non-editable factor of 40 ng-cm/µL)	Widely accepted value for RNA
	Oligo DNA with non-editable calculated factor in ng-cm/µL	Factor calculated from user-defined DNA base sequence. When selected, available DNA base units (i.e., G, A, T, C) appear as keys. Define sequence by selecting appropriate keys. Factor is calculated automatically based on widely accepted value for each base unit.

Setting	Available Options	Description
	Oligo RNA with non-editable calculated factor in ng-cm/µL	Factor calculated from user-defined RNA base sequence. When selected, available RNA base units (i.e., G, A, U, C) appear as keys. Define sequence by selecting appropriate keys. Factor is calculated automatically based on widely accepted value for each base unit.
	Custom (with user-specified factor in ng-cm/ $\mu$ L)	Enter factor between 15 ng-cm/µL and 150 ng-cm/µL
Dye 1/Dye 2 Type <sup>a</sup>	Cy3, 5, 3.5, or 5.5, Alexa Fluor 488, 546, 555, 594, 647, or 660	Select pre-defined dye(s) used to label sample material, or one that has been added using Dye Editor.
Dye 1/Dye 2 Unit	picomoles/microliter (pmol/µl), micromoles (uM), or millimoles (mM)	Select unit for reporting dye concentrations
Analysis Correction <sup>b</sup>	On or off Enter analysis correction wavelength in nm or use default value (340 nm)	Corrects sample absorbance measurement for any offset caused by light scattering particulates by subtracting absorbance value at specified analysis correction wavelength from absorbance value at analysis wavelength. Corrected value is used to calculate sample concentration.
		<b>Tip</b> : If the sample has a modification that absorbs light at 340 nm, select a different correction wavelength or turn off Analysis Correction.

<sup>a</sup> To add a custom dye or edit the list of available dyes, use the Dye/Chromophore Editor.


<sup>b</sup> The Analysis Correction affects the calculation for nucleic acid concentration only.

### **Dye/chromophore editor**

Use the Dye/Chromophore Editor to add a custom dye to the list of available dyes in Microarray Setup or Proteins & Labels Setup. You can also specify which dyes are available in that list.

To access the Dye/Chromophore Editor, from the Home screen, select **Settings > Dye Editor**.

#### **Dye Editor**

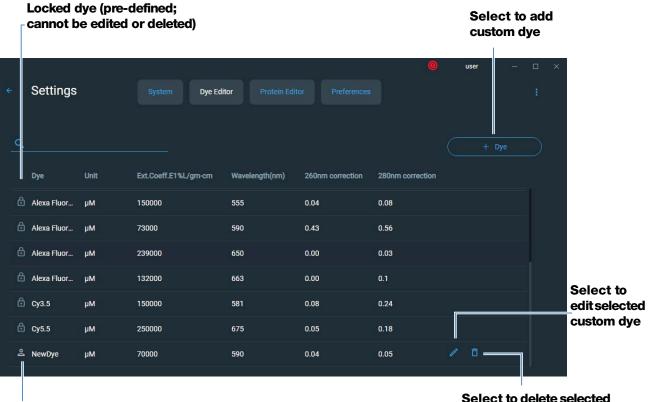

# Custom dye (user-defined; can be edited or deleted)



These operations are available from the Dye/Chromophore Editor:

#### Add custom dye

- tap + Dye to show Create Dye box
- enter unique Name for new dye
- select default Unit that will be used to display dye concentration
- enter dye's Extinction Coefficient (or molar absorptivity constant) in L/mole-cm (typically provided by dye manufacturer)
- specify Wavelength in nm (between 350 nm and 850 nm) that will be used to measure dye's absorbance
- specify dye's correction values at 260 nm and 280 nm
- select Save

**Note** To determine dye correction values (if not available from dye manufacturer):

- use instrument to measure pure dye and note absorbance at 260 nm,
   280 nm and at analysis wavelength for dye (see above)
- calculate ratio of A<sub>260</sub>/A<sub>dye wavelength</sub> and enter that value for 260 nm Correction
- calculate ratio of  $A_{280}/A_{dye\ wavelength}$  and enter that value for 280 nm Correction

When a custom dye is selected before a measurement, the dye's absorbance and concentration values are reported and the corrections are applied to the measured sample absorbance values, and to the resulting sample concentrations and purity ratios.

### Edit custom dye

**Tip** Dyes pre-defined in the software cannot be edited.

- select custom dye
- select edit 🧪
- edit any entries or settings
- click Save

#### **Delete custom dye**

**Tip** Dyes pre-defined in the software cannot be deleted.

- select custom dye
- click 🛅

**NOTICE** Deleting a custom dye permanently removes the dye and all associated information from the software.

### **Calculations for Microarray Measurements**

As with the other nucleic acid applications, the Microarray application uses a modification of the Beer-Lambert equation to calculate sample concentration where the extinction coefficient and pathlength are combined and referred to as a "factor." The Microarray application offers six options (shown at right) for selecting an appropriate factor for each measured sample, to be used in conjunction with Beer's Law to calculate sample concentration.

If the factor is known, choose the Custom Factor option and enter the factor in ng-cm/ $\mu$ L. Otherwise, choose the option that best matches the sample solution.

**Tip**: Ideally, the factor or extinction coefficient should be determined empirically using a solution of the study nucleic acid at a known concentration using the same buffer.

### **Available Options for Factors**

- **dsDNA** (factor = 50 ng-cm/µL)
- ssDNA (factor = 33 ng-cm/µL)
- **RNA** (factor = 40 ng-cm/µL)
- Oligo DNA (calculated from user entered DNA nucleotide sequence)
- Oligo RNA (calculated from user entered RNA nucleotide sequence)
- **Custom Factor** (user entered factor between 15 ng-cm/µL and 150 ng-cm/µL

Note: See Sample Type for more information.

Calculated nucleic acid concentrations are based on the absorbance value at 260 nm, the factor used and the sample pathlength. A single-point baseline correction (or analysis correction) may also be applied.

Concentration is reported in mass units. Calculators are available on the Internet to convert concentration from mass to molar units based on sample sequence.

Absorbance values at 260 nm, 280 nm and sometimes 230 nm are used to calculate purity ratios for the measured nucleic acid samples. Purity ratios are sensitive to the presence of contaminants in the sample, such as residual solvents and reagents typically used during sample purification.

### **Measured Values**

### A260 absorbance

**Note**: The absorbance value at 850 nm is subtracted from all wavelengths in the spectrum. As a result, the absorbance at 850 nm is zero in the displayed spectra. Also, for micro-volume absorbance measurements, the spectra are normalized to a 10 mm pathlength equivalent.

- Nucleic acid absorbance values for all Microarray sample types are measured at 260 nm using the 850-corrected and normalized spectrum.
- If Analysis Correction is selected, the absorbance value at the correction wavelength is subtracted from the absorbance at 260 nm.
- If one or more dyes are selected, the dye correction values at 260 nm are also subtracted from the absorbance at 260 nm.
- The final corrected absorbance at 260 nm is reported and used to calculate sample concentration.

### A280 absorbance

• Analysis-corrected and normalized absorbance value at 280 nm (minus the A280 dye correction) is used to calculate an A260/A280 ratio.

Dye concentrations are calculated from the absorbance value at the dye's analysis wavelength, the dye's extinction coefficient, and the sample pathlength. A sloped-line dye correction may also be used.

### **Dye absorbance**

- Dye absorbance values are measured at specific wavelengths. See Dye/Chromophore Editor for analysis wavelengths used.
- If Sloping Dye Correction is selected, a linear baseline is drawn between 400 nm and 850 nm and, for each dye, the absorbance value of the sloping baseline is subtracted from the absorbance value at each dye's analysis wavelength. Baseline-corrected dye absorbance values are reported and used to calculate dye concentrations.

### **Dye correction**

- Pre-defined dyes have known correction values for A260 and A280. See Dye/Chromophore Editor for correction values used.
- A260 dye corrections are subtracted from the A260 absorbance value used to calculate nucleic acid concentration, and from the A260 absorbance value used to calculate the A260/A280 purity ratio.

### **Sample Pathlength**

- For micro-volume measurements, the software selects the optimal pathlength (between 1.0 mm and 0.1 mm) based on sample absorbance at the analysis wavelength.
- Displayed spectra and absorbance values are normalized to a 10 mm pathlength equivalent.

### **Reported Values**

- Nucleic acid concentration. Reported in selected unit (i.e., ng/µL, µg/µl or µg/mL). Calculations are based on modified Beer's Law equation using corrected nucleic acid absorbance value.
- **A260/A280 purity ratio**. Ratio of corrected absorbance at 260 nm to corrected absorbance at 280 nm. An A260/A280 purity ratio of ~1.8 is generally accepted as "pure" for DNA (~2.0 for RNA). Acidic solutions may under represent the reported value by 0.2-0.3; the opposite is true for basic solutions.
- **Dye1/Dye2 concentration**. Reported in pmol/µL. Calculations are based on Beer's Law equation using (sloping) baseline-corrected dye absorbance value(s).

**Note**: Although purity ratios are important indicators of sample quality, the best indicator of DNA or RNA quality is functionality in the downstream application of interest (e.g., microarray).

### **Related Topics**

• Calculations for Nucleic Acid Measurements

#### **4** Nucleic Acid Applications Measure using a Custom Factor

## Measure using a Custom Factor

Measures the concentration of purified nucleic acids using a custom factor for the calculations.

Measure using Custom Factor

**Reported Results** 

Settings

**Detection Limits** 

Calculations



### **Measure Nucleic Acid using a Custom Factor**

Use the Custom Factor application to quantify purified DNA or RNA samples that absorb at 260 nm with a user-defined extinction coefficient or factor. The application reports nucleic acid concentration and two absorbance ratios (A260/A280 and A260/A230). A single-point baseline correction can also be used.

### To measure nucleic acid samples using a custom factor

### NOTICE

- Do not use a squirt or spray bottle on or near the instrument as liquids will flow into the instrument and may cause permanent damage.
- Do not use hydrofluoric acid (HF) on the pedestals. Fluoride ions will permanently damage the quartz fiber optic cables.

#### Before you begin...

Before taking pedestal measurements with the NanoDrop Eight instrument, lift the instrument arm and clean the upper and lower pedestals. At a minimum, wipe the pedestals with a new laboratory wipe. For more information, see Cleaning the Pedestals.

### To measure using a custom factor

1. From the Home screen, select the Nucleic Acids tab and select Custom Factor.

- Custom Factor Setup Factor 50 Done
- Enter the factor to be used for the calculations and specify a baseline correction if desired.


- 3. Pipette 1–2 µL blanking solution onto the lower pedestals and lower the arm.
- 4. Select **Blank** and wait for the measurement to complete.

**Tip**: If Auto-Blank is On, the blank measurement starts automatically after you lower the arm.

- 5. Lift the arm and clean all pedestals with a new laboratory wipe.
- 6. Pipette 1-2 µL sample solution onto the pedestals and lower the arm.
- 7. Start the sample measurement: If Auto-Measure is On, lower arm; if Auto-Measure is off, lower arm and select **Measure**.

When the sample measurement is completed, the spectrum and reported values are displayed (see the next section).

- 8. When you are finished measuring samples, select
- End Experiment.
- 9. Lift the arm and clean all pedestals with a new wipe.

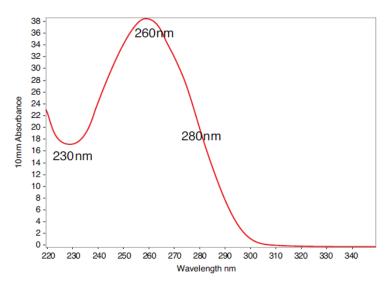

**Typical nucleic acid spectrum** 


#### 4 Nucleic Acid Applications Measure using a Custom Factor

#### **Related Topics**

- Measure a Micro-Volume Sample
- Best Practices for Micro-Volume Measurements
- Prepare Samples and Blanks
- Basic Instrument Operations

## **Custom Factor Reported Results**

For each measured sample, this application shows the absorbance spectrum and a summary of the results. Here is an example:

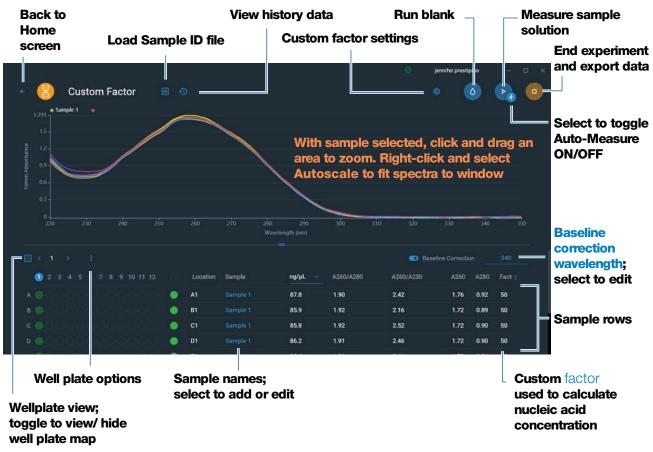

#### Tips:

Click sample row to select sample and update spectrum Shift-click multiple sample rows to overlay spectra Click a sample and hover locations on spectra to view measurement values

#### **Related Topics**

- Basic Instrument Operations
- Nucleic Acid Reported Results
- Nucleic Acid Calculations

## **Settings for Nucleic Acid Measurements using a Custom Factor**

To show the Custom Factor settings, from the Custom Factor measurement screen, select the settings icon to view the **Custom Factor Setup**.

Setting	Available Options	Description
Custom Factor	Enter an integer value between 15 ng-cm/µL and 150 ng-cm/µL	Constant used to calculate nucleic acid concentration in modified Beer's Law equation. Based on extinction coefficient and pathlength:
		$f = 1/(E_{260} * b))$
		where: f = factor E = molar extinction coefficient at 260 nm in ng-cm/µL $b = sample pathlength$ in cm (1 cm for nucleic acids measured with the NanoDrop Eight instruments)
Baseline Correction	On or off Enter baseline correction wavelength in nm or use default value (340 nm)	<b>Optional user-defined baseline correction</b> . Can be used to correct for any offset caused by light scattering particulates by subtracting measured absorbance at specified baseline correction wavelength from absorbance values at all wavelengths in sample spectrum. As a result, absorbance of sample spectrum is zero at specified baseline correction wavelength.
		NOTE: Baseline correction is selected from the measurement screen of the instrument software and is not shown in the Custom Factor Setup.

#### **Related Topics**

• Instrument Settings

# Detection Limits for Nucleic Acid Measurements using a Custom Factor

The lower detection limits and reproducibility specifications for nucleic acids are provided here. The upper detection limits are dependent on the upper absorbance limit of the instrument and the user-defined extinction coefficients.

#### To calculate upper detection limits for nucleic acid samples

To calculate upper detection limits in  $ng/\mu L$ , use the following equation:

(upper absorbance limit<sub>instrument</sub> \* extinction coefficient<sub>sample</sub>)

For example, for a sample measurement using an extinction coefficient of 31, the equation looks like this:

(200 AU \* 31 ng-cm/µL) = 6,200 ng/µL

#### **Related Topics**

• Detection Limits for All Applications

#### 4 Nucleic Acid Applications Measure Oligo DNA or Oligo RNA

## Measure Oligo DNA or Oligo RNA

Measures the concentration of purified ssDNA or RNA oligonucleotides that absorb at 260 nm.


Measure Oligo DNA or RNA

**Reported Results** 

Settings

**Detection Limits** 

Calculations

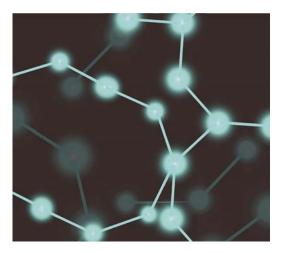

## **Measure Oligo DNA or Oligo RNA**

Use the Oligo DNA and Oligo RNA applications to quantify oligonucleotides that absorb at 260 nm. Molar extinction coefficients are calculated automatically based on the user-defined base sequence of the sample. These applications report nucleic acid concentration and two absorbance ratios (A260/A280 and A260/A230). A single-point baseline correction can also be used.

#### To measure Oligo DNA or Oligo RNA samples

#### NOTICE

- Do not use a squirt or spray bottle on or near the instrument as liquids will flow into the instrument and may cause permanent damage.
- Do not use hydrofluoric acid (HF) on the pedestals. Fluoride ions will permanently damage the quartz fiber optic cables.

#### Before you begin...

Before taking pedestal measurements with the NanoDrop Eight instrument, lift the instrument arm and clean the upper and lower pedestals. At a minimum, wipe the pedestals with a new laboratory wipe. For more information, see Cleaning the Pedestals.

#### To measure an oligonucleotide sample

- 1. From the Home screen, select the **Nucleic Acids** tab and select either **Oligo DNA** or **Oligo RNA**, as needed.
- 2. Specify the Oligo base sequence and a baseline correction if desired.


- 3. Pipette 1–2 µL blanking solution onto the lower pedestal and lower the arm.
- 4. Select **Blank** and wait for the measurement to complete.

**Tip**: If Auto-Blank is On, the blank measurement starts automatically after you lower the arm.

- 5. Lift the arm and clean all pedestals with a new laboratory wipe.
- 6. Pipette 1-2 µL sample solution onto the pedestal and lower the arm.
- 7. Start the sample measurement: If Auto-Measure is On, lower arm; if Auto-Measure is off, lower arm and select **Measure**.

When the sample measurement is completed, the spectrum and reported values are displayed (see the next section).

- 8. When you are finished measuring samples, tap **End Experiment**
- 9. Lift the arm and clean all pedestals with a new wipe.

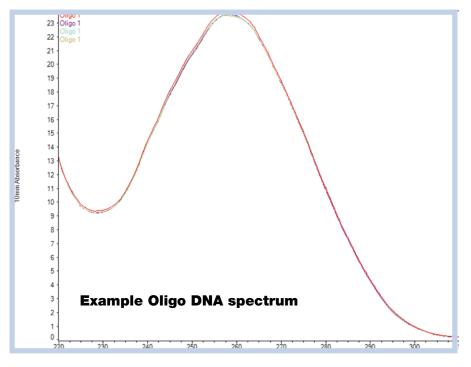

#### **Related Topics**


- Best Practices for Nucleic Acid Measurements
- Measure a Micro-Volume Sample
- Best Practices for Micro-Volume Measurements

- Prepare Samples and Blanks
- Basic Instrument Operations

## **Oligo Reported Results**

#### **Oligo DNA and RNA measurement screen**

For each measured sample, the Oligo DNA and Oligo RNA applications show the UV absorbance spectrum reported results. Here is an example:

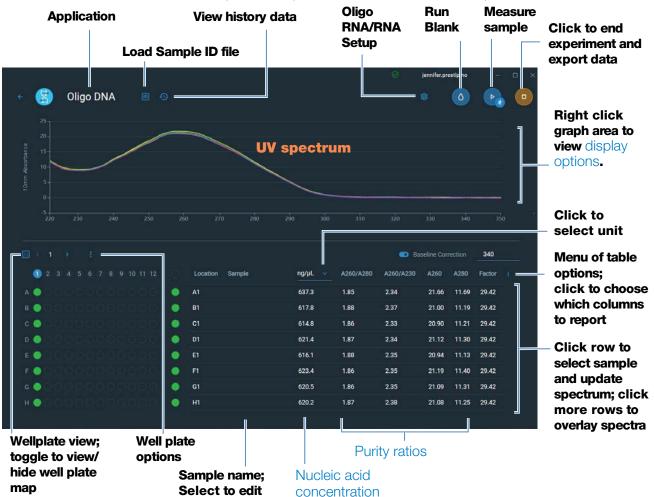

#### Oligo DNA and Oligo RNA reported values

• sample name

#### 4 Nucleic Acid Applications Measure Oligo DNA or Oligo RNA

- created on (date sample measurement was taken)
- nucleic acid concentration
- A260/A280
- A260/A230
- A260
- A280
- factor
- oligo sequence
- baseline correction

**Note** The five nucleotides that comprise DNA and RNA exhibit widely varying A260/A280 ratios. See Oligo Purity Ratios for more information.

#### **Related Topics**

- Basic Instrument Operations
- Oligo Calculations

## **Settings for Oligo DNA and Oligo RNA Measurements**

The Oligo setup screen appears after you select the Oligo DNA or Oligo RNA application from the Nucleic Acids tab on the Home screen. From the Oligo DNA or RNA measurement screen, select the settings icon to view **Oligo DNA Setup** or **Oligo RNA Setup**.

Setting	Available Options	Description			
Oligo Base Sequence for DNA: Use the G, A, T and C keys to specify the DNA base sequence	G, A, T and C keys to	Specify your DNA or RNA base sequence. click the corresponding keys:			
	Add guanine Add adenine Add thymine (DNA) or uracil (RNA)				
	for RNA: Use the G, A, U and C keys to specify the RNA base sequence	Cligo DNA Setup GAT CTA GAT CTA Factor Factor Vumber of bases Cancel Cancel Done Base sequence			
		You can also enter base sequence using the keyboard, or by copy and pasting a sequence from another application.			
		Each time a base is added to the sequence, the software			

• **Factor**. Constant used to calculate oligonucleotide concentration in modified Beer's Law equation. Based on extinction coefficient and pathlength:

$$f = 1/(\mathcal{E}_{260} * b)$$

calculates the following:

where:

- f= factor
- $\mathcal{E}$  = molar extinction coefficient at 260 nm in ng-cm/µL b = sample pathlength in cm (1 cm for nucleic acids measured with the NanoDrop Eight instrument)

#### 4 Nucleic Acid Applications Measure Oligo DNA or Oligo RNA

Setting	Available Options	Description
		• <b>Molecular Weight</b> of oligo calculated from user-defined base sequence.
		Number of Bases entered.
		• Molar Ext. Coefficient (260 nm). Molar extinction coefficient of oligo (in ng-cm/µL) at 260 nm calculated from entered base sequence.
		<ul> <li>%GC. Percentage of guanine and cytosine residues in total number of bases entered.</li> </ul>
Baseline Correction	On or off Enter baseline correction wavelength in nm or use default value (340 nm)	Corrects for any offset caused by light scattering particulates by subtracting measured absorbance at specified baseline correction wavelength from absorbance values at all wavelengths in sample spectrum. As a result, absorbance of sample spectrum is zero at specified baseline correction wavelength. <b>Tip</b> : If the sample has a modification that absorbs light at 340 nm, select a different correction wavelength or turn off Baseline Correction.

#### **Related Topics**

• Instrument Settings

## **Detection Limits for Oligo DNA and Oligo RNA Measurements**

The lower detection limits and reproducibility specifications for the oligonucleotide sample types (ssDNA and RNA) are provided here. The upper detection limits are dependent on the upper absorbance limit of the instrument and the extinction coefficients for the user-defined base sequences.

#### To calculate upper detection limits for nucleic acid samples

To calculate upper detection limits in  $ng/\mu L$ , use the following equation:

(upper absorbance limit<sub>instrument</sub> \* extinction coefficient<sub>sample</sub>)

For example, for a sample measurement using an extinction coefficient of 55, the equation looks like this:

 $(200 \text{ AU} * 55 \text{ ng-cm/}\mu\text{L}) = 11,000 \text{ ng/}\mu\text{L}$ 

## Calculations for Oligo DNA and Oligo RNA Measurements

As with the other nucleic acid applications, the Oligo applications use the Beer-Lambert equation to correlate absorbance with concentration based on the sample's extinction coefficient and pathlength. Because oligonucleotides are short, single-stranded molecules (or longer molecules of repeating sequences), their spectrum and extinction coefficient ( $\mathbf{E}$ ) are closely dependent on base composition and sequence.

(The generally accepted extinction coefficients and factors for single-stranded DNA and RNA provide a reasonable estimate for natural, essentially randomized, sequences but not for short, synthetic oligo sequences.) To ensure the most accurate results, we use the exact value of  $\mathcal{E}_{260}$  to calculate oligonucleotide concentration.

The NanoDrop software allows you to specify the base sequence of an oligonucleotide before it is measured. For any entered base sequence, the software uses the equation at the right to calculate the extinction coefficient.

**Tip**: The extinction coefficient is wavelength specific for each oligonucleotide and can be affected by buffer type, ionic strength and pH.

#### **Extinction Coefficients for Oligonucleotides**

The software uses the nearest neighbor method and the following formula to calculate molar extinction coefficients for specific oligonucleotide base sequences:

$$\varepsilon_{260} = \sum_{1}^{N-1} \varepsilon_1 - \sum_{2}^{N-1} \varepsilon_2 + \sum_{1}^{N} \varepsilon_3$$

where:

 $\mathcal{E}$  = molar extinction coefficient in L/mole-cm

 $\mathbf{\hat{E}}_1 = \mathbf{\hat{E}}_{nearest neighbor}$   $\mathbf{\hat{E}}_2 = \mathbf{\hat{E}}_{individual bases}$   $\mathbf{\hat{E}}_3 = \mathbf{\hat{E}}_{modifications, such as fluorescent dyes}$ 

Calculated nucleic acid concentrations are based on the absorbance value at 260 nm, the factor used and the sample pathlength. A single-point baseline correction (or analysis correction) may also be applied.

Concentration is reported in mass units. Calculators are available on the Internet to convert concentration from mass to molar units based on sample sequence.

Absorbance values at 260 nm, 280 nm and sometimes 230 nm are used to calculate purity ratios for the measured nucleic acid samples. Purity ratios are sensitive to the presence of contaminants in the sample, such as residual solvents and reagents typically used during sample purification.

#### **Measured Values**

#### A260 absorbance

**Note**: For micro-volume absorbance measurements, the spectra are normalized to a 10 mm pathlength equivalent.

- Nucleic acid absorbance values are measured at 260 nm using the normalized spectrum. This is the reported A260 value if Baseline Correction is not selected.
- If Baseline Correction is selected, the absorbance value at the correction wavelength is subtracted from the sample absorbance at 260 nm. The corrected absorbance at 260 nm is reported and used to calculate nucleic acid concentration.

#### A230, A280 absorbance

• Normalized absorbance values at 230 nm, 260 nm and 280 nm are used to calculate A260/A230 and A260/A280 ratios.

#### **Sample Pathlength**

- For micro-volume measurements, the software selects the optimal pathlength (between 1.0 mm and 0.1 mm) based on sample absorbance at the analysis wavelength.
- Displayed spectra and absorbance values are normalized to a 10 mm pathlength equivalent.

The five nucleotides that comprise DNA and RNA exhibit widely varying A260/A280 ratios. Estimated A260/A280 ratios for each independently measured nucleotide are provided below:

Guanine: 1.15 Adenine: 4.50 Cytosine: 1.51 Uracil: 4.00 Thymine: 1.47

The A260/A280 ratio for a specific nucleic acid sequence is approximately equal to the weighted average of the A260/A280 ratios for the four nucleotides present.

**Note**: RNA will typically have a higher 260/280 ratio due to the higher ratio of Uracil compared to that of Thymine.

#### **Reported Values**

- Nucleic acid concentration. Reported in selected unit (i.e., ng/µL, µg/µl or µg/mL). Calculations are based on modified Beer's Law equation using corrected nucleic acid absorbance value.
- **A260/A280 purity ratio**. Ratio of corrected absorbance at 260 nm to corrected absorbance at 280 nm.
- **A260/A230 purity ratio**. Ratio of corrected absorbance at 260 nm to corrected absorbance at 230 nm.

**Note**: The traditional purity ratios (A260/A280 and A260/A230), which are used as indicators of the presence of various contaminants in nucleic acid samples, do not apply for oligonucleotides because the shapes of their spectra are highly dependent on their base compositions. See side bar for more information.

## **Related Topics**

Calculations for Nucleic Acid Measurements

5

## **Protein Applications**

## **Measure Protein A280**

Measures the concentration of purified protein samples that absorb at 280 nm.


Measure A280 Proteins

Reported Results

Settings

**Detection Limits** 

Calculations

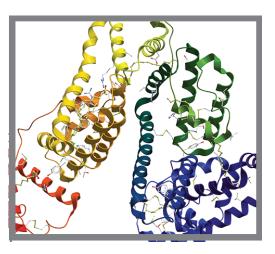

## **Measure Protein Concentration at A280**

Use the Protein A280 application to quantify purified protein samples that contain amino acids such as tryptophan or tyrosine, or cys-cys disulfide bonds, which exhibit absorbance at 280 nm. This application reports protein concentration measured at 280 nm and one absorbance ratio (A260/A280). A single-point baseline correction can also be used. This application does not require a standard curve.

#### To measure Protein A280 samples

#### NOTICE

- Do not use a squirt or spray bottle on or near the instrument as liquids will flow into the instrument and may cause permanent damage.
- Do not use hydrofluoric acid (HF) on the pedestals. Fluoride ions will permanently damage the quartz fiber optic cables.

#### Before you begin...


Before taking pedestal measurements with the NanoDrop Eight instrument, lift the instrument arm and clean the upper and lower pedestals. At a minimum, wipe the pedestals with a new laboratory wipe. For more information, see Cleaning the Pedestals.

#### To measure a Protein A280 sample

- 1. From the Home screen, select the Proteins tab and select Protein A280.
- 2. Specify a sample type and baseline correction if desired.
- 3. Pipette  $1-2 \mu$ L blanking solution onto the lower pedestal and lower the arm.
- 4. Click Blank and wait for the measurement to complete.

**Tip**: If Auto-Blank is On, the blank measurement starts automatically after you lower the arm.

- 5. Lift the arm and clean all pedestals with a new laboratory wipe.
- 6. Pipette 2 µL sample solution onto the pedestal and lower the arm.
- 7. Start the sample measurement:
  - If Auto-Measure is On, lower arm;
  - if Auto-Measure is off, lower arm and click

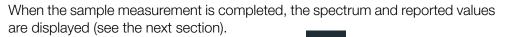


8. When you are finished measuring samples, click

End Experiment.


Measure.

9. Lift the arm and clean all pedestals with a new wipe.



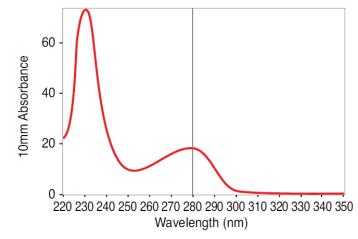

**High concentration BSA sample** 

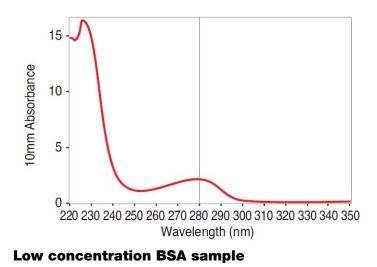

## Best practices for protein measurements

• Isolate and purify protein samples before measurement to remove impurities. Depending on the sample, impurities could include DNA, RNA and some buffer components. See Preparing Samples for more information.

**Note** Extraction reagents that contribute absorbance between 200 nm and 280 nm will affect measurement results if present in samples (even residual amounts).

• Ensure the sample absorbance is within the instrument's absorbance detection limits.

- Choosing a blank:
  - For the Protein A280 application, blank with the same buffer solution used to resuspend the analyte of interest. The blanking solution should be a similar pH and ionic strength as the analyte solution.
- Run a blanking cycle to assess the absorbance contribution of your buffer solution. If the buffer exhibits strong absorbance at or near the analysis wavelength (typically 280 nm or 205 nm), you may need to choose a different buffer or application. See Choosing and Measuring a Blank for more information.

**Note** Buffers such as Triton X, RIPA, and NDSB contribute significant absorbance and are not compatible with direct A280 or A205 measurements.

- For micro-volume measurements:
  - Ensure pedestal surfaces are properly cleaned and conditioned. (Proteins tend to stick to pedestal surfaces.)
  - Gently (but thoroughly) vortex samples before taking a measurement. Avoid introducing bubbles when mixing and pipetting.
  - Follow best practices for micro-volume measurements.
  - Use a 2 µL sample volume. See Recommended Sample Volumes for more information.

#### **Related Topics**


- Best practices for protein measurements
- Measure a Micro-Volume Sample
- Prepare Samples and Blanks
- Basic Instrument Operations



## **Protein A280 Reported Results**

#### Protein A280 measurement screen

For each measured sample, this application shows the absorbance spectrum and a measurement results. Here is an example:

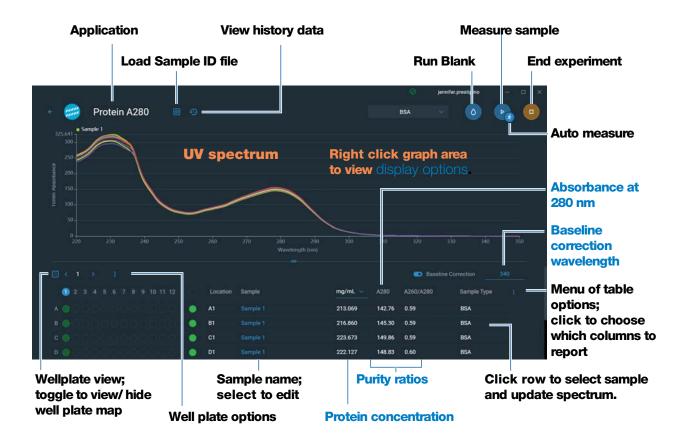

**Note** Micro-volume absorbance measurements are normalized to a 10.0 mm pathlength equivalent.

#### **Protein A280 reported values**

- Protein concentration
- Absorbance at 280 nm
- Purity ratio
- Sample type
- Baseline correction wavelength

#### **Thermo Scientific**

#### **5 Protein Applications** Measure Protein A280

- Baseline correction absorbance
- created on (date sample measurement was taken)
- monitor wavelength
- pathlength used
- Extinction Coefficient
- contaminant
- corrected

#### **Related Topics**


- Basic Instrument Operations
- Protein A280 Calculations

#### **Settings for Protein A280 Measurements**

To show the Protein A280 settings, from the Protein A280 measurement screen, select the settings icon **Protein A280 Setup**.

#### **Protein A280 settings**

The Protein A280 application provides a variety of sample type options for purified protein analysis.

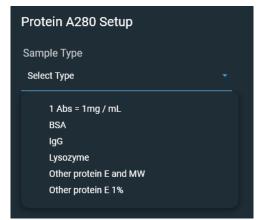

Each sample type applies a unique extinction coefficient to the protein calculations. If the extinction coefficient of the sample is known, choose the  $\mathcal{E}$  + MW (molar) or  $\mathcal{E}1\%$  (mass) option and enter the value. Otherwise, calculate the extinction coefficient or choose the option that best matches the sample solution. If you only need a rough estimate of protein concentration and the sample extinction coefficient is unknown, select the 1 Abs=1 mg/mL sample type option.

**Tip** Ideally, the extinction coefficient should be determined empirically using a solution of the study protein at a known concentration using the same buffer.

Setting	Available Options	Mass Ext. Coefficient (L/gm-cm)	Description
Sample type <sup>a</sup>	1 Abs = 1 mg/mL	General reference	Recommended when extinction coefficient is unknown and rough estimate of protein concentration is acceptable for a solution with no other interfering substances. Assumes 0.1% (1 mg/mL) protein solution produces 1.0A at 280 nm (where pathlength is 10 mm), i.e., $\varepsilon 1\% = 10$ .
	BSA	6.7	Calculates BSA (Bovine Serum Albumin) protein concentration using mass extinction coefficient (ε) of 6.7 L/gm-cm at 280 nm for 1% (i.e., 10 mg/mL) BSA solution. Assuming MW is 66,400 daltons (Da), molar extinction coefficient at 280 nm for BSA is approximately 43,824 M <sup>-1</sup> cm <sup>-1</sup> .
	IgG	13.7	Suitable for most mammalian antibodies (i.e., immunoglobulin G or IgG). Calculates protein concentration using mass extinction coefficient ( $\mathcal{E}$ ) of 13.7 L/gm-cm at 280 nm for 1% (i.e., 10 mg/mL) IgG solution. Assuming MW is 150,000 Da, molar extinction coefficient at 280 nm for IgG is approximately 210,000 M <sup>-1</sup> cm <sup>-1</sup> .
	Lysozyme	26.4	Calculates lysozyme protein concentration using mass extinction coefficient ( $\mathcal{E}$ ) of 26.4 L/gm-cm at 280 nm for 1% (i.e., 10 mg/mL) lysozyme solution. Assumes molar extinction coefficient for egg white lysozyme ranges between 36,000 M <sup>-1</sup> cm <sup>-1</sup> and 39,000 M <sup>-1</sup> cm <sup>-1</sup> .

Setting	Available Options	Mass Ext. Coefficient (L/gm-cm)	Description
	Other protein ( <b>E</b> + MW)	User entered molar extinction coefficient and molecular	Assumes protein has known molar extinction coefficient ( $\epsilon$ ) and molecular weight (MW), where:
		weight	$(\epsilon_{molar})^*10=(\epsilon_{percent})^*(MW_{protein})$
			Enter MW in kiloDaltons (kDa) and molar extinction coefficient ( $\epsilon$ ) in M <sup>-1</sup> cm <sup>-1</sup> divided by 1000 (i.e., $\epsilon$ /1000). For example, for protein with molar extinction coefficient of 210,000 M <sup>-1</sup> cm <sup>-1</sup> , enter 210.
	Other protein ( <b>E</b> 1%)	User entered mass extinction coefficient	Assumes protein has known mass extinction coefficient (£). Enter mass extinction coefficient in L/gm-cm for 10 mg/mL (£1%) protein solution.

<sup>a</sup> To add or edit a custom protein, use Protein Editor.

Setting	Available Options	Mass Ext. Coefficient (L/gm-cm)	Description
Baseline Correction	On or off Enter baseline correction wavelength in nm or use default value (340 nm)	N/A	Corrects for any offset caused by light scattering particulates by subtracting measured absorbance at specified baseline correction wavelength from absorbance values at all wavelengths in sample spectrum. As a result, absorbance of sample spectrum is zero at specified baseline correction wavelength. <b>Tip</b> : If the sample has a modification that absorbs light at 340 nm, select a different correction wavelength or turn off Baseline Correction.



## **Protein editor**

Use the Protein Editor to add a custom protein to the list of available protein sample types in Protein A280 Setup.

#### To access the Protein Editor:

• From the Home screen, select 😥 Settings > Protein Editor.

Protein E	ditor	Cli	ck to add c	ustom protein				
<ul> <li>← Settings</li> </ul>	System	Protein Editor	Preferences		Ø	operator	-3	
<u>_</u> Q						( + 1	Protein	$\supset$
Туре	Name	ɛ/1000	MW(kDA)	Ext.Coeff.E1%L/gm-cm	Description			
Mass	Protein AA			44.77		1	ī s'	Ð
list in Prote	oteins <sup>-</sup> in Sample T in A280 Setu is & Labels S	ype p Cli		elected custom	-			

These operations are available from the Protein Editor:

#### Add custom protein

- 1. In Protein Editor, select + Protein to show the **New Protein Type** box.
- 2. Enter a unique Name for the new protein.
- 3. Enter a **Description** for the new protein.
- 4. Specify whether to enter **Molar Extinction** coefficient or **Mass Extinction** coefficient for custom protein.
  - If Mass Extinction coefficient is selected, enter mass extinction coefficient in L/gm-cm for 10 mg/mL (£1%) protein solution.

New Protein Typ	e		×
• Molar Extinction	O Mass Extinction		
Name		Description	
ε/1000	MW(kDA)		
0.01			

#### - if **Molar Extinction** is selected,

- enter molar extinction coefficient (E) in M<sup>-1</sup>cm<sup>-1</sup> divided by 1000 (that is, E/1000). For example, for protein with molar extinction coefficient of 210,000 M<sup>-1</sup>cm<sup>-1</sup>, enter 210.
- Enter molecular weight (MW) in kiloDaltons (kDa)
- 5. Select Save to close the New Protein Type box.

The new custom protein appears in the **Type** list in Protein A280 Setup and Proteins & Labels Setup.

#### **Edit custom protein**

- 1. In Protein Editor, click to select custom protein
- 2. Click **m** to show the **Edit Protein Type** box
- 3. Edit any entries or settings
- 4. Click OK

#### **Delete custom protein**

- 1. In Protein Editor, click to select a custom protein to delete
- 2. click 📋

**Note** Deleting a custom protein permanently removes the protein and all associated information from the software.



## **Detection Limits for Protein A280 Measurements**

Detection limits and reproducibility specifications for purified BSA proteins are provided here. The BSA lower detection limit and reproducibility values apply to any protein sample type. The upper detection limits are dependent on the upper absorbance limit of the instrument and the sample's extinction coefficient.

#### To calculate upper detection limits for other (non-BSA) protein sample types

To calculate upper detection limits in mg/mL for proteins, use the following equation:

(upper absorbance limit<sub>instrument</sub> /mass extinction coefficient<sub>sample</sub>) \* 10

For example, if the sample's mass extinction coefficient at 280 nm is 6.7 for a 1% (10 mg/mL) solution, the equation looks like this:

(200 / 6.7) \* 10 = 298.5 (or ~300)

## **Calculations for Protein A280 Measurements**

The Protein A280 application uses the Beer-Lambert equation to correlate absorbance with concentration. Solving Beer's law for concentration yields the equation at the right.

#### **Beer-Lambert Equation (solved for concentration)**

$$c = A / (\varepsilon * b)$$

where:

A = UV absorbance in absorbance units (AU)

 $\epsilon$  = wavelength-dependent molar absorptivity coefficient (or extinction coefficient) in liter/mol-cm

b = pathlength in cm

c = analyte concentration in moles/liter or molarity (M)

**Note**: Dividing the measured absorbance of a sample solution by its molar extinction coefficient yields the molar concentration of the sample. See Published Extinction Coefficients for more information regarding molar vs. mass concentration values.

#### **5** Protein Applications Measure Protein A280

The extinction coefficient of a peptide or protein is related to its tryptophan (W), tyrosin (Y) and cysteine (C) amino acid composition.

**Tip**: The extinction coefficient is can be affected by buffer type, ionic strength and pH.

wavelength specific for each protein and

where:

equation:

 $\boldsymbol{\varepsilon}$  = molar extinction coefficient **n** = number of each amino acid residue

**Extinction Coefficients for Proteins** 

5500, 1490 and 125 = amino acid molar absorptivities at 280 nm

At 280 nm, the extinction coefficient is approximated by the

weighted sum of the 280 nm molar extinction coefficients of the three constituent amino acids, as described in this

 $\mathcal{E} = (nW * 5500) + (nY * 1490) + (nC * 125)$ 

#### **Available Options for Extinction Coefficient**

- 1 Abs = 1 mg/mL, where sample type and/or ext. coefficient is unknown (produces rough estimate of protein concentration)
- **BSA** (Bovine Serum Albumin, 6.7 L/gm-cm)
- **IgG** (any mammalian antibody, 13.7 L/gm-cm)
- Lysozyme (egg white lysozyme, 26.4 L/gm-cm)
- Other protein (E + MW), user-specified molar ext. coefficient
- Other protein (£1%), user-specified mass ext. coefficient

**Note**: See Sample Type for details.

This application offers six options (shown at right) for selecting an appropriate extinction coefficient for each measured sample, to be used in conjunction with Beer's Law to calculate sample concentration.

If the extinction coefficient of the sample is known, choose the  $\epsilon$  + MW (molar) or ε1% (mass) option and enter the value. Otherwise, calculate the extinction coefficient or choose the option that best matches the sample solution.

**Tip**: Ideally, the extinction coefficient should be determined empirically using a solution of the study protein at a known concentration using the same buffer.



Most sources report extinction coefficients for proteins measured at or near 280 nm in phosphate or other physiologic buffer. These values provide sufficient accuracy for routine assessments of protein concentration.

The equation at the right shows the relationship between molar extinction coefficient ( $\epsilon_{molar}$ ) and percent extinction coefficient ( $\epsilon_{0}$ ).

To determine concentration (c) of a sample in mg/mL, use the equation at the right and a conversion factor of 10.

**Tip**: The NanoDrop Eight software includes the conversion factor when reporting protein concentrations.

#### **Published Extinction Coefficients**

Published extinction coefficients for proteins may be reported as:

- wavelength-dependent molar absorptivity (or extinction) coefficient (ε) with units of M<sup>-1</sup>cm<sup>-1</sup>
- percent solution extinction coefficient (£1%) with units of (g/100 mL)<sup>-1</sup>cm<sup>-1</sup> (i.e., 1% or 1 g/100 mL solution measured in a 1 cm cuvette)
- protein absorbance values for 0.1% (i.e., 1 mg/mL) solutions

**Tip**: Assess published values carefully to ensure unit of measure is applied correctly.

#### Conversions Between $\epsilon_{\text{molar}}$ and $\epsilon_{\text{1\%}}$

$$(\varepsilon_{molar}) * 10 = (\varepsilon 1\%) * (MW_{protein})$$

Example: To determine percent solution extinction coefficient ( $\pounds$ 1%) for a protein that has a molar extinction coefficient of 43,824 M<sup>-1</sup>cm<sup>-1</sup> and a molecular weight (MW) of 66,400 daltons (Da), rearrange and solve the above equation as follows:

 $\varepsilon 1\% = (\varepsilon_{molar} * 10) / (MW_{protein})$ 

£1% = (43,824 \* 10) / 66,400 Da)

## $\epsilon$ 1% = 6.6 g/100 mL

Conversions Between g/100 mL and mg/mL

 $C_{\text{protein}}$  in mg/mL = (A /  $\varepsilon$ 1%) \* 10

Example: If measured absorbance for a protein sample at 280 nm relative to the reference is 5.8 A, protein concentration can be calculated as:

 $C_{protein} = (A / \epsilon 1\%) * 10$ 

$$C_{\text{protein}} = (5.8/6.6 \text{ g}/100 \text{ mL}) * 10$$

 $C_{\text{protein}} = 8.79 \text{ mg/mL}$ 

Calculated protein concentrations are based on the absorbance value at 280 nm, the selected (or entered) extinction coefficient and the sample pathlength. A single-point baseline correction (or analysis correction) may be applied.

Concentration is reported in mass units. Calculators are available on the Internet to convert concentration from mass to molar units based on sample sequence.

Absorbance values at 260 nm and 280 nm are used to calculate purity ratios for the measured protein samples.

Purity ratios are sensitive to the presence of contaminants in the sample, such as residual solvents and reagents typically used during sample purification.

#### **Measured Values**

#### A280 absorbance

**Note**: For micro-volume absorbance measurements, the spectra are normalized to a 10 mm pathlength equivalent.

- Protein absorbance values are measured at 280 nm using the normalized spectrum. If Baseline Correction is not selected, this is the reported A280 value and the value used to calculate protein concentration.
- If Baseline Correction is selected, the normalized and baseline-corrected absorbance value at 280 nm is reported and used to calculate protein concentration.

#### Sample Pathlength

- For micro-volume measurements, the software selects the optimal pathlength (between 1.0 mm and 0.1 mm) based on sample absorbance at the analysis wavelength.
- Displayed spectra and absorbance values are normalized to a 10 mm pathlength equivalent.

#### **Reported Values**

- Protein concentration. Reported in selected unit (mg/mL or µg/mL). Calculations are based on Beer-Lambert equation using corrected protein absorbance value.
- **A260/A280 purity ratio**. Ratio of corrected absorbance at 260 nm to corrected absorbance at 280 nm. An A260/A280 purity ratio of ~0.57 is generally accepted as "pure" for proteins.

**Note**: Although purity ratios are important indicators of sample quality, the best indicator of protein quality is functionality in the downstream application of interest (e.g., real-time PCR).

- **Contaminant** Displays contaminant identified by the Acclaro software, if one is available.
- **Corrected** Display the corrected analyte concentration determined using the Acclaro software, if one is available.

Protein Applications Measure Protein A205

## **Measure Protein A205**

Measures the concentration of purified protein populations that absorb at 205 nm.


Measure A205 Proteins

**Reported Results** 

Settings

**Detection Limits** 

Calculations

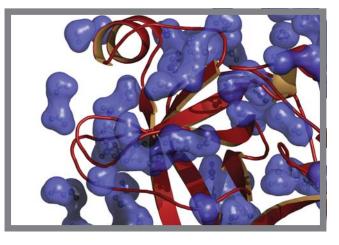

## **Measure Protein Concentration at A205**

Use the Protein A205 application to quantify purified peptides and other proteins that contain peptide bonds, which exhibit absorbance at 205 nm. This application reports protein concentration and two absorbance values (A205 and A280). A single-point baseline correction can also be used. This application does not require a standard curve.

**Note** If your samples contain mainly amino acids such as tryptophan or tyrosine, or cys-cys disulfide bonds, use the Protein A280 application instead of Protein A205.

#### To measure Protein A205 samples

#### NOTICE

- Do not use a squirt or spray bottle on or near the instrument as liquids will flow into the instrument and may cause permanent damage.
- Do not use hydrofluoric acid (HF) on the pedestals. Fluoride ions will permanently damage the quartz fiber optic cables.

#### Before you begin...

Before taking pedestal measurements with the NanoDrop Eight instrument, lift the instrument arm and clean the upper and lower pedestals. At a minimum, wipe the pedestals with a new laboratory wipe. For more information, see Cleaning the Pedestals.

#### To measure a Protein A205 sample

- 1. From the Home screen, select the Proteins tab and select Protein A205.
- 2. Specify a sample type and baseline correction if desired.
- 3. Pipette  $1-2 \mu$ L blanking solution onto the lower pedestal and lower the arm.
- 4. Click G Blank and wait for the measurement to complete.

**Tip**: If Auto-Blank is On, the blank measurement starts automatically after you lower the arm.

- 5. Lift the arm and clean all pedestals with a new laboratory wipe.
- 6. Pipette 2 µL sample solution onto the pedestal and lower the arm.
- 7. Start the sample measurement:
  - If Auto-Measure is On, lower arm;
  - if Auto-Measure is off, lower arm and click

When the sample measurement is completed, the spectrum and reported values are displayed (see the next section).

8. When you are finished measuring samples, click

End Experiment.

Measure.

9. Lift the arm and clean all pedestals with a new wipe.

#### **Related Topics**


- Best Practices for Protein Measurements
- Measure a Micro-Volume Sample
- Prepare Samples and Blanks
- Basic Instrument Operations



## **Protein A205 Reported Results**

#### Protein A205 measurement screen

For each measured sample, this application shows the absorbance spectrum and a summary of the results. Here is an example:

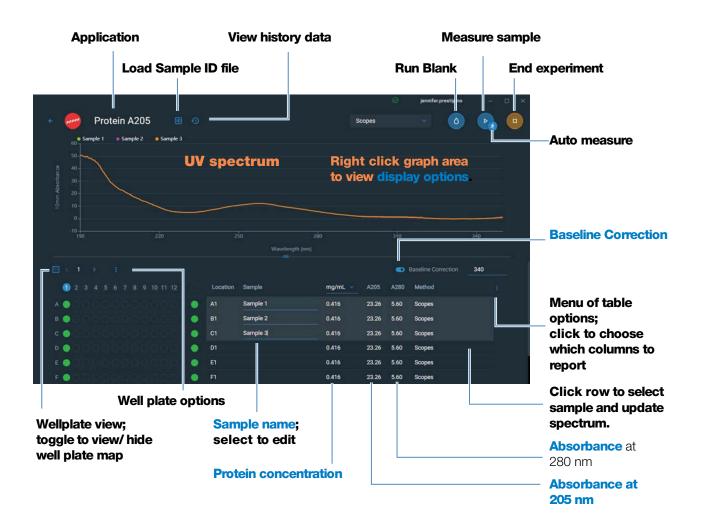

#### **Protein A205 reported values**

- Protein concentration
- Absorbance at 205 nm
- Absorbance at 280 nm
- Sample type
- Baseline correction

- created on (date sample measurement was taken)
- monitor wavelength
- pathlength used
- Extinction Coefficient

#### **Related Topics**

- Basic Instrument Operations
- Protein A205 Calculations

## **Settings for Protein A205 Measurements**

To show the Protein A205 settings, from the Protein A205 measurement screen, select Protein A205 Setup

Setting	Available Options	Mass Ext. Coefficient (L/gm-cm)	Description
Sample type	31	31	Assumes $\boldsymbol{\epsilon}$ 0.1% (1 mg/mL) at 205 nm = 31
	Scopes	27 + 120 * (A280/A205)	Assumes <b>&amp;</b> 0.1% (1 mg/mL) at 205 nm = 27 + 120 * (A280/A205)
	Other protein ( <b>ɛ</b> 1%)	User entered mass extinction coefficient	Assumes protein has known mass extinction coefficient (£). Enter mass extinction coefficient in L/gm-cm for 1 mg/mL (£0.1%) protein solution.
Baseline Correction	On or off Enter baseline correction wavelength in nm or use default value (340 nm)	N/A	Corrects for any offset caused by light scattering particulates by subtracting measured absorbance at specified baseline correction wavelength from absorbance values at all wavelengths in sample spectrum. As a result, absorbance of sample spectrum is zero at specified baseline correction wavelength.
			<b>Tip</b> : If the sample has a modification that absorbs light at 340 nm, select a different correction wavelength or turn off Baseline Correction.

#### **Related Topics**

• Instrument Settings



## **Calculations for Protein A205 Measurements**

As with the other protein applications, Proteins A205 uses the Beer-Lambert equation to correlate absorbance with concentration based on the sample's extinction coefficient and pathlength.

This application offers three options (shown at right) for selecting an appropriate extinction coefficient for each measured sample, to be used in conjunction with Beer's Law to calculate sample concentration.

If the extinction coefficient of the sample is known, choose the  $\mathcal{E}1\%$  (mass) option and enter the value. Otherwise, calculate the extinction coefficient or choose the option that best matches the sample solution.

**Tip**: Ideally, the extinction coefficient should be determined empirically using a solution of the study protein at a known concentration using the same buffer.

Calculated protein concentrations are based on the absorbance value at 205 nm, the selected (or entered) extinction coefficient and the sample pathlength. A single-point baseline correction may also be applied.

Concentration is reported in mass units. Calculators are available on the Internet to convert concentration from mass to molar units based on the sample sequence.

#### **Available Options for Extinction Coefficient**

- **31**, assumes £0.1% (1 mg/mL) at 205 nm = 31
- **Scopes**, assumes £0.1% (1 mg/mL) at 205 nm = 27 + 120 \* (A280/A205)
- **Other protein**, enter mass extinction coefficient in L/gm-cm for 1 mg/mL (£0.1%) protein solution

**Note**: See Sample Type for details.

#### **Measured Values**

#### A205 absorbance

**Note**: For micro-volume absorbance measurements, the spectra are normalized to a 10 mm pathlength equivalent.

- Protein absorbance values are measured at 205 nm using the normalized spectrum. If Baseline Correction is not selected, this is the reported A205 value and the value used to calculate protein concentration.
- If Baseline Correction is selected, the normalized and baseline-corrected absorbance value at 205 nm is reported and used to calculate protein concentration.

#### A280 absorbance

• Normalized and baseline-corrected (if selected) absorbance value at 280 nm is also reported.

#### **Sample Pathlength**

- For micro-volume measurements, the software selects the optimal pathlength (between 1.0 mm and 0.1 mm) based on sample absorbance at the analysis wavelength.
- Displayed spectra and absorbance values are normalized to a 10 mm pathlength equivalent.

#### **Reported Values**

• **Protein concentration.** Reported in selected unit (mg/mL or µg/mL). Calculations are based on Beer-Lambert equation using corrected protein absorbance value.

#### **Related Topics**

- Beer-Lambert Equation
- Protein A280 Calculations

## **Measure Proteins and Labels**

Measures the concentration of purified proteins that have been labeled with up to two fluorescent dyes.

Measure Labeled Proteins

**Reported Results** 

Settings

**Detection Limits** 

Calculations



## **Measure Labeled Protein Samples**

Use the Proteins and Labels application to quantify proteins and fluorescent dyes for protein array conjugates, as well as metalloproteins such as hemoglobin, using wavelength ratios. This application reports protein concentration measured at 280 nm, an A269/A280 absorbance ratio, and the concentrations and measured absorbance values of the dyes, allowing detection of dye concentrations as low as 0.2 picomole per microliter. This information is useful for evaluating protein/dye conjugation (degree of labeling) for use in downstream applications.

#### To measure labeled protein samples

#### NOTICE

- Do not use a squirt or spray bottle on or near the instrument as liquids will flow into the instrument and may cause permanent damage.
- Do not use hydrofluoric acid (HF) on the pedestals. Fluoride ions will permanently damage the quartz fiber optic cables.

#### Before you begin...

Before taking pedestal measurements with the NanoDrop Eight instrument, lift the instrument arm and clean the upper and lower pedestals. At a minimum, wipe the pedestals with a new laboratory wipe. For more information, see Cleaning the Pedestals.

#### To measure a labeled protein sample

1. From the Home screen, select the **Proteins** tab and then select **Protein & Labels**.

2. Specify the sample type and the type of dye(s) used.

**Tip**: Select a dye from the pre-defined list or add a custom dye using the Dye/Chromophore Editor.


- Pipette 1–2 μL of the blanking solution onto the lower pedestal and lower the arm.
- 4. Tap **Blank** and wait for the measurement to complete.

**Tip**: If Auto-Blank is On, the blank measurement starts automatically after you lower the arm.

- 5. Lift the arm and clean all pedestals with a new laboratory wipe.
- 6. Pipette 2 µL sample solution onto the pedestal and lower the arm.
- 7. Start the sample measurement:
  - If Auto-Measure is On, lower arm;
  - if Auto-Measure is off, lower arm and select **Measure**.

When the sample measurement is completed, the spectrum and reported values are displayed (see the next section).

- 8. When you are finished measuring samples, select **End Experiment**.
- 9. Lift the arm and clean all pedestals with a new wipe.

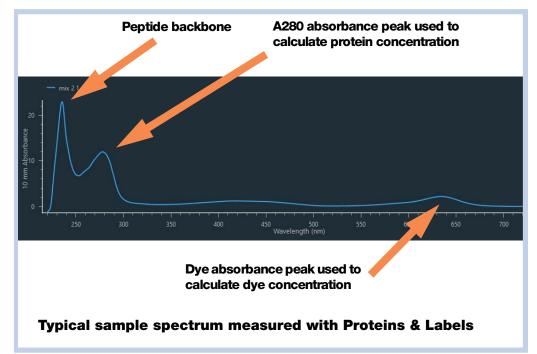


#### **Related Topics**

- Best practices for protein measurements
- Measure a Micro-Volume Sample
- Prepare Samples and Blanks
- Basic Instrument Operations

## **Proteins & Labels Reported Results**

#### **Proteins & Labels measurement screen**

For each measured sample, this application shows the absorbance spectrum and a summary of the results. Below is an example of the measurement screen:

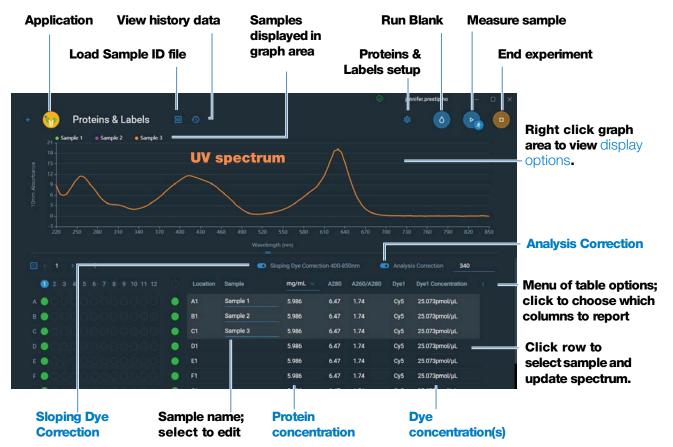

#### Note

- A baseline correction is performed at 340 nm (absorbance value at 340 nm is subtracted from absorbance values at all wavelengths in sample spectrum).
- Micro-volume absorbance measurements are normalized to a 10.0 mm pathlength equivalent.

#### **Proteins & Labels reported values**

The initial screen that appears after each measurement (see previous image) shows a summary of the reported values. To view all reported values, press and hold the sample row. Here is an example:

#### **Reported values for Proteins & Labels application**

- Sample Name
- Creation date
- Protein
- A280
- Sample Type
- Dye 1/Dye 2
- Sloping Dye Correction
- Analysis Correction

#### **Related Topics**

- Basic Instrument Operations
- Proteins & Labels calculations

## **Settings for Proteins and Labels Measurements**

To show the Proteins & Labels settings, from the Proteins & Labels measurement screen, select roteins & Labels Setup.

#### **5** Protein Applications Measure Proteins and Labels

Setting	Available Options	Mass Ext. Coefficient (L/gm-cm)	Description
Sample	1 Abs = 1 mg/mL	General reference	Select Sample type for detailed description of
type <sup>a</sup>	BSA	6.7	each available setting.
	lgG	13.7	Each sample type applies a unique extinction coefficient to the protein calculations. If the
	Lysozyme	26.4	extinction coefficient of the sample is known, choose the $\mathcal{E}$ + MW (molar) or $\mathcal{E}1\%$ (mass)
	Other protein ( $\varepsilon$ + MW) Other protein ( $\varepsilon$ 1%)	user-entered molar extinction coefficient/ molecular weight User entered mass extinction coefficient	option and enter the value. Otherwise, calculate the extinction coefficient or choose the option that best matches the sample solution. If you only need a rough estimate of protein concentration and the sample extinction coefficient is unknown, select the 1 Abs=1 mg/mL sample type option.
		Coemcient	<b>Tip</b> : Ideally, the extinction coefficient should be determined empirically using a solution of the study protein at a known concentration using the same buffer.
Analysis Correction <sup>b</sup>	On or off Enter analysis correction wavelength in nm or use default value (340 nm)	N/A	Corrects sample absorbance measurement for any offset caused by light scattering particulates by subtracting absorbance value at specified analysis correction wavelength from absorbance value at analysis wavelength. Corrected value is used to calculate sample concentration.
			<b>Tip</b> : If the sample has a modification that absorbs light at 340 nm, select a different correction wavelength or turn off Analysis Correction.
Dye 1/Dye 2 Type <sup>c</sup>	Cy3, 5, 3.5, or 5.5, Alexa Fluor 488, 546, 555, 594, 647, or 660	See Dye/Chromophore Editor for specific values for each dye	Select pre-defined dye used to label sample material, or one that has been added using Dye/Chrom. Editor.
Dye 1/Dye 2 Unit	picomoles/microliter (pmol/µl), micromoles (uM), or millimoles (mM)	not applicable	Select unit for reporting dye concentrations.

#### **5 Protein Applications** Measure Proteins and Labels

Setting	Available Options	Mass Ext. Coefficient (L/gm-cm)	Description
Sloping Dye Correction <sup>d</sup>	On or off		Corrects dye absorbance measurements for any offset caused by light scattering particulates by subtracting absorbance value of a sloping baseline from 400 nm to 850 nm from absorbance value at dye's analysis wavelength.

<sup>a</sup> To add or edit a custom protein, use Protein Editor.

<sup>b</sup> Analysis Correction affects calculation for protein concentration only.

<sup>c</sup> To add custom dye or edit list of available dyes, use Dye/Chromophore Editor.

<sup>d</sup> Sloping Dye Correction affects calculations for dye concentration only.

#### **Related Topics**

- Instrument Settings
- Protein Editor
- Dye/Chromophore Editor

### **Detection Limits for Proteins and Labels Measurements**

Detection limits and reproducibility specifications for purified BSA proteins and dyes that are pre-defined in the software are provided here. The BSA lower detection limit and reproducibility values apply to any protein sample type. The upper detection limits are dependent on the upper absorbance limit of the instrument and the sample's extinction coefficient.

#### To calculate upper detection limits for other (non-BSA) protein sample types

To calculate upper detection limits in mg/mL for proteins, use the following equation:

(upper absorbance limit<sub>instrument</sub> /mass extinction coefficient<sub>sample</sub>) \* 10

For example, if the sample's mass extinction coefficient at 280 nm is 6.7 for a 1% (10 mg/mL) solution, the equation looks like this:

(200 / 6.7) \* 10 = 298.5 (or ~300)

#### **Related Topics**

• Detection Limits for All Applications

### **Calculations for Proteins and Labels Measurements**

As with the other protein applications, Proteins & Labels uses the Beer-Lambert equation to correlate absorbance with concentration based on the sample's extinction coefficient and pathlength.

This application offers six options (shown at right) for selecting an appropriate extinction coefficient for each measured sample, to be used in conjunction with Beer's Law to calculate sample concentration.

If the extinction coefficient of the sample is known, choose the  $\epsilon$  + MW (molar) or  $\epsilon$ 1% (mass) option and enter the value. Otherwise, calculate the extinction coefficient or choose the option that best matches the sample solution.

**Tip**: Ideally, the extinction coefficient should be determined empirically using a solution of the study protein at a known concentration using the same buffer.

Calculated protein concentrations are based on the absorbance value at 280 nm, the selected (or entered) extinction coefficient and the sample pathlength. A single-point baseline correction (or analysis correction) may be applied.

Concentration is reported in mass units. Calculators are available on the Internet to convert concentration from mass to molar units based on sample sequence.

#### Available Options for Extinction Coefficient

- **1** Abs = **1** mg/mL, where sample type and/or ext. coefficient is unknown (produces rough estimate of protein concentration)
- BSA (Bovine Serum Albumin, 6.7 L/gm-cm)
- IgG (any mammalian antibody, 13.7 L/gm-cm)
- Lysozyme (egg white lysozyme, 26.4 L/gm-cm)
- **Other protein** (ε + MW), user-specified molar ext. coefficient
- **Other protein** (£1%), user-specified mass ext. coefficient

Note: See Sample Type for details.

#### **Measured Values**

#### A280 absorbance

**Note**: The absorbance value at 850 nm is subtracted from all wavelengths in the spectrum. As a result, the absorbance at 850 nm is zero in the displayed spectra. Also, for micro-volume absorbance measurements, the spectra are normalized to a 10 mm pathlength equivalent.

- Protein absorbance values are measured at 280 nm using the Analysis-corrected and normalized spectrum. If Analysis Correction and Dye Correction are not selected, this is the reported A280 value and the value used to calculate protein concentration.
- If Analysis Correction is selected, the normalized and analysis-corrected absorbance value at 280 nm is reported and used to calculate protein concentration.
- If a Dye is used, the normalized, analysis-corrected and dye-corrected absorbance value at 280 nm is reported and used to calculate protein concentration.

Dye concentrations are calculated from the absorbance value at the dye's analysis wavelength, the dye's extinction coefficient, and the sample pathlength. A sloped-line dye correction may also be used.

#### Dye absorbance

- Dye absorbance values are measured at specific wavelengths. See Dye/Chromophore Editor for analysis wavelengths used.
- If Sloping Dye Correction is selected, a linear baseline is drawn between 400 nm and 850 nm and, for each dye, the absorbance value of the sloping baseline is subtracted from the absorbance value at each dye's analysis wavelength. Baseline-corrected dye absorbance values are reported and used to calculate dye concentrations.

#### **Dye correction**

- Pre-defined dyes have known correction values for A260 and A280. See Dye/Chromophore Editor for correction values used.
- A280 dye correction is subtracted from A280 absorbance value used to calculate protein concentration.

#### **Sample Pathlength**

- For micro-volume measurements, the software selects the optimal pathlength (between 1.0 mm and 0.1 mm) based on sample absorbance at the analysis wavelength.
- Displayed spectra and absorbance values are normalized to a 10 mm pathlength equivalent.

#### **Reported Values**

- **Protein concentration.** Reported in selected unit (mg/mL or µg/mL). Calculations are based on Beer-Lambert equation using corrected protein absorbance value.
- **Dye1/Dye2 concentration**. Reported in pmol/µL. Calculations are based on Beer's Law equation using (sloping) baseline-corrected dye absorbance value(s).

#### **Related Topics**

- Beer-Lambert Equation
- Protein A280 Calculations

6

# **Custom Applications**

Use the NanoDrop Eight to perform UV-Vis measurements.

The UV-Vis application allows the instrument to function as a conventional spectrophotometer. Up to 40 wavelengths from 190 nm to 850 nm can be monitored and reported.

- Measure UV-Vis ..... page 82
- Measure Custom ..... page 87

6 Custom Applications Measure UV-Vis

### **Measure UV-Vis**

Measures the absorbance of any sample at up to 40 wavelengths across the ultra-violet (UV) and visible regions of the spectrum.

Measure UV-Vis

**Reported Results** 

Settings

**Detection Limits** 



### **Measure UV-Vis**

The UV-Vis application allows the instrument to function as a conventional spectrophotometer. Sample absorbance is displayed on the screen from 190 nm to 850 nm. Up to 40 wavelengths can be designated for absorbance monitoring and inclusion in the report. Automatic pathlength adjustment and a single-point baseline correction can also be used.

#### To make UV-Vis measurements

#### NOTICE

- Do not use a squirt or spray bottle on or near the instrument as liquids will flow into the instrument and may cause permanent damage.
- Do not use hydrofluoric acid (HF) on the pedestals. Fluoride ions will permanently damage the quartz fiber optic cables.

#### Before you begin...

Before taking pedestal measurements with the NanoDrop Eight instrument, lift the instrument arm and clean the upper and lower pedestals. At a minimum, wipe the pedestals with a new laboratory wipe. For more information, see Cleaning the Pedestals.

#### To measure a sample using the UV-Vis application

- 1. From the Home screen, from the Custom tab, select UV-Vis.
- 2. Specify up to 40 wavelengths to monitor (or you can specify them later if desired) and whether automated pathlength adjustment, analysis wavelength, and baseline correction will be used.
- 3. Pipette 1–2 µL blanking solution onto the lower pedestal and lower the arm.

#### 6 Custom Applications Measure UV-Vis

4. Select **Blank** and wait for the measurement to complete.

**Tip**: If Auto-Blank is On, the blank measurement starts automatically after you lower the arm.

- 5. Lift the arm and clean all pedestals with a new laboratory wipe, .
- 6. Pipette 1-2 µL sample solution onto the pedestal and lower the arm.
- 7. Start the sample measurement: If Auto-Measure is On, lower arm; if Auto-Measure is off, lower arm and select **Measure**.

When the sample measurement is completed, the spectrum and reported values are displayed (see the next section).

- 8. When you are finished measuring samples, select End Experiment
- 9. Lift the arm and clean all pedestals with a new wipe.

### **Best practices for UV-Vis measurements**

- Ensure the sample absorbance is within the instrument's absorbance detection limits.
- Blank with the same buffer solution used to re-suspend the analyte of interest. The blanking solution should be a similar pH and ionic strength as the analyte solution.
- Run a blanking cycle to assess the absorbance contribution of your buffer solution. If the buffer exhibits strong absorbance at or near an analysis wavelength, you may need to choose a different buffer or application. See Choosing and Measuring a Blank for more information.
- For micro-volume measurements:
  - Ensure pedestal surfaces are properly cleaned and conditioned.
  - Ensure samples are homogeneous before taking a measurement. Avoid introducing bubbles when mixing and pipetting.
  - Follow best practices for micro-volume measurements.
  - Use a 1-2 µL sample volume. See Recommended Sample Volumes for more information.

### **UV-Vis Reported Results**

#### **UV-Vis measurement screen**


For each measured sample, the absorbance spectrum and results are shown:



Shift-click multiple sample rows to overlay spectra Click a sample and hover locations on spectra to view measurement values

**Note** Micro-volume absorbance measurements are normalized to a 10.0 mm pathlength equivalent.

#### **UV-Vis reported values**

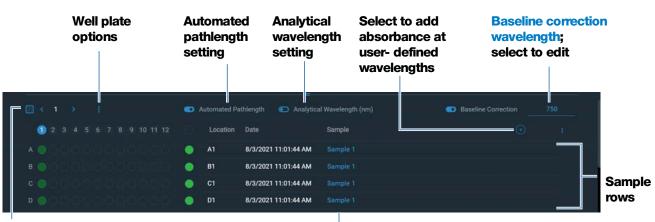

The lower half of the measurement screen shows the reported values:

Wellplate view; toggle to view/ hide well plate map

# Sample names; select to edit

#### Tips:

Click sample row to select sample and update spectrum Shift-click multiple sample rows to overlay spectra Click a sample and hover locations on spectra to view measurement values

### **Settings for UV-Vis Measurements**

To show the UV-Vis settings, from the Home screen, from the **Custom** tab, select **UV-Vis**.

Setting	Available Options	Description
Monitored wavelengths	Enter up to 40 wavelengths between 190 nm and 850 nm	<b>User-defined wavelengths to be measured and</b> <b>reported at run time</b> . Absorbance values for the first three entered wavelengths are displayed in the measurement screen. To see absorbance values for 8 monitored wavelengths, swipe left in the measurement screen to show the Data table. To see all monitored wavelengths, press and hold a sample row to show the Sample Details screen (scroll up to display absorbance values for any additional user-defined wavelengths).
		<b>Note</b> : If Baseline Correction is selected, all displayed absorbance values are the corrected values.
Analytical Wavelength	Any wavelength between 190 nm and 850 nm	This is the wavelength the software will use to determine the pathlength selection.
Automated Pathlength	On or Off (affects pedestal measurements only)	<b>Optional automated pathlength selection</b> . Allows the software to use the optimal (shorter) pedestal pathlength for high concentration samples to help prevent detector saturation (see Detection Limits for details).
		<ul> <li>When selected, the shorter pathlength is used when any wavelength has 10 mm equivalent absorbance value of 12.5 or higher.</li> </ul>
		• When deselected, the pedestal pathlength is restricted to 1.0 mm across all wavelengths.
		<b>Note</b> : In either case, displayed absorbance values have been normalized to a 10 mm pathlength equivalent.
Baseline Correction	On or off Enter baseline correction wavelength in nm or use default value (750 nm)	<b>Optional user-defined baseline correction</b> . Can be used to correct for any offset caused by light scattering particulates by subtracting measured absorbance at specified baseline correction wavelength from absorbance values at all wavelengths in sample spectrum. As a result, absorbance of sample spectrum is zero at specified baseline correction wavelength.

### **Measure Custom**

Runs a custom measurement method created using NanoDrop Eight software.

Measure Custom Method

Delete Custom Method

Reported Results



### Measure using a Custom Method

Use the Custom application to run a user-defined method created using the NanoDrop Eight software. For more information, see "Create Custom Method" on page 90.

Custom methods can only be created on a personal computer running the NanoDrop Eight software.

#### To measure using a custom method

#### NOTICE

- Do not use a squirt or spray bottle on or near the instrument as liquids will flow into the instrument and may cause permanent damage.
- Do not use hydrofluoric acid (HF) on the pedestals. Fluoride ions will permanently damage the quartz fiber optic cables.

#### Before you begin...

Before taking measurements with the NanoDrop Eight instrument, lift the instrument arm and clean the upper and lower pedestals. At a minimum, wipe the pedestals with a new laboratory wipe. For more information, see Cleaning the Pedestals.

#### To measure a sample using a custom method

1. From the Home screen, from the **Custom** tab, select **Custom Methods**.

NANODROP EIGHT	thermoscie	ntific		8
Ouston	Nucleic Acid	Protein OD60	0 Custom	
Custom Tab				
Custom ——— Methods Icon	f(x)			
	Custom Method	s	UV-Vis	

2. In the method selection pane, select the method to run.

Method custom 01 12/2/2021 12:51:33 PM operator  Inputs Formula Standards  Result Name Concentration ng/µL Measurement Range UV range 190-350 (nm) Analysis wavelength 340 nm Correction for Analysis Wavelength None Miscellanous Baseline Correction (nm)	۹	Method Name 🔱	Method custom 01		► Run
Measurement Range UV range 190-350 (nm) Analysis wavelength 340 nm Correction for Analysis Wavelength None					
Analysis wavelength 340 nm Correction for Analysis Wavelength None			Result Name	Concentration	ng/µL
Correction for Analysis Wavelength None			Measurement Range	UV range	190-350 (nm)
			Analysis wavelength	340	nm
Miscellanous Baseline Correction (nm)			Correction for Analysis Wavelength	None	
			Miscellanous	Baseline Correction (nm)	
Automated Pathlength				Automated Pathlength	

Information about the selected method appears in the method details pane.

3. Select ▷ Run

4. Follow the on-screen instructions to measure a sample.

### **Delete Custom Method**

- From Home screen, from the Custom tab, select Custom Methods.
- In Select Method box, select a method to delete


From the drop-down menu



### **Custom Method Reported Results**

#### **Custom method measurement screen**

For each measured sample, this application shows the absorbance spectrum and results details. Here is an example:

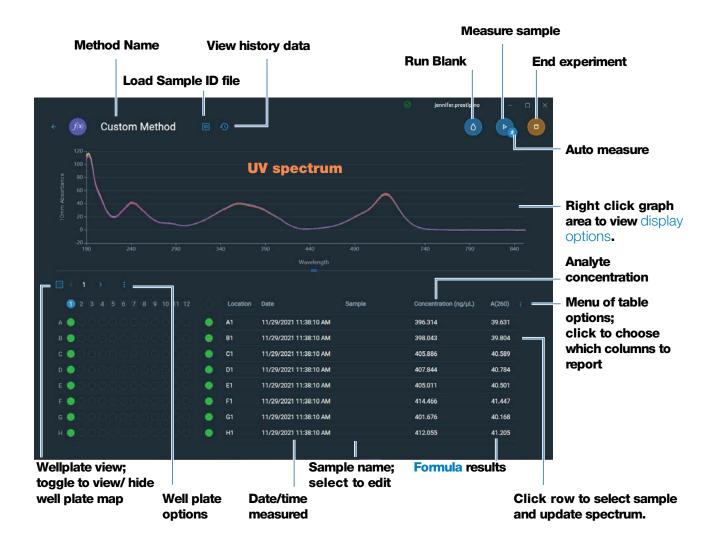

#### **Custom method reported values**

- Result name
- Measurement Range
- Analysis wavelength correction

- Factor or Extinction coefficient
- Standards (Standard curve methods only)
- Baseline correction

### **Manage Custom Methods**

The NanoDrop Eight software is your tool for creating and managing custom methods, which contain user-defined settings that can be used to acquire data with the instrument. Custom methods can be made with or without standards.

#### **Create Custom Method**

Create method to be used for sample measurements with user-defined settings.

#### **Create new custom method**

 From the NanoDrop Home screen, select the Custom tab and select Custom Methods



- Formula (if your method will not have standards)
- Standard Curve (if your method will have standards)
- In the setup window, enter Method Name
- Enter detailed **Description** of method, if desired
- Specify how to calculate and report the method results:
  - if method does not have standards, specify factor or extinction coefficient of analyte (enter "1" to report absorbance measurements only)



- if method has standards, enter name and concentration of each standard and select the curve fit type.

Inputs Standards Formula			
		Linear	
Standard ID	Concentration (ng/µL)		
Standard 1	396		
Standard 2	405.11		
Standard 3			
Standard 4			


Alternatively, load a standard curve.

- Enter or choose remaining custom settings as needed
- Choose Save

Method Description					Save	~
Inputs Formula						
Results			Analysis wavelength	340		
Concentration	ng/µL		Measurement Range			
			UV range 🗸	190 <b>nm</b> -		nr
Correction for Analysis Wavelength			Choose Factor/Extinction Coefficie			
• None			• User defined factor	10	ng/µL	
O Single point		nm	O Extinction coefficient			
O Sloping baseline	nm -	nm	Molecular weight			
Miscellanous		aseline Correc	tion 🕕 💽	Automated Pa		

**Note** Any errors in the method will be listed in red text at the bottom of the method editor screen. Errors must be addressed before the method can be saved.

#### Load standard curve

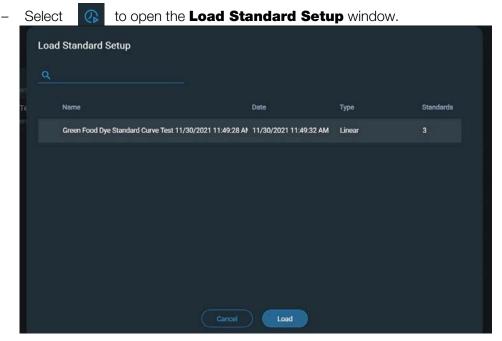

- Select a standard curve and click Load
- The measurement screen opens with the standard curve loaded. You can begin measuring samples



#### View or edit custom method

- Select **Custom Method** (existing methods are listed in Select Method box along with their type (formula or standards) and Description

select Edit

- From the Custom Method Management screen, select the method you would like to edit from the list of loaded methods.
- From the drop-down menu



- View and adjust the method settings as desired
- Select Save

#### **Custom method settings**

These settings are available for creating custom methods.

Setting	Available Options
Result name	Enter descriptive name for calculated concentration result (for example, "MTT Assay") and use adjacent drop down list to select appropriate unit. Result name appears as column heading for reported concentration value.
Measurement range	Select spectral range in which method will acquire data. Available options:
	<ul> <li>Ultra-violet only (190 nm - 350 nm)</li> </ul>
	<ul> <li>Visible only (350 nm - 850 nm)</li> </ul>
	<ul> <li>Ultra-violet and visible (190 nm - 850 nm)</li> </ul>
	<ul> <li>Custom (specify starting and ending point in nanometers)</li> </ul>
	Notes
	• If a Baseline correction and/or Analysis wavelength correction are used, make sure your selected spectral range includes your specified baseline correction and/or analysis correction wavelength.
	• For micro-volume absorbance measurements, the spectra are normalized to a 10 mm pathlength equivalent.
Analysis wavelength correction	Use this option to specify absorbance correction at analysis wavelength only. Available options:
	• <b>None.</b> No correction at analysis wavelength.
	• <b>Single point</b> . Enter wavelength for analysis correction. (Absorbance value at specified analysis correction wavelength is subtracted from absorbance value at analysis wavelength. Corrected value is used to calculate sample concentration.)
	• <b>Sloping baseline</b> . Enter two wavelengths that define sloping baseline for analysis correction. (Absorbance value of sloping baseline at analysis wavelength is subtracted from absorbance value at analysis wavelength. Corrected value is used to calculate sample concentration.)

Setting	Available Options
Factor or Extinction coefficient at 1 cm pathlength	Specify whether to use factor or extinction coefficient to calculate concentra- tion result:
(Formula methods only)	• <b>User-defined factor</b> . Enter <b>factor</b> for 1 cm pathlength and use adjacent drop down list to select appropriate <b>unit</b> . Equation below shows how factor is used to calculate sample concentration:
	c = (A * f) / b
	where: c = analyte concentration A = absorbance in absorbance units (A) $f = factor (typically 1/\epsilon, where \epsilon = wavelength-dependent molarabsorptivity coefficient, or extinction coefficient)b = pathlength$ in cm (determined at measurement time, then normalized to 10 mm (1 cm) pathlength equivalent)
	• Extinction coefficient and molecular weight. Enter extinction coefficient for 1 cm pathlength and use adjacent drop down list to select appropriate unit. Equation below shows how extinction coefficient is used to calculate sample concentration:
	$\mathbf{c} = \mathbf{A} / (\varepsilon * \mathbf{b})$
	where: c = analyte concentration A = absorbance in absorbance units (A) $\varepsilon = wavelength-dependent molar absorptivity coefficient (or extinction coefficient) b = pathlength in cm (determined at measurement time, then normalizedto 10 mm (1 cm) pathlength equivalent)$
	Notes:
	<ul> <li>Refer to product literature for information about factors and extinction coefficients for specific materials.</li> </ul>
	• To set up a method that reports absorbance measurements only, select Factor or Extinction Coefficient with the factor or extinction coefficient set to "1".
	<ul> <li>If specified unit for factor or extinction coefficient is based on mass (such as mg/mL) and specified unit for calculated result is based on molarity (such as pmol(ul) or vise verse, onter melecular weight and use</li> </ul>

(such as  $pmol/\mu L$ ) or vice versa, enter **molecular weight** and use

adjacent drop down list to select appropriate **unit**.

Setting	Available Options
Standards (Standard curve methods only)	Define the standards:
	• Enter name and analyte concentration of each standard and a reference, if desired:
	<ul> <li>Depending on the Curve Type setting, a standard curve can be generated using two or more standards. (The software allows a reference and up to 7 standards.)</li> </ul>
	<ul> <li>All reference and standards solutions should be in the same buffer used to resuspend the samples plus the same volume of reagent added to the samples.</li> </ul>
	<ul> <li>First standard can be a reference measurement. The reference solution should contain none of the analyte of interest. (The reference measurement is not the same as a blank measurement.)</li> </ul>
	<ul> <li>Concentration values for standards can be entered in any order but the standards must be measured in the order in which they were entered; however, best practice dictates that standards be measured from the lowest concentration of the standard analyte stock to the highest.</li> </ul>
	<ul> <li>Concentration range of the standards must cover the dynamic range of the assay and the expected range of the unknown samples. Sample analyte concentrations are not extrapolated beyond the concentration of the highest standard.</li> </ul>
	Select curve fit type.
	Specify type of equation used to create standard curve from standard concentration values. Available options:
	<ul> <li>Linear: Draws the linear least squares line through all measured standards (requires reference measurement and at least one standard)</li> </ul>
	<ul> <li>Interpolation: Draws a series of straight lines to connect all measured standards (requires reference measurement ans at least one standard)</li> </ul>
	<ul> <li>2<sup>nd</sup> order polynomial: Draws the 2<sup>nd</sup> order least squares polynomial using all measured standards (requires reference measurement and at least standards)</li> </ul>
	<ul> <li>3<sup>rd</sup> order polynomial: Draws the 3<sup>rd</sup> order least squares polynomial using all measured standards (requires reference measurement and at least three standards)</li> </ul>

Setting	Available Options
Analysis wavelength (Standard curve methods	Monitor absorbance at specified wavelength (enter the wavelength in nano- meters).
only)	<b>Note</b> : The specified wavelength must fall within the selected measurement range.
	The measurement results or the concentration will be calculated automatically using the absorbance value at the specified wavelength and applying the selected method type (factor or standard curve).
Baseline correction	Select this option to correct offset caused by light scattering particulates by subtracting the absorbance at a specified baseline point. Then specify wavelength for baseline correction.
	<b>Note</b> : Software subtracts absorbance value at specified baseline correction wavelength from absorbance values at all wavelengths in sample spectrum. As a result, absorbance of sample spectrum is zero at specified baseline correction wavelength.
Automated pathlength	Affects micro-volume measurements only.
	• When Automated Pathlength is selected, software selects the optimal pathlength (between 1.0 mm and 0.1 mm) based on sample absorbance at the analysis wavelength. For example, when sample absorbance at the analysis wavelength is less than or equal to 12.5 (10 mm pathlength equivalent), the optimal longer pathlength is used. When sample absorbance is greater than 12.5, the optimal shorter pathlength is used. Recommended for samples that are highly absorbing at the analysis wavelength. (This option may cause reduced sensitivity when the sample spectra have a large absorbance peak that is not at the analysis wavelength.)
	<b>Note</b> : When the analysis wavelength is between 190 nm and 219 nm, the optimal longer pathlength is used when sample absorbance is less than or equal to 10 (10 mm pathlength equivalent), and the optimal shorter pathlength is used when sample absorbance is greater than 10.
	<ul> <li>When Automated Pathlength is deselected, the software uses a 1 mm pathlength regardless of the sample absorbance. This can cause detector saturation (resulting in jagged peaks) for highly absorbing samples (e.g., ~15 A at 10 mm pathlength equivalent).</li> </ul>

Setting	Available Options				
Formula table (optional)	Use the Formula table to specify additional reported results, such as a purity ratio, for each sample.				
	Available options:				
	• <b>Predefined</b> . Select from a list of predefined formulas, which can be used as is or edited, and choose <b>Add</b> . The predefined formula is listed in the Formula Table.				
	Add. Create formula for current method. Available options:				
	• <b>Formula Name</b> . Enter a name for the formula. After a measurement, the name is reported in Data Table and Sample Details screens.				
	• <b>Formula</b> . Enter valid formula (see below for rules and examples). After a measurement, the measured or calculated value is reported in Data Table and Sample Details screens.				
	• <b>Unit</b> . Enter unit for reported result. After a measurement, the unit is reported in Data Table and Sample Details screens.				
	• Edit. Edit selected formula for current method.				
	Delete. Delete selected formula from current method.				
Formula rules	Custom formulas can include the following operators and functions:				
	Path(). Returns sample pathlength in cm.				
	• <b>A(nm)</b> . Returns sample absorbance at specified wavelength (for example, enter A(650) to add the measured absorbance at 650 nm to your equation).				
	• <b>Operators</b> : + (add), - (subtract), * (multiply), / (divide).				
	• Functions: Log(x), Pow(x,y).				
	Notes: Follow these additional rules for all languages:				
	<ul> <li>Use period "." decimal separators for floating point and double-floating point numbers.</li> </ul>				
	• Use comma "," list separators (for example, "POW(2,8)").				
	• Do not use comma "," group separators for large numbers (for example, enter 1000 rather than 1,000).				

#### **Copy Custom Method**

To create a custom method that is similar to an existing one, open the existing method, make your changes, then select **Save As** and enter a new name.

#### **Copy custom method**

- From the Custom Methods screen, select a custom method
- From the drop-down menu
   choose Edit
- Enter new Method name and Description
- Select Save As
- Enter a filename for the method and click Save

You can now select the saved method and edit the **Description** and settings.

#### **Run Custom Method**

To run a custom method, first create or import a custom method.

#### **Run custom method**

- From the Custom Methods screen, select a custom method
- Select ( ▷ Run
- Follow the on-screen instructions to measure a sample.

#### **Export Custom Method**

Export a custom method in order to run it and store the measurement results on another PC for use with another NanoDrop Eight instrument.

- From the Custom Methods screen, select a custom method
- From the drop-down menu error message is displayed; errors must be fixed before method can be exported)
- Choose Save (method is exported to method file (\*.method filename extension) in proprietary format)

#### Import custom method

You can import an existing custom method file to edit the method settings.

- From the Custom Methods screen, choose Import
- Locate and select ".method" file
- Choose Open (imported method is added to end of Select Method list)

#### **Edit custom method**

Edit a custom method in order to change the method settings.

 From Custom Methods screen, select a custom method from the list of available methods

•

- From the drop-down menu
- , choose **Edit**
- Edit method settings as desired
- Choose Save

#### **Delete custom method**

- From Custom Methods screen, select a custom method from the list of available methods
- From the drop-down menu , choose **Delete**
- After the confirmation message, choose Yes

7


# **Learning Center**


#### Contents

- Micro-Volume Sampling-How it Works 102
- Set Up the Instrument 104
- Measure a Micro-Volume Sample 106
- Sample Modes 110
- Prepare Samples and Blanks 110
- Basic Instrument Operations 116
- Acclaro Sample Intelligence 141
- Instrument Settings 147
- Measurement Screen Display Options 148

# Micro-Volume Sampling – How it Works

Surface Tension Absorbance Spectrum Sample Absorbance Sample Concentration Baseline Correction

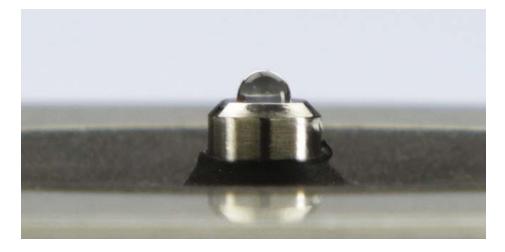


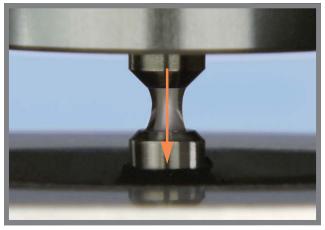

#### **Surface Tension**

The NanoDrop Eight spectrophotometer uses surface tension to hold a small volume of sample between two pedestals (upper and lower). The patented sample retention system enables the measurement of highly concentrated samples without the need for dilutions.

A fiber optic cable embedded in the upper pedestal leads to a xenon light source. A second cable embedded in the lower pedestal leads to a detector. When the instrument arm is down, the sample forms a liquid column, essentially bridging the gap between the two fiber optic cables.

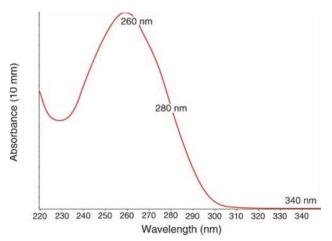

Absorbance = 
$$-\log\left[\frac{\text{intensity}_{\text{sample}}}{\text{intensity}_{\text{blank}}}\right]$$

#### **Beer-Lambert equation**

$$\mathbf{A} = \mathbf{\varepsilon} \cdot \mathbf{b} \cdot \mathbf{c}$$

where:

 $\mathbf{A}$  = absorbance in absorbance units (A)

 $\epsilon$  = wavelength-dependent molar absorptivity coefficient (or extinction coefficient) in liter/mol-cm

 $\mathbf{b}$  = pathlength in cm

**c** = analyte concentration in moles/liter or molarity (M)

#### Absorbance Spectrum

The light passes through the liquid column to the detector, which generates a spectrum of absorbance versus wavelength. The spectrum shows the amount of light absorbed by the molecules of the sample at each measured wavelength.

**Note**: To prevent evaporation, which affects measurement accuracy, close the arm quickly after you finish loading a sample or blank.

The example at the left shows a typical absorbance spectrum taken of a nucleic acid sample. The spectrum is measured from 190 nm to 850 nm. The displayed range may vary for each application.

#### Sample Absorbance

When the instrument is blanked, a reference spectrum is taken of the blanking solution and stored in memory. For each sample measurement, the sample intensities along with the blank intensities are used to calculate the total absorbance of the sample according to the equation at the left.

#### **Sample Concentration**

The Beer-Lambert equation (Beer's law) shown at the left is used to correlate sample absorbance with concentration.

The pathlength is the distance between the two pedestals, which varies in real time during each measurement. This auto-ranging pathlength technique produces accurate concentration results over a wide dynamic range.

#### **Baseline Correction**

For some applications, the instrument can be set up to apply a baseline correction to each measurement to minimize any offset caused by light scattering particulates in the sample spectra. The correction subtracts the absorbance value at a reference wavelength that is close to zero from the absorbance value at each wavelength across the spectrum, essentially "anchoring" the spectrum to zero absorbance units at the reference wavelength.

### Set Up the Instrument



### **Connect Power**




**CAUTION** Avoid shock hazard. Each wall outlet used must be equipped with a ground. The ground must be a noncurrent-carrying wire connected to earth ground at the main distribution box.

Connect the provided power cord to a grounded wall outlet. See "Power Cords" on page 172 for more information.

### **Connect to a Computer**

Connect the USB cable to the NanoDrop Eight and to an available USB port on your PC.

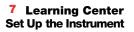

### **Operating Specifications**

The instrument operates reliably when the room environment meets these specifications:

- operating temperatures: 5 °C 35 °C (41 °F 95 °F)
- relative humidity (non-condensing): 20-80%


Locate the instrument away from air vents and exhaust fans to minimize evaporation.

**Note** If operating the instrument at the low end of the recommended humidity range, use adequate sample volume to avoid evaporation.

After the instrument is installed, you can leave it turned on.

# Measure a Micro-Volume Sample

The NanoDrop Eight spectrophotometer uses surface tension to hold a small volume of sample between two pedestals. The patented sample retention system enables the measurement of highly concentrated samples without the need for dilutions. See "Micro-Volume Sampling—How it Works" on page 102 for details.

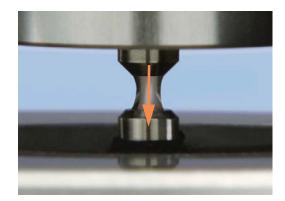

#### Supplies needed

- NanoDrop Eight spectrophotometer
- lint-free laboratory wipes
- calibrated precision pipettor
- sample material resuspended in appropriate buffer solution (see Preparing Samples)
- pure buffer solution for blanking instrument (see Choosing and Measuring a Blank.

### Best practices for micro-volume measurements

#### **Cleaning pedestals for daily operation**

- Before first measurement, clean all pedestals with a new laboratory wipe.
- Run a blanking cycle to verify pedestals are clean.
- After each measurement, clean all pedestals with new wipe to prevent carryover.
- After each set of measurements, clean pedestals with DI H2O (see Clean pedestals between users)
- Recondition pedestals periodically to maintain their hydrophobic property.

#### **Pipetting Samples**

- Position the instrument at an angle for optimal use of the pipette guide.
- Use calibrated precision pipettor (0–2 μL volume range) with well-fitting, low-retention precision tips to apply sample material to instrument for measurement. If using low accuracy (0-10 μL) pipettor, use 2 μL sample volumes.

- Use recommended sample volumes to ensure proper liquid column formation.
- Use new tip for each blank and sample aliquot.
- Use new aliquot of sample for each measurement.
- When the measurement is complete, open the sampling arm and wipe the samples from both the upper and lower pedestals using a soft laboratory wipe.

If solvents are used, make sure they are compatible with the pedestals. (see "Compatible Solvents" in Hazardous Materials).

### **Recommended sample volumes**

Application	Sample Volume
Nucleic acid (aqueous solution)	1 µL <sup>a</sup>
Purified protein	2 µL
Other protein applications such as Bradford or BCA	2 µL
Microbial cell suspensions	2 µL

<sup>a</sup> Use 2 µL for samples that contain materials that may reduce surface tension such as a surfactant.

#### To measure a micro-volume sample

#### NOTICE

- Do not use a squirt or spray bottle on or near the instrument as liquids will flow into the instrument and may cause permanent damage.
- Do not use hydrofluoric acid (HF) on the pedestals. Fluoride ions will permanently damage the quartz fiber optic cables.



 From the Home screen, select an application from one of the application categories, such as UV-Vis, or Custom Methods.

#### 7 Learning Center Measure a Micro-Volume Sample

2. Lift the instrument arm and clean the upper and lower pedestals with new laboratory wipe.





- 3. Measure a blank:
  - Pipette 1–2 µL blanking solution onto the lower pedestal and quickly lower the arm
  - Click Blank and wait for the measurement to complete


**Tip**: If Auto-Blank is On, blank measurement starts automatically after you lower the arm.


- Lift the arm and clean all pedestals with a new laboratory wipe
- 4. Measure the first sample:
  - Pipette 1-2 µL sample solution onto the pedestal and quickly lower the arm (see Recommended Sample Volumes for more information)
  - Start the sample measurement:
    - if Auto-Measure is On, lower arm
    - if Auto-Measure is off, lower arm and click



 When the sample measurement is completed, the spectra and reported values are displayed.

#### 7 Learning Center Measure a Micro-Volume Sample

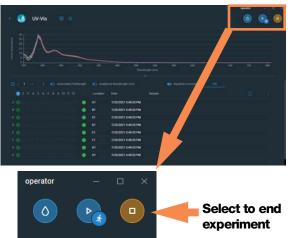

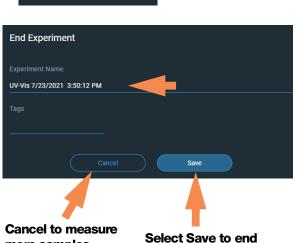

more samples

and save experiment

- 5. To measure another sample:
  - Lift the arm
  - Clean all pedestals with new wipe
  - Load the next sample(s) and quickly lower the arm
  - Start the sample measurement
  - Wait for the measurement to complete

The new spectrum replaces the previous one on the spectral display and the new reported values replace the previous ones in the table.


- 6. When you are finished measuring samples:
  - Select the End Experiment icon (see previous image)
  - Enter an experiment name or leave the default experiment name
  - Select Save
  - Lift the arm and clean all pedestals with a new wipe

If finished with the instrument for the day, clean the pedestals with DI H2O (see Clean pedestals between users)

Acquired data are automatically saved in an experiment with the entered name. In the default configuration, experiments are stored in a database according to acquisition date, experiment name, application used and any assigned labels (see Manage identifiers).

# **Sample Modes**

The NanoDrop Eight can be used in either Single sample or an Eight sample mode. Select the desired mode from the drop-down menu on the Home screen. In Single Sample mode, measurements will only be taken using Channel A.

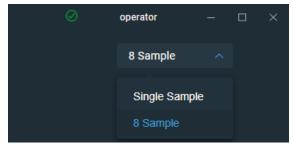

### Module startup

Instrument self-test and preparation begins upon selection of an application after first launching the software. The message 'Please wait - Initializing Spectrometer' will appear. When this message disappears, the instrument will be ready for use and you can prepare blanks. All data taken will automatically be logged in the appropriate archive file.

### Sample Plate Map

When running in Eight Sample mode, a plate map will display in the sample area. This on-screen sample plate will populate with Sample ID information from an imported sample ID file. You can also manually enter Sample ID or other identifying information. The sample status color code displayed on the plate map is determined by plate configuration at set up. See "Sample Position Illuminator" on page 126 for details.

## **Prepare Samples and Blanks**

Before making a sample measurement, a blank must be measured and stored. All eight channels are blanked with each blanking command when using the 8 sample mode. Only channel A is blanked for the single position mode.

**Note:** When using the 8 Sample mode, the software initiates each blank and measurement cycle on the first position to be read. The user will, therefore, hear one less position increment than expected. After making an initial blank measurement, a straight line will appear on the individual graphs. Subsequent blanks will clear any sample spectrum and again display straight baselines.

#### 7 Learning Center Prepare Samples and Blanks

### **Preparing Samples**

• Isolate and purify samples before measuring them with the instrument. Commercial sample isolation kits are available for these purposes, or use an in-house protocol. After purification, analyte of interest is typically dissolved in aqueous buffer solution before it is measured.

**Tip**: Any molecule that absorbs light at analysis wavelength will contribute to total absorbance value used to calculate sample concentration.

- Ensure final analyte concentration is within instrument's absorbance detection limits.
- For micro-volume measurements, gently (but thoroughly) vortex each sample before taking a measurement.


Avoid introducing bubbles when mixing and pipetting.

#### **Choosing and Measuring a Blank**

The buffer used to resuspend a sample analyte can contribute absorbance. Blanking minimizes any absorbance contribution due to the buffer components from the sample measurement. The resulting sample spectrum represents the absorbance of only the analyte of interest.

#### For best results:

- For most applications, blank with the same buffer solution used to resuspend the analyte of interest. The blanking solution should be a similar pH and ionic strength as the analyte solution. For details, see "To measure samples" in the application used.
- Measure new blank before each set of samples. It is not necessary to blank the instrument before each sample measurement unless the samples are dissolved in different buffer solutions.
- It is recommended to measure a new blank every 30 minutes.

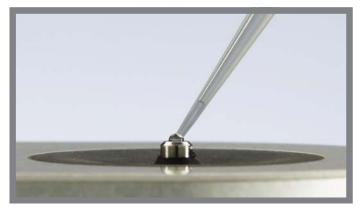

#### 7 Learning Center Prepare Samples and Blanks

 Run a blanking cycle to assess the suitability of your blanking solution before using it to perform sample measurements. For a quick demonstration, watch the multimedia training Evaluating a Blanking Solution for Suitability.

The resulting spectrum should vary no more than 0.04 A (10 mm equivalent) across the spectrum, especially at the analysis wavelength as in the example at the right.


If the resulting spectrum is greater than 0.04 A around the analysis wavelength, that buffer solution may interfere with the sample analyses, especially for low concentration samples. See below for details.

#### **Problems associated with blanking**

- Residual sample was left on pedestal before blank measurement was performed. (Resulting sample spectra may exhibit negative absorbance values, indicating blank had more absorbance than sample in that region of spectrum.)
- Blank measurement exhibits higher absorbance than unknown sample at analysis wavelength. (If buffer used as blank differs in composition from that used to resuspend sample, measurement results will be incorrect.)
- Sample was inadvertently used to blank instrument. (Resulting sample spectra may exhibit negative absorbance values or, in some cases, resemble a mirror image of a typical pure nucleic acid or protein spectrum.)



Good blanking buffer (measured abs < 0.04)

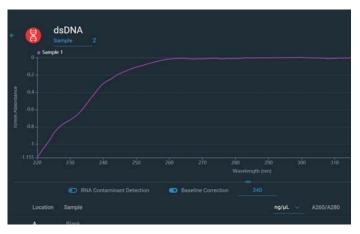

Solution containing salt used to blank instrument results in "mirror image" spectrum



#### **Solutions for blanking problems**

- Thoroughly clean and/or recondition both pedestals and then:
  - rerun blanking cycle, or
  - measure new blank using new aliquot of appropriate buffer solution, then measure new aliquot of unknown sample
- For most applications, blank with the same buffer solution used to resuspend the analyte of interest. The blanking solution should be a similar pH and ionic strength as the analyte solution. For details, see "To measure samples" in the application used.

## **Run a Blanking Cycle**

Run a blanking cycle to verify the following:

- instrument is operating normally (with flat baseline)
- pedestals are clean (i.e., no dried-down sample material on pedestals)
- absorbance contribution of buffer solution you plan to use for sample analyses

### **Supplies needed**

- lint-free laboratory wipes
- calibrated precision pipettor (0-2 μL)
- buffer solution for evaluation

## To run a blanking cycle

For quick demonstration, watch multimedia training Evaluating a Blanking Solution for Suitability.

#### NOTICE

- Do not use a squirt or spray bottle on or near the instrument as liquids will flow into the instrument and may cause permanent damage.
- Do not use hydrofluoric acid (HF) on the pedestals. Fluoride ions will permanently damage the quartz fiber optic cables.

#### 7 Learning Center Prepare Samples and Blanks


- 1. From the Home screen, select an application name.
- 2. Lift the instrument arm and clean the upper and lower pedestals with new laboratory wipe.
- 3. Measure a water blank:
  - Pipette exactly 1 µL deionized water (DI H<sub>2</sub>O) onto the lower pedestal and lower the arm.
  - Select **Blank** and wait for the measurement to complete.
  - Lift the arm and clean all pedestals with new laboratory wipe.
- 4. Measure the buffer solution:
  - Pipette 1-2 µL buffer solution onto the pedestal and lower the arm.
  - Start the sample measurement:
    - if Auto-Measure is On, lower arm
    - if Auto-Measure is off, lower arm and select **Measure**
  - Wait for measurement to complete.

The resulting spectrum should vary no more than 0.04 A from the baseline at the analysis wavelength.

If your spectrum does not meet these criteria, repeat steps 2–4.

If spectrum is still outside specifications, see Solutions for Blanking Problems.

- When you are finished with the blanking cycle, select
   End Experiment.
- 6. Lift the arm and clean all pedestals with a new wipe.

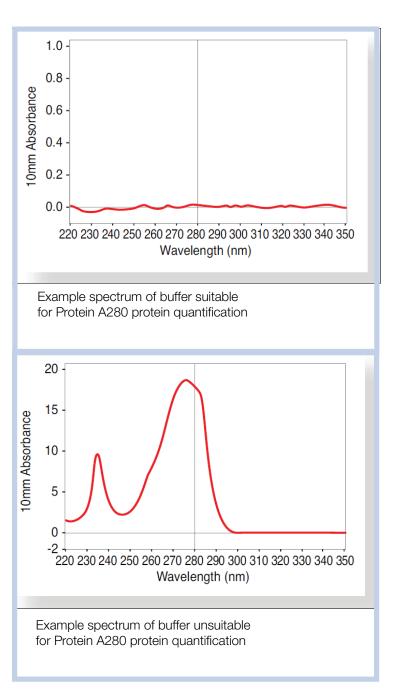


# **Basic Instrument Operations**

- NanoDrop Eight Home Screen
- NanoDrop Eight Measurement Screens
- View History
- NanoDrop Eight General Operations

# NanoDrop Eight Home Screen

These operations are available from the NanoDrop Eight Home screen.

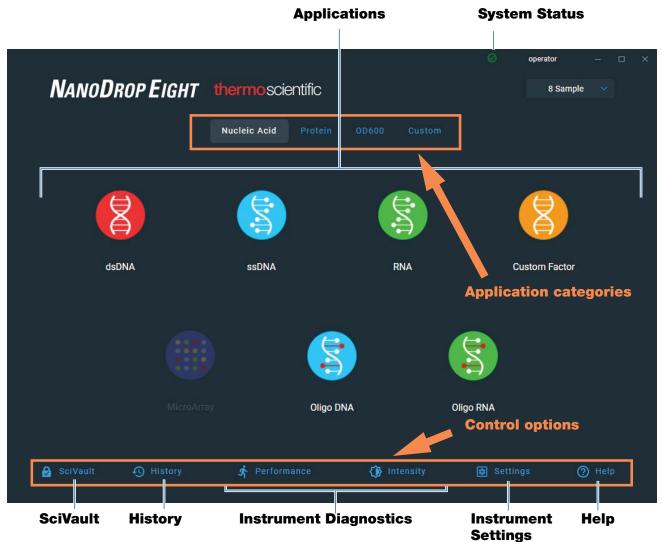

## **Applications**

The NanoDrop Eight software offers several configurable applications, which gives users full control of the measurement. See "Custom Applications" on page 81 for detailed information about each available application.

## **System Status**

Select the O icon on the Home screen to open the system status box.

thermo scientific	
NANODROP EIGHT	
Serial Number ND8_0011	
Status	Connected
Services Status	Connected
Data Storage	Local
Firmware	34
Build Number	
Software Version	1.0.0
Copyright © 2021 Thermo Fisher Scientific, Inc. All rights reserved.	
Close	

The available information is described below.

Serial number	Instrument serial number
Status	Current status of the instrument
Data storage	Indicates location of database set where instrument is currently storing data.
Firmware	Version of instrument firmware installed
Build Number	Current Software build
Software Version	Version of instrument operating software installed

### **Control options**

History:	View data stored locally. Filter by date or application.
Performance:	Performance verification process using PV-1 solution. See "Performance Verification" on page 162
Intensity:	Run an intensity check for the pedestal. See "Intensity Check" on page 160.

Settings:	Set security server location and path if desired.
Help:	View help

#### History

Select O History on the Home screen to view any data acquired earlier today, last week, last month, last six months, last year or in a specific date range. See "View History" on page 128 for more information about the History feature.

#### **Instrument Settings**

Select Settings on the Home screen to access instrument settings for software updates, Protein Editor, and more. See "Instrument Settings" on page 147 for detailed information about all available instrument settings.

#### **Instrument Diagnostics**

Instrument diagnostics (Performance and Intensity check) should be run periodically according to the recommended maintenance schedule. See "Instrument Diagnostics" on page 159 for information about how to run the available instrument diagnostics.

#### SciVault Software


The Thermo Scientific<sup>™</sup> SciVault<sup>™</sup> software is available as an optional add-on. This companion software allows users to operate their NanoDrop Eight instrument in a manner that complies with US FDA 21 CFR Part 11. When purchased, the SciVault software is shipped to you on a USB stick and integrates directly into the NanoDrop Eight software user interface. For more information, visit thermofisher.com/nanodrop.

#### Help

Select **O** Help to launch the user guide, email a question to tech support, or find the software logs.



## NanoDrop Eight Measurement Screens

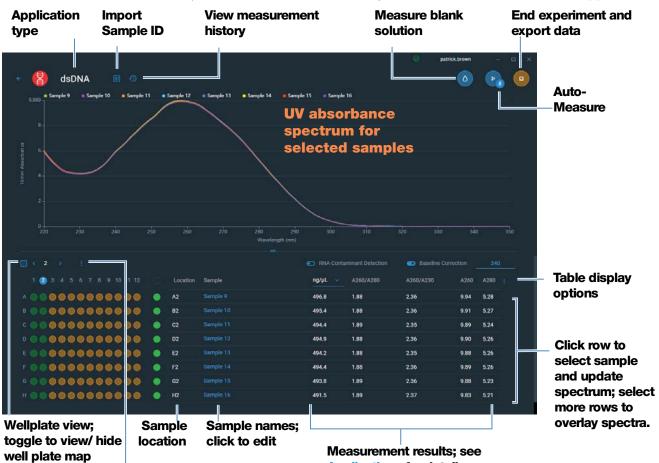

These operations are available from any measurement screen within an Application.

Well plate options

Measurement results; see Applications for details

## **Measurement Screen Display Options**


When performing measurements, right-click the graph on the measurement screen to bring up the following display options:

Overlay Mode	Display multiple samples in overlay.
✓ Show CrossHairs	Hover spectrum to see plot data
✓ Show Legend	Display graph legend
Find Peaks	Calculate peaks for specified range
Autoscale	Scale axes to fit spectra measurement
Format X-axis	Select to manually enter x-axis range
Format Y-axis	Select to manually enter y-axis range

#### Display options- Right click graph to view

## **Find Peaks**

Select **Find Peaks** to view calculated peaks for specified range. You can enter the range by dragging the color-coded limit lines, or enter values into the fields at the top of the spectrum. Found peaks for the defined range are listed in the table below the spectrum.

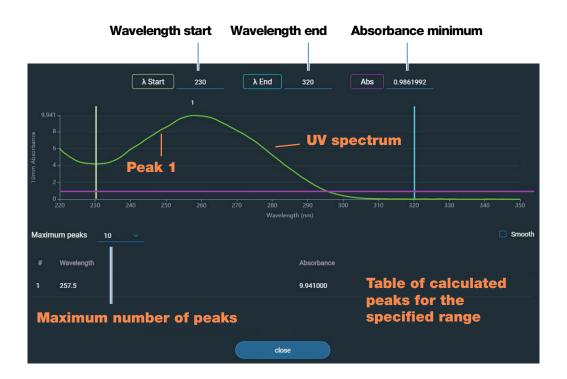



#### **Sample Name**

Click the Sample Name field in any measurement screen to edit the sample name.

In Single Sample mode, each sample is given a default base name "sample" followed by the number sample in the sequence. For example, the first sample would be named "Sample 1" followed by "Sample 2," etc. You can edit the default base name and overwrite any sample name.

In 8 Sample mode, each sample is automatically labeled by the sample location. Add sample names by clicking the sample name field, or by importing Sample ID file.

**Note** If you edit the sample base name during an experiment when Auto-Naming is selected, the assigned sample ID numbers restart.

#### Edit sample base name

After you measure a blank and before the first sample is measured:

- click Sample Name field
- enter new base name
- press Enter key

#### **Edit sample name**

- from Home screen, click 🔨 History to open History
- select experiment
- select name field of sample
- enter new sample name
- press Enter key

#### **Measurement Results**

The types of results that appear in the measurement screens depend on the selected application. For details, see the reported results section of that application in this guide:

Applications > [application group] > Measure [application name] > Reported Results

## Absorbance Spectrum

For each measured sample, each application shows the UV or UV-Visible absorbance spectrum and a summary of the results. The vertical axis shows absorbance in absorbance units (A). The horizontal axis shows wavelength in nm. Here is an example for a UV-Vis method.



#### **Sample Pathlength**

All applications display the sample pathlength along the spectrum's vertical axis. Micro-volume absorbance measurements are normalized to a 10.0 mm pathlength equivalent.

#### **Measurement Alerts**

The Acclaro Sample Intelligence technology built into the NanoDrop Eight instrument provides important features to help you assess sample integrity. click a Sample Intelligence icon in the software to view its associated information.



contaminant analysis is available to help qualify a sample before use in downstream applications



on-demand technical support is available for measurements that are atypical or very low concentrations

#### **Blank Button**

Select Blank to measure a blank for the selected experiment.





A blank must be measured before each group of similar samples. The blank solution is typically the pure buffer that was used to resuspend the sample. For more information, see Choosing and Measuring a Blank.

#### **Measure Button**

Select **Measure** to measure a sample for the selected experiment.



Samples must be properly isolated and prepared before they can be measured with the instrument and the concentration must be within the instrument's absorbance detection limits. For more information, see Preparing Samples, Measure a Micro-Volume Sample, and Absorbance Detection Limits.

**Note** The **Measure** button is enabled after a valid blank measurement is completed.

#### Auto-Measure and Auto-Blank Options

Speed up sample analysis with the NanoDrop Eight Auto-Measure and Auto-Blank features, which cause the instrument to start the measurement immediately after you lower the instrument arm. These options eliminate the need for repetitive Measure or Blank operations for large batches of samples.

**Note** Auto-Measure and Auto-Blank are available for micro-volume measurements only.

#### **Auto-Measure**

To select or deselect Auto-Measure, from any sample measurement screen, click the **On** or **Off** button at the right of the Measure button.



**Auto-Blank** 

To select or deselect Auto-Blank, from any blank measurement screen, click the **On** or **Off** button at the right of the Blank button.



## **End Experiment Button**

Select **End Experiment** when you are ready to name and save your experiment, add a label to help you locate the experiment later or export the data. Depending on the administrative settings, you may be prompted to sign the experiment upon ending the experiment.

**Note** The **End Experiment** button is enabled after the first sample measurement is completed.

After you select End Experiment, the End Experiment dialog box is displayed:

End Experime	ent	
Experiment Name		
UV-Vis 6/19/2021	3:30:09 PM	
Tags		
	Cancel Save	
	End Experiment ×	
	Experiment was saved successfully.	
	Print 👔 Export	

Click Save and you will have the option to Print or Export the experiment data.

Available options:

Experiment Name	Enter a name for this group of measurements. The measurement results are saved in the selected database location using the entered experiment name.
Tags	Enter a descriptive label to help you find this experiment later or to associate it with another experiment (see Manage identifiers on the instrument for details).

Export Select an available location for exporting the measurements in this experiment. Experiments can be exported to a USB device connected to the PC, any USB port, or to a network location. Available export file formats: comma-separated values spreadsheet (.csv) file • tab-separated values spreadsheet (.tsv) file NanoDrop Eight software (.n8db) file The filename is the entered experiment name (see above). The file is stored in a folder named "NanodropEight" followed by the instrument serial number. (Use System Status to view your instrument serial number.) Cancel (Return To Close the End Experiment box and display the results for Experiment) the most recent measurement. From there you can add measurements to the current experiment and save it later. Print button Print measurement results for current experiment Save (End Experiment) End the experiment and save the measurement results using the entered experiment name. The experiment is saved in the selected database location.

## **Sample Details**

Sample details are shown in the **sample row** in any measurement screen or History. This information includes all available measurement results and associated details for the selected sample. To add or remove results shown in the sample row, click the menu in the upper right corner of the data table. Information about the measured values displayed in Sample Details is provided in this Help system, under the application used to acquire the data.

**Note** You can also edit the sample name by clicking in the name field of the sample.

## **Data Table**

The data table for the current experiment is shown below the spectrum graph. The data table contains the measurement results for all samples in the experiment. The image below highlights the available features.

Well plate map Active channel indicator;	clie	mple name; ck to edit click ro	Ар	asurement plications f	-	e	Menu o options click to open
H • • • • • • • • • • • • • • • • • • •	<u>н</u> 2	Sample 16	491.5	1.89	2.37	9.83	5.21
G 000000000000000000000000000000000000	<b>G</b> 2	Sample 15	493.8	1.89	2.36	9.88	5.23
F • • • • • • • • • • • • • • • • • • •	F2	Sample 14	494.4	1.88	2.36	9.89	5.26
E 00000000000	🔵 E2	Sample 13	494.2	1.88	2.35	9.88	5.26
	D2	Sample 12	494.9	1.88	2.36	9.90	5.26
c •••••••••••••	O C2	Sample 11	494.4	1.89	2.35	9.89	5.24
B • • • • • • • • • • • • • • • • • • •	🔵 В2	Sample 10	495.4	1.88	2.36	9.91	5.27
A 00000000000	A2	Sample 9	496.8	1.88	2.36	9.94	5.28
1 2 3 4 5 6 7 8 9 10 11 12	Locatio	n Sample	ng/µL 🔍	A260/A280	A260/A230	A260	A280
□ < 2 > ፤			RNA Cont	aminant Detection	Baseline Co	rrection	


Data Table is in the bottom half of the measurement screen

## **Sample Position Illuminator**


The NanoDrop Eight Sample Position Illuminator and accompanying software user interface are designed to help users keep track of which samples have already been measured and which samples should be measured next. There are three different templates designed for different sample holding formats. You can choose between 96-well plate, 1.5 mL tubes, and 0.5 mL tubes by clicking the 3-dot menu above the plate matrix user interface in the bottom left of the measurement acquisition screen.

- 96-well plate: allows users to keep track of sample measurements for samples stored in a standard 96-well microtiter plate
- 1.5 mL tubes: allows users to keep track of samples measurements for samples stored in the 1.5 mL-sized spaces in the NanoDrop Eight tube rack accessory
- 0.5 mL tubes: allows users to keep track of samples measurements for samples stored in the 0.5 mL-sized spaces in the NanoDrop Eight tube rack accessory

**NOTE:** This feature works most efficiently when a user imports sample names from a file.

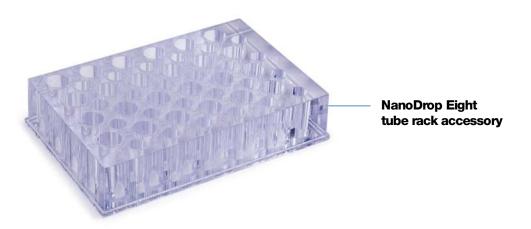

The wells/tubes which should be measured next are filled with bright orange in the software user interface. They correspond to the wells/tubes illuminated by the green LEDs on the top of the instrument. Wells/tubes that have sample IDs pre-populated will be filled with light orange. Wells/tubes that have already been measured will either be filled bright green (indicating no sample alerts) or with an A in a circle (indicating an Acclaro alert is present for that sample).

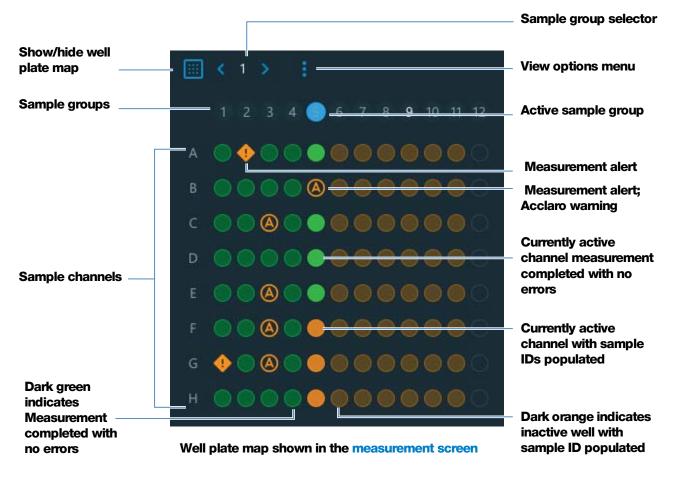


**Thermo Scientific** 

After a set of measurements is complete, users can manually advance the sample names and well illumination by clicking the sample and sarrows above the well plate map in the bottom left. This will prepare the instrument software for the next set of measurements. Alternatively, users can click the 3-dot menu sabove the well plate map and enable Auto Advance. Auto Advance sets the software to automatically advance the columns a few seconds after the last measurement is made.

The entire well plate map interface can be hidden by clicking on the plate icon

### Manual Sample ID Entry

Users can manually enter sample names using a keyboard or barcode scanner. When sample names are entered manually, the user needs to activate each channel they wish to measure. To activate/inactive all channels at once, click the circle in the column header area next to Sample. Alternatively, users can activate/inactivate individual channels by clicking the circles within each row.

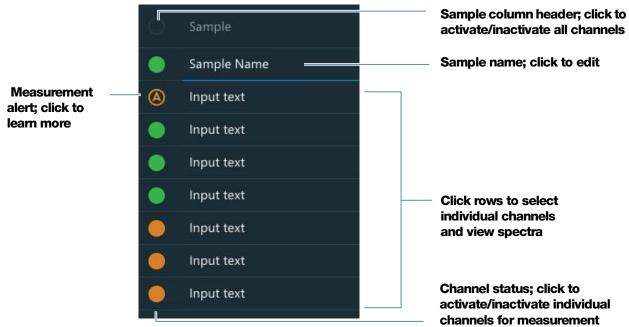



## **View History**

Whether you collect one sample or many in a row, after you choose End Experiment, the acquired data are automatically saved in an experiment with an experiment name. In the default configuration, experiments are stored in the NanoDrop Eight database according to acquisition date, experiment name, application used and any assigned labels.

Use the History feature to open the database in order to view acquired spectra and associated data from any experiment at any time.



#### **Open instrument database of measurement results**

- to open NanoDrop Eight database, from the Home screen, select History.

### **History Menu**

Available options:

Home	Return to NanoDrop Eight Home screen
Import	Import data from a USB flash drive or folder on your computer or network hard drive
Export	Export the experiment


#### Measurement Menu

When viewing measurement data from History, Click the menu in the upper right of the measurement screen to access available menu options.

Export	Export the experiment
Print	Print selected measurement results
Tags	Manage experiment tags
Details	Experiment details such as Name, Application type, Creation date, and number of measurements

### **Search Experiment Database**

Use the search feature in History to search the selected database for an experiment or to change the time range or other search filters. The database is filtered using the current settings in the Search box. Filters include time range, application type and any user-defined labels (see Manage Identifiers for information about adding and deleting labels). Here is an example:

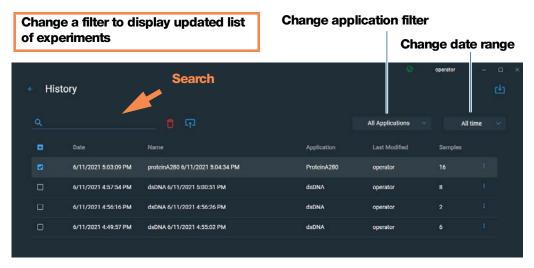

#### **Export Selected Experiments**

Use **Select** in History to select experiments to be exported.

#### **Export selected experiments**


- Open History. You can filter or use the Search feature to find experiment

+ His	tory	Search		0	operator
٩		Ō 中		All Applications	<ul> <li>✓ All tin</li> </ul>
	Date	Name	Application	Last Modified	Samples
	6/11/2021 5:03:09 PM	proteinA280 6/11/2021 5:04:34 PM	ProteinA280	operator	16
	6/11/2021 4:57:54 PM	dsDNA 6/11/2021 5:00:51 PM	dsDNA	operator	8)

- click the checkbox to select row or rows to be exported
- click Export



select one or more formats to export to (see "Export Selected Experiments")

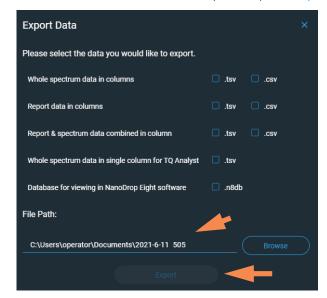

- enter or browse to an available export location USB drive, PC storage, or a network location) and select Export
- after "Export Success" message, select OK

#### **Delete Selected Experiments**

Select experiments to be deleted.

#### **Delete selected experiments**

- Filter experiments as needed, use Search feature to find desired experiment
- Click the checkbox of the **row** or rows to select one or more experiments to delete (click again to deselect an experiment)
- Click **Delete** and **OK**

Note Deleted data cannot be recovered.

#### **Open Experiment and View Associated Data**

Use History to locate and open any experiment to see the measurement data it contains.

#### **Open an experiment**


- In History, if you don't see the experiment you want to open, you can use the Search feature to find the desired experiment
- click the experiment name to open the experiment

The History provides measurement data as spectral data and data tables, similar to what you see after you complete a measurement.

**Note** The data shown are dependent upon the application used to measure the samples (nucleic acids in these examples). For more information, see the application details.

#### Spectral data -

After you open an experiment, the software shows the UV or UV-visible absorbance spectrum and associated data for the selected experiment, much like it appears during a measurement. The image below describes the available features.

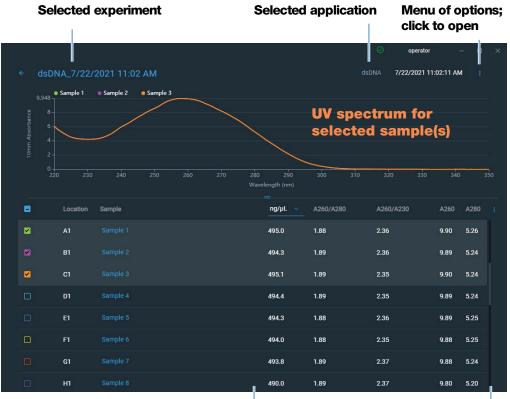

Click row to select sample and display spectrum. Select multiple rows to overlay spectra

Measurement results; see Applications for details

#### Data Table -

The lower area of the measurement history screen (shown above) shows the data table for the current experiment. The data table contains the measurement results for all samples in the experiment. The image below describes the available features.

#### Menu

Select the menu in the upper right corner from any Spectral Data or Data Table screen to see the available menu options.

Home	Return to NanoDrop Eight Home screen
Manage Identifiers	Add or delete labels for selected experiment to make it easier to find (see Manage identifiers on the instrument)
Export	Export selected experiments
Print	Print plot or data table for selected measurement results; if no results are selected, prints all results in data table
Settings	View or change instrument settings

## **NanoDrop Eight General Operations**

These operations are available from any measurement screen or from the History.

#### Manage Identifiers

You can add one or more "tags" (i.e., labels or metadata tags) to an experiment to make the experiment easier to find.

Use the History to add labels to experiments, assign existing labels, view assigned labels and remove or delete labels on the instrument. You can filter the list of experiments in the History based on one or more user-defined labels.

#### Label new experiment when you save it



End Experiment.

- after the last sample has been measured, select

- End Experiment
  Experiment Name
  UV-Vis 7/23/2021 3:50:12 PM
  Tags
  Cancel Save
- in End Experiment box, enter user defined labels in the Tags field

- click Save

### Label experiment in History

- from the Home screen, select History
- Click the menu of the experiment row and select Tags
- in Manage Tags box, enter a tag as an Identifier
- Click Save to add the tag. Add additional tags as needed.
- Click OK

#### **Remove a label**

- from Home screen, select History
- Click the menu of the experiment row and select Tags
- in Manage Tags box, click the X for any tag you wish to delete
- Click **OK**.

#### **Edit Experiment Name**

You can edit the experiment name when you save the experiment or afterwards from the History.

#### Edit experiment name at end of experiment

- when finished measuring samples, select End Experiment



End Experiment				
Experiment Name				
UV-Vis 7/23/2021 3:50:12 PM				
 Tags				
Cancel Save				

- enter a name for this group of measurements in the Experiment Name box

click Save

#### **Edit experiment name from History**

- from Home screen, open History
- use Search feature to find experiment
- Double click experiment to open experiment
- Click the **Experiment Name** field
- enter new experiment name
- the new name is saved automatically

#### **Export Selected Experiments**

You can export measurement data when you save the experiment or afterwards from the History

**Note** Data exported during a save are still saved to a database (local or remote, depending on the Data Storage setting.

Measurement data can be exported in three formats:

- as comma-separated values (.csv) files containing:
  - whole spectrum data in columns
  - report data in columns
  - both report and spectrum data combined in columns.
- as tab-separated values (.tsv) files containing:
  - whole spectrum data in columns
  - report data in columns
  - both report and spectrum data combined in columns.

- whole spectrum data in single column for TQ Analyst.
- as NanoDrop Eight software (.n8db) files containing spectra and measurement results for each exported experiment

Use any spreadsheet or word processing application to open a CSV or TSV file. Here is an example of several sample measurement results in CSV format:

1	Α	В	С	D	E	F	G	н	I I	J
1	Application:	dsDNA								
2	Serial Number:	ND8_0014								
3	User Name:	p.brown								
4	Sample Id	Sample Name	Date & Time	Location	ng/µL	A260/A280	A260/A230	A260	A280	
5	ef545555-6982-4cf1-	Sample 1	7/22/2021 11:03	A1	494.9728	1.8826	2.3554	9.8995	5.2584	
6	175af623-c7b2-4e94	Sample 2	7/22/2021 11:03	B1	494.3488	1.8854	2.3578	9.887	5.2441	
7	ef763938-31ba-4eb9	Sample 3	7/22/2021 11:03	C1	495.1052	1.8885	2.3474	9.9021	5.2432	
8	16031e24-e1d6-4628	Sample 4	7/22/2021 11:03	D1	494.3507	1.8852	2.349	9.887	5.2446	
9	a08fc6f9-bd2e-4f28-	Sample 5	7/22/2021 11:03	E1	494.3455	1.8823	2.3554	9.8869	5.2525	
10	0c88c4ce-0fdd-489e	Sample 6	7/22/2021 11:03	F1	493.9864	1.8827	2.3517	9.8797	5.2476	
11	69e3aac4-50da-4244	Sample 7	7/22/2021 11:03	G1	493.788	1.8861	2.3715	9.8758	5.2361	

**Note** The types of data exported are dependent upon the application used to measure the samples (nucleic acids in this example). For more information, see the application details.

Data can be exported to any file path connected to the PC or to a network location. If you select multiple experiments for export, each exported experiment has a corresponding file. The filenames are the same as the experiment names. The files are stored in a folder named "NanodropEight" followed by the instrument serial number. (Use System Status to view your instrument serial number.)


#### Export data at end of experiment

when finished measuring samples, click



End Experiment

- from End Experiment box, enter Tags and edit experiment name if desired, and clock Save
- Once the experiment is saved, you can click Export

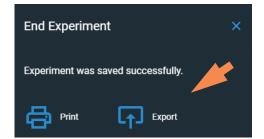

- select one or more formats to export to (see above for details)

Export Data	×	
Please select the data you would like to export.		
Whole spectrum data in columns	🗌 .tsv 🔲 .csv	
Report data in columns	🗋 .tsv 📄 .csv	
Report & spectrum data combined in column	🗋 .tsv 📋 .csv	
Whole spectrum data in single column for TQ Analyst	🗋 .tsv	
Database for viewing in NanoDrop Eight software	.n8db	
File Path:		
C:\Users\operator\Documents\2021-8-5 1221		
Export		

#### - Set File Path: to an available export path

- click **Export** 

## **Export data from History**

- From Home screen, open History
- Click the checkbox for the experiment(s) you want to export. You can use the Search feature to find experiment
- Click Export
- Select one or more formats to export to (see above for details)
- Set File Path: to an available export path and click Export

Export Data		×
Please select the data you would like to export.		
Whole spectrum data in columns	🗌 .tsv 🗌 .csv	
Report data in columns	🗌 .tsv 🔲 .csv	
Report & spectrum data combined in column	🗌 .tsv 🔲 .csv	
Whole spectrum data in single column for TQ Analyst	🗌 .tsv	
Database for viewing in NanoDrop Eight software	.n8db	
File Path:		
C:\Users\operator\Documents\2021-6-11 505		
Export		

- after "Export Success" message, click OK

#### **Delete Selected Measurements**

You can delete selected sample measurements from any experiment, or all the measurements in the database.

Note Deleted data cannot be recovered.

#### Delete data from any measurement screen

- click a sample measurement row
- click 📋

#### **Delete data from History**


- from Home screen, open History
- select **row** in History. You can filter experiments by date or type, or use Search feature to find desired experiment
- click to select one or more experiments to export
- click Delete

#### **Print Selected Measurements**

Connect a compatible printer to the PC to quickly print measurement results, including spectral data, standard curves, data tables, sample details and diagnostic results.

#### Print data from any measurement screen

- after you have measured a sample, display the measurement results to be printed such as the spectral data, standard curve, data table or sample details (see NanoDrop Eight Measurement Screens)
- from the options menu, select **Print**
- In Print Review, select the data you want to include in the data tables.

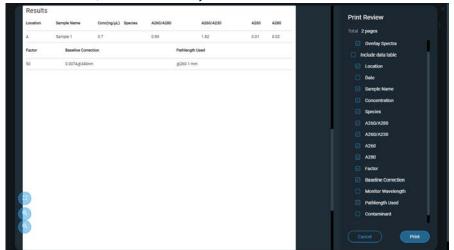

- choose Print to confirm
- in the Print Preview window, make sure the correct printer is selected and set other print options as desired such as paper size and orientation ("Auto" setting is recommended), margin and alignment to adjust the image in the preview window

**Note** The software saves the print settings each time you print.

choose Print

The selected measurement screen is printed for each selected measurement.

#### **Print data from History**

- from Home screen, open History
- click a row in History or use Search feature to find desired experiment
- click the experiment name to open the experiment
- click the menu in the upper right corner of the screen and choose Print

Location	Sample Name	Conc(ng/µL) Species	A260/A290	A260/A230	A260	A280	and the statements
A.	Sample 1	0.7	0.89	1.82	0.01	0.02	Total 2 pages
Factor	Baseline Com	ection	Puthi	ength Used			Overlay Spectra
							Include data table
50	0.007Ag340	nm .	@260	1 mitti			C Location
							Date
							Sample Name
							Concentration
							Species
							A260/A280
							A260/A230
							A260
							A280
							E Factor
							Baseline Correction
							Monitor Wavelength
9							Pathlength Used
							Contaminant
5							

- In Print Review, select the data you want to include in the data tables.

- choose **Print** to confirm
- in the Print Preview window, make sure the correct printer is selected and set other print options as desired such as paper size and orientation ("Auto" setting is recommended), margin and alignment to adjust the image in the preview window

**Note** The software saves the print settings each time you print.

- choose Print

The selected measurement screen is printed for each selected measurement.

#### 7 Learning Center Acclaro Sample Intelligence

# **Acclaro Sample Intelligence**

The Thermo Scientific<sup>™</sup> Acclaro<sup>™</sup> Sample Intelligence technology built into the NanoDrop Eight instrument provides these exclusive features to help you assess sample integrity:



contaminant analysis to help qualify a sample before use in downstream applications



on-demand technical support for measurements that are atypical or very low concentration



Use these embedded resources to quickly troubleshoot possible problem measurements and make informed decisions on whether to use, re-purify or take other actions with an atypical sample result. The Sample Intelligence feature also serves as a resource for further study and a learning tool for new or novice users.

## **Activate Detection**


From the measurement screen, click the activate contaminant detection option.

## RNA Contaminant Detection

When enabled, RNA Contaminant Detection/DNA Contaminant detection will apply mathematical models to predict the amount of RNA contaminant in dsDNA or dsDNA in RNA. These models are specific to the source of the nucleic acid. The mouse icon is representative of all mammalian sources of nucleic acid. If you are measuring nucleic acid from a source for which we do not have mathematical model, leave the selections unchecked.

## **View Acclaro Sample Intelligence Information**

Measurements that include a contaminant analysis or technical information are flagged automatically (see examples below). Select the icon to review the associated data or information.

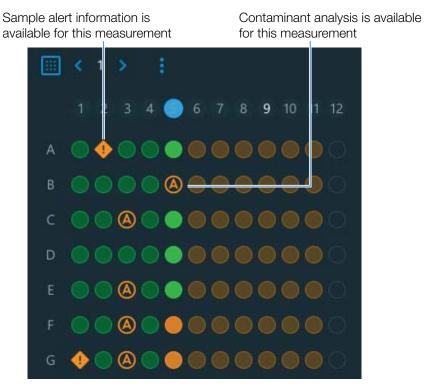

The icons appear in the Sample Plate Map, the data table, and in History (see below).

		Location	Sample	ng/µL 💊
		G	Blank	
		н	Blank	
M	٠	A1	1000 ng/µl	-0.0
	٠	B1	500 ng/µl	-0.2
	٠	C1	250 ng/µl	-0.5
		D1	125 ng/µl	-0.6

The icons are active in all three places; the information remains with the data indefinitely, even after it has been exported.

## **Contaminant Analysis**

For the dsDNA, RNA and Protein A280 applications, the NanoDrop Eight software automatically initiates a spectral analysis for several known contaminants during the measurement. Examples of known contaminants include:


- for dsDNA and RNA measurements:
  - in the analysis region: protein and phenol
  - detects presence of guanidine HCl and guanidinium isothiocyanate
  - detects species-specific dsDNA contamination in the RNA application and detects species-specific RNA contamination is in the dsDNA application
- for protein measurements:
  - in the analysis region: nucleic acids and phenol

If contaminants are identified in a sample, the "Contaminant Analysis" icon appears to the left of the measurement results.



Click the icon to view the contaminant analysis and associated information.

The following is an example of results from a nucleic acid contaminant analysis that contains enough protein contaminant to affect the measurement results.

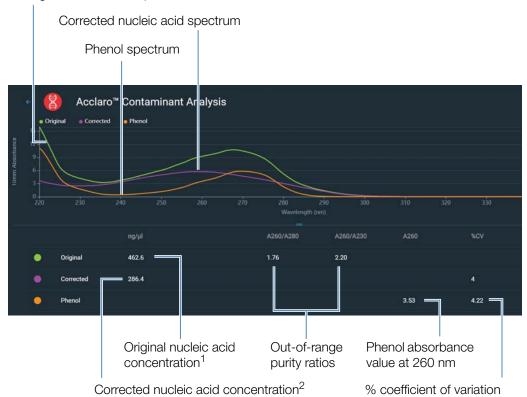

Original nucleic acid spectrum

<sup>1</sup>Based on total sample absorbance (sample plus contaminant)

<sup>2</sup>Based on corrected sample absorbance (sample minus contaminant)

Since phenols absorb light near the analysis wavelengths for nucleic acid (230 nm, 260 nm, and 280 nm), the presence of phenol in the nucleic acid sample shown above has pushed the A260/A280 and A260/A230 ratios out of range and caused the reported nucleic acid concentration to be higher than the real value. The software identifies the impurity (phenol), and reports the following:

- baseline-corrected absorbance due to phenol (2.53) at the analysis wavelength (260 nm)
- % coefficient of variation for the measurement result (uncertainty x 100/measurement result = 4.22%; a high %CV indicates the measurement result is close to the instrument detection limit or there is an interfering component)
- original nucleic acid concentration (462.6 ng/µL), which is based on the total baseline-corrected absorbance (sample plus contaminant) at the analysis wavelength
- corrected nucleic acid concentration (286.4 ng/µL), which is based on the corrected absorbance (sample minus contaminant) at the analysis wavelength

#### Theory behind contaminant analysis

UV and UV-visible absorbance measurements are used to quantify nucleic acid and protein samples at 260 nm and 280 nm, respectively. The analysis is based on the fact that the total absorbance of a mixture solution at a given wavelength is the sum of the absorbance values of each component in the mixture.

An ongoing challenge of this method is that a number of materials used in the extraction process can absorb in various regions across the spectrum. When these contaminants are present in a sample, they can interfere with the analysis by artificially inflating the absorbance at the wavelength of interest, which causes the analyte concentration to be overestimated.

Traditionally, purity ratios are used to detect the presence of contaminants that could affect downstream applications. However, purity ratios do not always provide a complete picture of possible contamination. When a purity ratio falls outside the expected range, the spectral profile is often examined qualitatively.

Our Acclaro technology applies a quantitative approach to contaminant analysis. Through sophisticated mathematical algorithms, Acclaro analyzes the spectral data to identify probable contaminants in a sample and removes any contribution due to the contaminant from the sample result. This results in a more accurate concentration value of the analyte of interest and a more quantitative analysis of the level of contamination.

Since the spectrum of a pure compound is unique to that compound, a mixture spectrum of mostly known materials that have few interactions can be mathematically broken down into its component spectra and the components identified. The contaminant analysis algorithm uses a narrow spectral region (220-285 nm) around the analysis wavelength (260 nm for nucleic acids, 280 nm for proteins) to determine any absorbance contribution from possible known contaminants (protein or nucleic acid, and phenol) that absorb in that region. The entire spectrum is analyzed to determine the presence of other possible contaminants such as guanidine HCl and/or guanidinium isothiocyanate, which are common reagents used for nucleic acid purification.

**Note** Achieving consistent, high quality contaminant analysis results is dependent on the quality of the measured sample spectra, which is dependent on the maintenance status of the instrument. For more information, see Maintenance Schedule.

## **On-Demand Technical Support**

For the dsDNA and Protein A280 applications, the NanoDrop Eight software monitors all sample measurements for the presence of contaminants or other anomalies that may affect the measurement. Examples of monitored characteristics include:


• absorbance ratios, which indicate the presence of compounds that may interfere with sample measurements (also referred to as "purity ratios"). For more information, watch the multimedia training What is a Purity Ratio?.

If sample alert information is available, where appears to the left of the measurement results.



Click the icon to view the information.

Here are results from a nucleic acid analysis for which measured purity ratios are above the expected value

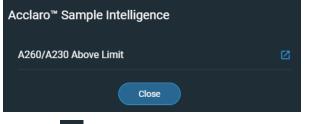

Click the **More Information** icon for the next level of information



# **Instrument Settings**

From the Home screen, select Settings. Available options include System Setings, Protein Editor, and Preferences.


## **System Settings**

These options are available:

<ul> <li>Settings System</li> </ul>	Protein Fisitor Preferences		<ul> <li>operator</li> </ul>	- ¤ ×
Database				
NanoDrop Eight Built-in Database				
Security		Language		
	Save	English		
Language	Select language f	or displaying NanoDro	op Eight so	ftware.
	<b>Notice</b> : Changin restart.	g the language require	es a softwa	are
Database	The NanoDrop Ei	ght uses a built-in dat	abase	
Security	software is installed populated for sing	ptional Thermo Scient ed. The server location gle computer installation for different compute	n is automa ons, but m	atically ust be
Auto-Export		for exported files and e end of experiments	enable/disa	able
LIMS Integration Support		orting data into LIMS ecksum into the expo		

7 Learning Center Measurement Screen Display Options

## Preferences

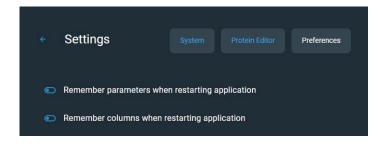

Set your preferences for remembering certain settings when restarting the application

## **Protein Editor**

See "Protein editor" on page 61 for information on using the protein editor.

# **Measurement Screen Display Options**


When performing measurements using the PC Control software, right-click the graph on the measurement screen to bring up the following display options:

Overlay Mode	Display multiple samples in overlay.
<ul> <li>Show CrossHairs</li> </ul>	Hover spectrum to see plot data
<ul> <li>Show Legend</li> </ul>	Display graph legend
Find Peaks	Calculate peaks for specified range
Autoscale	Scale axes to fit spectra measurement
Format X-axis	Select to manually enter x-axis range
Format Y-axis	Select to manually enter y-axis range

Display options- Right click graph to view

### **Find Peaks**

Select **Find Peaks** to view calculated peaks for specified range. You can enter the range by dragging the color-coded limit lines, or enter values into the fields at the top of the spectrum. Found peaks for the defined range are listed in the table below the spectrum.

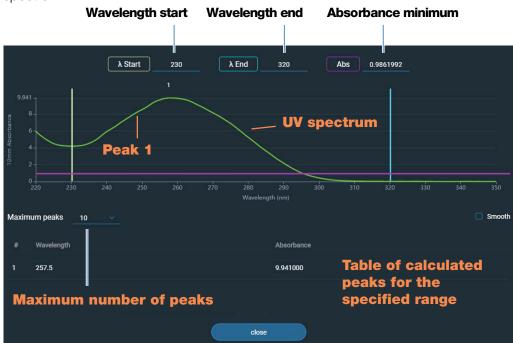

8

# **Maintenance**

- Maintenance Schedule 152
- Maintaining the Pedestals 153
- Decontaminating the Instrument 157
- Instrument Diagnostics 159

# **Maintenance Schedule**

## **Daily Maintenance**

• Clean pedestals with deionized water

# **Periodic Maintenance**

- Clean pedestals with 0.5M HCl
- Recondition pedestals



# **Every 6 Months**

- Recondition pedestals
- Run Intensity Check
- Run Performance Verification

If you are experiencing an issue with your system, refer to the troubleshooting information. If the issue persists, contact us. If you are outside the U.S.A. and Canada, please contact your local distributor.

If your instrument requires maintenance or repair, contact us or your local distributor.

# **Maintaining the Pedestals**

The pedestals require periodic maintenance to maintain measurement integrity. Time lines and procedures for cleaning and reconditioning the pedestals are provided below.

## **Cleaning the Pedestals**

To avoid carryover and cross contamination, clean the pedestals before the first blank or sample measurement and at the end of each measurement. Additional cleaning (see below) or reconditioning may be required for periodic maintenance.

#### Note

- Do not use hydrofluoric acid (HF) on the pedestals. Fluoride ions will permanently damage the quartz fiber optic cables.
- To prevent damage from spills, keep containers of liquids away from the instrument.
- Do not use a squirt or spray bottle on or near the instrument as liquids may flow into the instrument and may cause permanent damage.
- Do not attempt to remove the diaphragm around the lower pedestals as it is permanently affixed to the instrument.
- Do not allow HCI, alcohol, bleach, acetone or any other solvent to remain on the diaphragm for more than one minute or it may loosen the seals. If the diaphragm becomes loose, contact us.
- •

**Note** Solutions containing detergent or isopropyl alcohol may uncondition the pedestals. If these are required for sample analyses, follow immediately with 3– 5  $\mu$ L DI H<sub>2</sub>O.

#### 8 Maintenance Maintaining the Pedestals

#### **Supplies needed**

- lint-free laboratory wipes
- deionized water (DI H<sub>2</sub>O)
- for thorough cleaning: PR-1 kit or 0.5M HCl

# To clean the pedestals between measurements

Lift the instrument arm and clean the upper and lower pedestals with a new laboratory wipe.

# To clean the pedestals between users

- 1. Lift the arm and clean all pedestals with a new laboratory wipe.
- Pipette 3–5 μL DI H<sub>2</sub>O onto the lower pedestal.
- 3. Lower the arm and wait 2–3 minutes.
- 4. Lift the arm and clean all pedestals with a new wipe.


**Tip**: When thorough cleaning is required (for example, to remove dried sample left on the pedestals), substitute 0.5M HCl for the DI H2O in the procedure above and follow with 3-5  $\mu$ L DI H2O. You can also recondition the pedestals using PR-1 compound.




# **Reconditioning the Pedestals**

The pedestal surfaces may lose their "conditioned" properties over time, especially after measurements with isopropyl alcohol or solutions that contain surfactants or detergents such as the Bradford reagent. An unconditioned pedestal causes droplets on the lower pedestal to "flatten out," preventing proper formation of the liquid column when the arm is lowered. The resulting spectrum may look "rough" or "jagged."

If samples flatten out on the pedestal (rather than "beading up" or forming a rounded droplet) or the liquid column breaks during a measurement, recondition the pedestals.

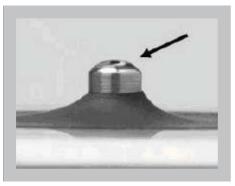

Unconditioned pedestal (droplet flattens out)

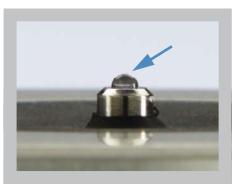

**Properly conditioned pedestal** (droplet beads up)


#### **Supplies needed**

- lint-free laboratory wipes
- PR-1 pedestal reconditioning kit (available from us or a local distributor)
- calibrated precision pipettor (0-2 μL)
- canned air

# CENTROL CENTRO

To recondition the pedestals



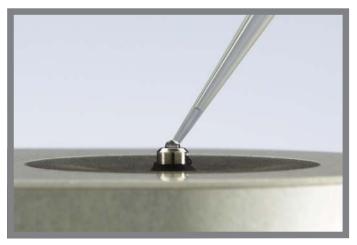

- 1. Open the container of PR-1 compound and use the provided applicator to remove a pin-head sized amount of the compound.
- 2. Apply a thin, even layer of reconditioning compound to the surface of the upper and lower pedestal.

Wait 30 seconds for the PR-1 compound to dry.


3. Fold a clean laboratory wipe into quarters and use it to vigorously buff the surface of each pedestal.

**Notice**: Support the instrument arm with one hand while you buff the upper pedestal to avoid damaging the arm.

**Tip**: Black residue on the wipe is normal.

- 4. Repeat step 3 with a new folded wipe until all residue is removed and the pedestals buff clean.
- 5. Use canned air to remove any paper residue from the pedestals.
- Pipette 1 μL DI H<sub>2</sub>O onto the lower pedestals.

The DI H<sub>2</sub>O should "bead up" or form a rounded droplet.

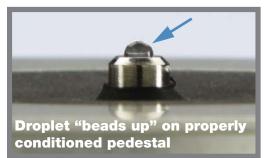

**Tip** The PR-1 pedestal reconditioning compound is the easiest way to recondition the pedestals. If you don't have a PR-1 kit, follow these steps:

- 1. Lift the instrument arm and pipette 3  $\mu$ L 0.5M HCl onto the lower pedestals.
- 2. Lower the arm and wait 2–3 minutes.
- 3. Lift the arm and clean all pedestals with a new laboratory wipe.
- 4. Pipette 3  $\mu$ L DI H<sub>2</sub>O onto the lower pedestals.
- 5. Lower the arm and wait 2–3 minutes.
- 6. Lift the arm and clean all pedestals with a new wipe.

NOTICE: Support the instrument arm with one hand while you buff the upper pedestal to avoid damaging the arm.

- 7. Fold a clean laboratory wipe into quarters and use it to vigorously buff the surface of each pedestal at least 50 times.
- 8. Use canned air to remove any paper residue from the pedestals.

# **Decontaminating the Instrument**

Decontaminate the instrument after measurements with samples that contain hazardous materials and before returning the instrument to us for maintenance or repair.

**Note** If your instrument requires maintenance or repair, contact us or your local distributor.

#### Note

- Do not use hydrofluoric acid (HF) on the pedestals. Fluoride ions will permanently damage the quartz fiber optic cables.
- To prevent damage from spills, keep containers of liquids away from the instrument.
- Do not use a squirt or spray bottle on or near the instrument as liquids may flow into the instrument and may cause permanent damage.
- Do not attempt to remove the diaphragm around the lower pedestals as it is permanently affixed to the instrument.
- Do not allow HCl, alcohol, bleach, acetone or any other solvent to remain on the diaphragm for more than one minute or it may loosen the seals. If the diaphragm becomes loose, contact us.

#### 8 Maintenance Decontaminating the Instrument

#### **Supplies needed**

- lint-free laboratory wipes
- deionized water (DI H<sub>2</sub>O)
- 0.5% sodium hypochlorite solution (1:10 dilution of commercial bleach, freshly prepared)
- pipettor

#### To decontaminate the pedestals

- 1. Lift the instrument arm and clean the upper and lower pedestal with a new laboratory wipe.
- 2. Pipette 2–3 µL diluted bleach solution (see Supplies needed) onto the lower pedestal.
- 3. Lower the arm and wait 2–3 minutes.
- 4. Lift the arm and clean all pedestals with a new wipe.
- 5. Pipette 3–5  $\mu$ L DI H<sub>2</sub>O onto the lower pedestal.
- 6. Lower the arm and wait 2–3 minutes.
- 7. Lift the arm and clean all pedestals with a new wipe.



#### To decontaminate the instrument surfaces

- 1. Dampen a clean, soft cloth or laboratory wipe with the diluted bleach solution (see Supplies needed) and use it to gently wipe the outside surfaces of the instrument.
- 2. Use a clean cloth or wipe dampened with DI  $H_2O$  to remove the bleach solution.



Diagnostics can be performed using the NanoDrop Eight software. **Intensity Check** and **Performance Verification**, are all accessible from the software Home screen:



#### Figure 1. Control options

History:	View data stored locally. Filter by date or application.
Performance:	Performance verification process using PV-1 solution
Intensity:	Run an intensity check for the pedestal
Settings:	Set security server location and path if desired
Help:	View help

### **Intensity Check**


Run Intensity Check every 6 months to verify operation of the instrument's internal components. The test measures the intensity of light from the xenon source through the instrument to verify that throughput, wavelength accuracy, and bias are within specifications. The test is automatically performed using the pedestal paths.

#### **Supplies needed**

• lint-free laboratory wipes

#### To run intensity check

- 1. From the Home screen, select **Intensity**.
- 2. Lift the instrument arm and clean the upper and lower pedestal with a new laboratory wipe.
- 3. Click **Ready** and lower the arm.
- 4. If **Auto-Measure** feature is OFF, click **Measure** to begin the measurement. If **Auto-Measure** feature is ON, the measurement will begin automatically after the arm is lowered.

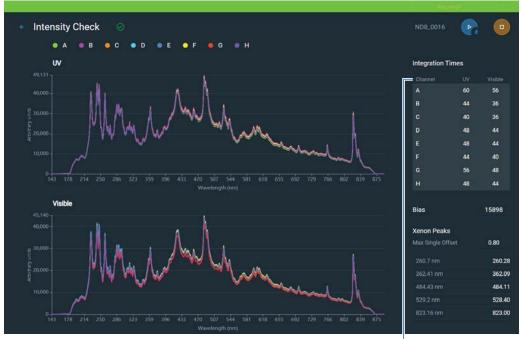

Here is an example of a typical intensity check result screen.

**Channel table** 

- 5. To rerun the intensity check, select **Run**.
- 6. When finished, select End Experiment.

After the test is completed, the results are available from the History. See Manage identifiers on the instrument for details.

#### To interpret intensity check results

The results will display the lamp intensity using the UV and Visible integration times in two separate graphs (see image above).

Users can see the intensity graphs of a single channel by clicking a single row from the Channel table or Shift + Click to highlight multiple rows (see image to the right).

A green check mark and banner will be displayed at the top if the UV intensity, Visible intensity, bias, and maximum single offset are all within specification. If one or more is not within specification, a message will be displayed at the top indicating what parameter(s) failed. If any parameter has failed, clean the pedestals with deionized water and then repeat the Intensity Check.

If a yellow triangle appears next to the Bias indicator, make sure the room is within the temperature specifications for the instrument.

If the Intensity Check fails again, contact us.

## **Performance Verification**

Run Performance Verification every 6 months to confirm pathlength accuracy is within specifications.

A CHEM-PV-8 kit containing two vials of PV-1 {aqueous nicotinic acid ( $C_6H_5NO_2$ ), potassium nitrate (KNO<sub>3</sub>)} is required to verify the performance of the Thermo Scientific<sup>TM</sup> NanoDrop<sup>TM</sup> Eight microvolume UV-Vis spectrophotometer.


#### **Supplies needed**

- lint-free laboratory wipes
- deionized water (DI H<sub>2</sub>O)
- Calibrated Precision 8-channel Pipettor (0.5 10 μL)
- CHEM-PV-8 kit containing two ampoules of PV-1, two 8-well PCR strip tubes, and transfer pipettes
- laboratory gloves

**Note** The PV-1 solution comes in a single-use ampoule. Before you open the ampoule, shake it vigorously and then allow the liquid to collect in the bottom portion of the ampoule. After the ampoule is opened, its contents must be used within one hour. Pipette directly from the ampoule; do not transfer the solution.

#### Before you begin

First make sure the pedestals are properly conditioned. To test pedestal conditioning, clean the upper and lower pedestal surfaces with a dry, lint free laboratory wipe, then pipette  $1.5 \ \mu$ L DI H<sub>2</sub>O onto the eight lower pedestal surfaces. The droplets should "bead up" as shown below. If it does not, recondition the pedestals. Remove the water sample from the upper and lower pedestal surfaces with a dry laboratory wipe.

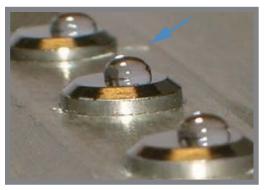

# Droplets "bead up" on properly conditioned pedestals

#### Sample Loading Hints

- Ensure the NanoDrop Eight instrument is not situated near an air vent or an exhaust fan from a nearby instrument.
- Use a small-volume (0.5 10  $\mu L)$  8-channel pipettor to load the diH\_20 and the PV-1 aliquots.

#### **Discharge and Touch Sample Delivery Method Practice**

It is very important to deliver the PV-1 aliquots to all eight pedestals in a single motion using a multichannel pipettor. It is highly recommended that the steps below be practiced using water before starting the calibration check procedure.

- 1. When the pipette tips are close to the measurement pedestals, discharge the fluid and allow the drops to hang on the end of the tips.
- 2. Gently touch the droplets to the pedestals and allow them to be pulled off the tips and onto the pedestals by surface tension.

**Tip:** It is important to use tips with a rigid structure to ensure the tips do not splay or skew when delivering the samples.

**Tip:** It is suggested that this technique be practiced until consistent delivery to all eight positions is routine and quick.

#### To run performance verification

- 1. Click on **Performance** at the bottom of the screen.
- 2. Enter Target Absorbance Value #1 (found on PV-1 ampoule label) into corresponding **Target #1 Abs** entry box.
- 3. Repeat step 2 for the **Target #2 Abs** entry box, using Target Absorbance Value #2 found on PV-1 ampoule label.

**Note** Target Abs. values are lot specific and must be entered into the correct, corresponding entry box.

- 4. Once the target values have been entered, click **Continue**.
- 5. Add 55  $\mu$ L of diH<sub>2</sub>0 to each well of one of the 8-well strips provided. Do **NOT** aliquot the PV-1 at this time.
- 6. Use an 8-channel pipettor to pipette a 1.5 iL aliquot of diH2O onto each of the lower pedestals, lower the arm, and click **Blank**.

**Tip:** It is important to use tips with a rigid structure to ensure the tips do not splay or skew when delivering the samples.

7. Remove water from upper and lower pedestal surfaces using a clean, dry laboratory wipe.

#### 8 Maintenance Instrument Diagnostics

- 8. Ensure PV-1 solution is thoroughly mixed by vigorously shaking both ampoules. Allow the solution to collect in the bottom portion of the ampoule, if needed gently tap the ampoule.
- 9. Carefully snap off top portion of each ampoule using an ampoule cracker, discard top along with ampoule cracker (use proper safety precautions for disposal).
- 10. Use the provided transfer pipette to dispense the contents of both PV-1 ampoules into each well of the 8-well PCR strip tube.
- Using an 8-channel pipettor, withdraw 1.5 µL of PV-1 solution from each of the 8- wells of the PCR strip tube and pipette onto each of the 8 lower pedestals. Lower arm.

**Note** If Auto-Measure feature is **OFF**, click Measure to begin the measurement. If Auto-Measure feature is **ON**, the measurement will begin automatically after the arm is lowered.

- 12. After the measurement is complete, remove sample from both upper and lower pedestals using a dry laboratory wipe.
- 13. Repeat steps 11 and 12 to measure five additional individual replicates of the PV-1 solution.
  - a. Always use fresh pipette tips to remove each aliquot and use a fresh aliquot of PV-1 for each measurement.
  - b. In between each measurement, remove PV-1 solution from all 16 pedestals, upper and lower, using a clean, dry laboratory wipe.
- 14. After each measurement is complete, the individual results will be displayed on screen and subsequently added to the existing results.
- 15. After six replicates have been measured, the **Status** columns for each pathlength will indicate whether the channel passed the Performance Verification.

#### To interpret performance verification results

- 1. Results will display  $\checkmark$  for Pass,  $\triangle$  for Conditional Pass, and  $\triangle$  for Fail.
  - a. If results are not within specifications, repeat the procedure using 2  $\mu L$  aliquots of PV-1.
  - b. If results fail to meet specifications using 2 µL aliquots, contact support or local distributor for assistance.
- 2. Click End Experiment when done.
  - a. Results can be exported and printed at this time or later from the History.
  - b. The Experiment name can be changed at this time and Identifiers can be added.

- 3. Click **End Experiment** when done.
- 4. To review results from a previous Performance Verification check, select **History** from the Home screen and locate the Performance Verification check results from the list of experiments.