Thermo Scientific

Heratherm Advanced Protocol and Advanced Protocol Security Microbiological Incubators

IMH 60/100/180 IMH 60-S/100-S/180-S/400-S/750-S

Operating Instructions

50125547 D

07.08.12







Contents

Chapter 1	Safety Notes	1-1
	Basic Operating Precautions	1-1
	Operational Safety Rules	1-2
	Warranty	1-2
	Explanation of Safety Information and Symbols	1-3
	Safety Notes and Symbols Used Throughout These Operating Instructions	1-3
	Additional Symbols for Safety Information	
	Symbols on the Incubator	1-5
	Intended Purpose of the Incubator	1-6
	Intended Purpose of the Incubator	1-6
	Incorrect Use	1-6
	Standards and Directives	1-6
Chapter 2	Delivery of the Incubator	2-1
•	Packaging	
	Acceptance Inspection	
	Scope of Supply	
Chapter 3	Installation	3-1
•	Ambient Conditions	3-1
	Location Requirements	
	Intermediate Storage	
	Room Ventilation	
	Space Requirement	
	Installation Built-in Units	
	Table-top incubators	
	Floor Stand Incubators	
	Transport	
	Stacking kit	
	Installing the Stacking Feet	
	Installing the Stacking Adapter	
	Installing the Anti-tilt Anchor	
	Floor stand unit spacers	
Chapter 4	Product Description	4-1
	Heratherm IMH Series Incubator Overview	
	Heratherm IMH-S Incubator Overview	
	Safety Devices	
	Work Space Atmosphere	
	Door Switch	
	Sensing and Control System	4-12

	Data Communications & Alarm Interface	4-13
	RS-232 Interface	4-13
	Alarm Contact	4-13
	AC Power Socket	4-13
	Fuses	4-13
	Work Space Components	4-14
	Inner Chamber	4-14
	Connecting Nozzles for Fresh-air Filters (Accessory Items)	4-14
	Access Port	
	Tube Access Ports	
Chapter 5	Installation procedures	5-1
•	Table-top incubators	
	Installing the Shelf System	5-1
	Initial Installation	5-1
	Installing the Perforated Shelves	
	Preparing the useable space	
	Installation or Removal of the Support Rails	
	Installing and Uninstalling the Air Baffle	
	Levelling the Incubator	
	Commissioning of floor stand incubators	
	Installing/Removing air baffles (IMH-S Series)	
	Commissioning, general	
	Installing the Shelf Support Brackets	
	Installing the Perforated Shelves	
	Mains connection	
	Connecting to the Power Supply Source	
	Connecting the RS-232 Interface	
	Interconnecting the Incubator with a Computer	
	Wiring the Alarm Contact	
	Functional Description	
	Alarm Relay Specifications	
	Connection Example	
Chapter 6	Operation	6-1
	Preparing the Incubator	
	Starting Operation	
Chapter 7	Handling and Control	7-1
•	Powering Up	
	Switching the Incubator Off / Powering Down	
	Temperature Set Value	7-7
	Timer	7-8
	Setting a Fixed-time On or Off Timer	7-10
	Setting a Weekly Timer	7-11
	Stopping a Timer	7-13
	Power Outlet (table-top incubators only)	7-15
	Decontamination	
	Fan	
	Table-top incubators	7-19

	Floor stand incubators	7-20
	Settings	7-22
	Error Log	7-22
	Calibration	
	Date and Time	7-24
	Temperature Display Unit	7-27
Chapter 8	Shut-down	8-1
	Shutting the Incubator Down	8-1
Chapter 9	Cleaning and Disinfection	9-1
·	Cleaning	
	Cleaning Exterior Surfaces	9-1
	Wipe / Spray Disinfection	9-1
	Preparing the Manual Wipe/Spray Disinfection	9-2
	Predisinfection	9-3
	Cleaning	9-4
	Final Disinfection	9-4
	Starting the Decontamination Routine	9-4
Chapter 10) Maintenance	10-1
·	Inspections and Checks	10-1
	Regular Checks	10-1
	Semi-annual Inspection	10-1
	Service Intervals	10-2
	3-monthly service	10-2
	Annual Service	10-2
	Preparing Temperature Calibration	10-3
	Comparison Measurement Procedure	10-3
	Temperature Calibration Procedure	10-4
	Replacing the Door Seal	
	Replacing the Power Cord	
	Returns for Repair	10-5
Chapter 11	l Disposal	
	Overview of Materials Used	11-1
Chapter 12	2 Error Codes	12-1
Chapter 13	3 Technical Data	13-1
Chapter 14	Spare Parts and Accessories	14-1
Chapter 15	5 Device Log	15-1
Chapter 16	6 Contact	16-1

Delivery of the Incubator

Packaging

Heratherm incubators are delivered in a rugged packaging box. All packaging materials can be separated and are reusable:

Packaging materials

Packaging carton: Recycled paper

Foam elements: Styrofoam (CFC-free)

Pallet: Chemically untreated wood

Packaging film: Polyethylene

Packaging ribbons: Polypropylene

Acceptance Inspection

After the incubator has been delivered, check the delivery immediately for:

- completeness,
- possible damage.

If components are missing or damage is found on the incubator or the packaging, in particular damage caused by humidity and/or water, please notify the carrier as well as Thermo Scientific Technical Support immediately.



Risk of injury

Should sharp edges have formed in damaged areas or elsewhere on the device, take all necessary precautions to protect personnel handling the incubator. For example, have them wear protective gloves and other personal protection equipment.

Scope of Supply

Incubators

Quantity of components supplied (pieces)	IMH Series IMH-S Series
Perforated shelves	2
Support rail for shelf table-top incubators	2
Shelf support	4
Power cord	1
Connector, potential-free contact	1
Clip springs for table-top incubators	2
Plug	1
Anti-tilt anchor	1
Operating manual	1
Short reference guide	1

2-2 Heratherm Thermo Scientific

Installation

Ambient Conditions

Location Requirements

Built-in incubators can, heating and drying ovens must be operated with an air exhaust system and exhaust hose (only original Thermo accessory should be used).

For safety reasons, the installation space should be made of non-combustible materials, according to DIN 4102.



During installation of built-in units, ensure that the escaping air will be safely discharged out of the installation space.

The incubator must only be operated in a location that meets all of the ambient condition requirements listed below:

- Installation location indoors in dry areas free from drafts.
- The dust burden may not exceed the contamination category 2 based on EN 61010-1. Using the incubator in an atmosphere with electrically conductive dust is prohibited.
- The minimal distance to adjacent surfaces must be observed on all sides (see Section "Space Requirement" on page 3-3)
- The operating room must be equipped with appropriate ventilation.
- Solid, level, fire-proof surface; no flammable materials opposite to the rear panel of the incubator.
- Vibration-proof substructure (floor stand, lab table) capable of bearing the dead weight of the incubator and its accessories (particularly if two devices are stacked).
- The electrical circuitry of the incubator has been designed for an operating height of up to 2000 m above sea level.
- Relative humidity up to 80% (maximum; preferably 60-70%), non condensing.
- Should condensation exist, wait until the moisture has evaporated completely before connecting the incubator to a power source and powering up.
- The ambient temperature must be within a range of +18 °C to +32 °C (64.4 °F to 89.6 °F).
- Avoid direct exposure to sunlight.

- Devices that produce excessive amounts of heat must not be placed near the incubator.
- Place the incubator on a floor stand (option; to be ordered separately), never on the lab floor. This prevents the penetration of dust or dirt into the device.
- Power line voltage variations must not exceed ±10 % of the nominal voltage.
- Transient surges must lie within the range of levels that normally occur in the power supply system. The impulse withstand voltage based on surge category II of IEC 60364-4-443 shall be applied at the nominal voltage level.
- Consider installing one dedicated upstream circuit breaker per incubator to avoid multiple device failures in case of an electrical fault.





Do not place the incubator directly on the lab floor, but mount it on the floor stand or on a lab work surface (option; to be ordered separately). Contaminants, such as bacteria, viruses, fungi, prions, and other biological substances may use the open door to migrate easily from the floor into the incubator's work space.

Intermediate Storage

When the incubator is placed in intermediate storage, which is permissible for a maximum of four weeks, make sure that the ambient temperature is between 20 °C to 60 °C (68 °F to 140 °F) and the maximum relative humidity does not exceed 90%, non-condensing.

Room Ventilation


Heat dissipating from the incubator during continuous operation may cause a change in the room climate.

- Therefore, the incubator must only be installed in rooms with sufficient ventilation.
- Do not install the incubator in room recesses without ventilation.
- When several devices are to be placed in the same room, additional ventilation may have to be provided as necessary.
- To avoid any impact of the heat dissipated by the incubator on the ambient climate the room must be vented by means of a laboratory-grade ventilation system that complies with applicable local and national health and safety regulations and has sufficient capacity.
- If excessive temperatures tend to occur in the operating room, be sure to provide a thermal protection means that cuts out the power supply to mitigate the impact of overtemperature scenarios.

3-2 Heratherm **Thermo Scientific**

Space Requirement

For built-in units following clearances should be kept:

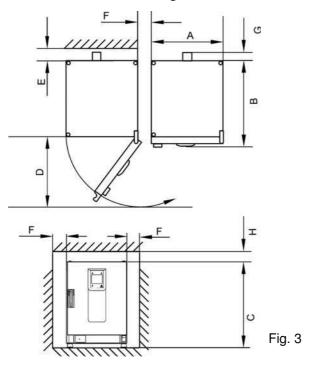

A, B, C and D see Operating Instructions of the unit.

E (mm/inch)	F (mm/inch)	G (mm/inch)	H (mm/inch)
100 / 4	50 / 2	80 / 3.2	30 / 1.2

Installation Built-in Units

IGS	Clearances without exhaust hose.
IMH/IMH-S	Clearances, close air port with plug, supplied with the unit.
OGS/OMS	Clearances, with additional space to operate to air slide, with air exhaust system and exhaust hose Ø 40 mm (1.58 in)/1.5 m (59 in) (original Thermo accessory), shortened to required length.
OGH/OGH-S OMH/OMH-S	Clearances, with air exhaust system and exhaust hose Ø 40 mm (1.58 in)/1.5 m (59 in) (original Thermo accessory), shortened to required length. Operating temperature up to max. 250 °C (482 °F).



After connecting the built-in unit to the electrical power supply, avoid damaging the power cord, then sliding into the installation position.

When installing the incubator, make sure that the installation and supply connections remain freely accessible.

The specified side clearances represent minimum distances.

Table-top incubators

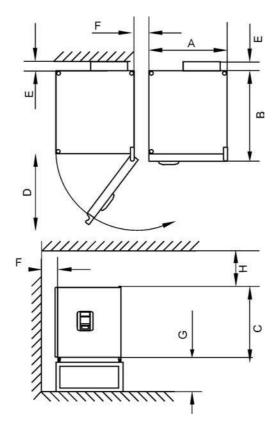


Figure 3-1 Table-top incubators, dimensions and required clearances

Table 3-1 Incubator Dimensions

Model	A (mm)	B (mm)	C (mm)	D (mm)
IMH 60	530/20.1	565/22.2	720/28.3	540/21.3
IMH 100	640/25.2	565/22.2	820/32.3	650/25.6
IMH 180	640/25.2	738/29.1	920/36.2	650/25.6
IMH 60-S	530/20.1	565/22.2	720/28.3	540/21.3
IMH 100-S	640/25.2	565/22.2	820/32.3	650/25.6
IMH 180-S	640/25.2	738/29.1	920/36.2	650/25.6

^{*} Depth of handle /display (66 mm/2.6 in) not included in overall depth specified; height of adjustable feet (36 mm/1.4 in) not included in overall height specified.

Table 3-2 Required Clearances

E (mm)	F (mm)	G (mm)	H (mm)
80/3.1	50/2.0	200/7.9	300/11.8

3-4 Heratherm Thermo Scientific

Floor Stand Incubators

400 liter units

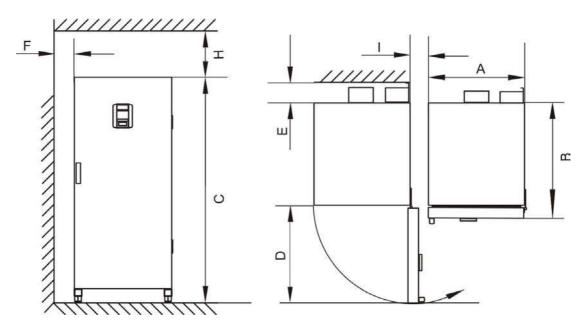


Figure 3-2 Floor stand incubators, dimensions and required clearances

Table 3-3 Incubator Dimensions

Model	A (mm/inch)	B (mm/inch)	C (mm/inch)	D (mm/inch)
IMH 400-S	755/29.7	770/30.3	1655/65.2	810/31.9

^{*} Depth of handle /display (66 mm/2.6 in) not included in overall depth specified. Width of hinge (23 mm) not included in overall width.

Table 3-4 Required Clearances

E (mm/inch)	F (mm/inch)	H (mm/inch)	I (mm/inch)
120/4.7	50 / 2	200 / 8	200/7.9

750 liter units

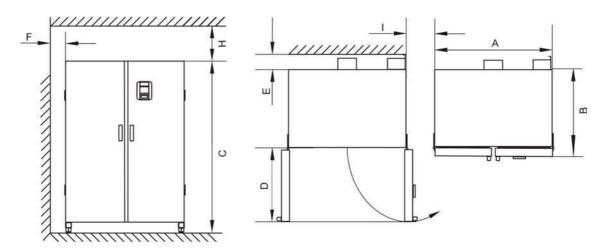


Figure 3-3 Floor stand incubators, dimensions and required clearances

Table 3-5 Incubator Dimensions

Model	A (mm/inch)	B (mm/inch)	C (mm/inch)	D (mm/inch)
IMH 750-S	1215 / 47.8	770/30.3	1655 / 65.2	670/26.4

^{*} Depth of handle /display (66 mm/2.6 in) not included in overall depth specified. Width of hinge (23 mm) not included in overall width.

Table 3-6 Required Clearances

E (mm/inch)	F (mm/inch)	H (mm/inch)	I (mm/inch)
120/4.7	50/2.0	200/7.9	350/13.8

3-6 Heratherm Thermo Scientific

Transport

Table-top incubators

For transport, do not lift the incubator using the doors or components attached to the incubator as lift points.

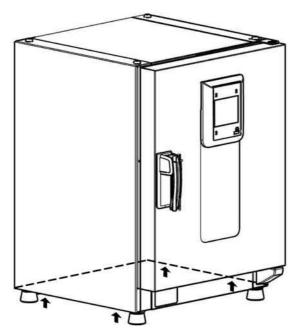


Figure 3-4 Lift Points



Heavy loads! Lift with care!

To avoid injury through physical strain, such as strain traumata and slipped discs, do not attempt to lift the incubator alone! To avoid injury through dropped loads, be sure to wear Personal Protection Equipment, such as safety shoes, when lifting the incubator.

To avoid crushing your fingers or hands (particularly in a closing door) or damaging the incubator, do not use any other lift points than those indicated in the illustration above.

Floor stand incubators



The floor stand incubators come equipped with four (4) casters. The lever for releasing the caster is located above the locking lever. After positioning the unit in its installation location ensure that the locking levers are pressed down on the casters.

To ensure the degree of stability specified by safety requirements the front casters must be turned so that they are facing forward after the unit has been positioned in its installation location and the locking levers pressed down on these casters.





Danger of tipping when moving!

Before moving the unit, ensure that it has been unplugged.

Move the Heratherm floor stand incubators with caution.

Quick starts and stops can result in tipping!

Always ensure that the doors are closed when moving the unit.

3-8 Heratherm Thermo Scientific

3-9

Stacking kit

The stacking adapter is available only for table-top incubators.

Material Number	Description
50126665	Stacking adapter Heratherm 60L
50126666	Stacking adapter Heratherm 100L
50126667	Stacking adapter Heratherm 180L

Scope of delivery:

- 1 Stacking adapter
- 1 Anti-tilt anchor
- 1 Plastic bag with 2 stacking feet and 2 M4x16 Torx screws.

Required Tools:


Slotted screwdriver 5,5x100 or Torx screwdriver 20x100.

Installing the Stacking Feet

Remove the left and right blank plugs at the top.

Install the stacking feet with the enclosed screws using a slotted or Torx screwdriver.

Installing the Stacking Adapter

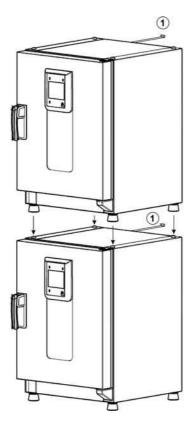

Anti-tilt anchor

Stack as follows when using a stacking adapter (numbers denote incubator volumes in liters):

- for 60 l or 100 l on 180 l use stacking adapter Heratherm 180L,
- for 60 l on 100 l use stacking adapter Heratherm100L.

To prevent the top device from slipping and dropping down, the following requirements must be fulfilled before devices may be stacked:

- Only two units may be stacked together. When stacking devices with the same type of enclosure, yet with a different footprint the device with the smaller footprint must be placed on top of the one with the larger footprint at all times.
- The bottom device must be correctly levelled.
- Place the corresponding stacking adapter.
- The levelling feet on the top device must be screwed in all the way.
- The levelling feet of the top device must be aligned with and placed exactly on the stacking pads of the stacking adapter.
- The anti-tilt anchor must be installed at the top device.


3-10 Heratherm Thermo Scientific

Installing the Anti-tilt Anchor

Table-top incubators

The anti-tilt anchor secures the top device in a stack to a solid part of a building. The anti-tilt anchor is to be mounted on the side opposite of the door hinges.

Bend the fixing tabs of the anti-tilt anchor up on one side and down on the other by an angle of approx. 90°.

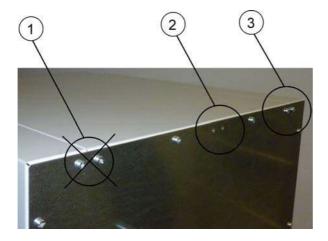

- 1. Do not use this position if the door is hinged on this side. Right-hand hinges represent the standard configuration.
- 2. Preferred position.
- 3. Alternative position. Do not use if the door is hinged on this side.

Remove the bracket screws. Use the preferred position, if possible.

Fix the anti-tilt anchor with the bracket side down to the unit.

Position the unit with the anti-tilt anchor to in an angle of approx. 90° +/- 20%.

Take care that the stacking feet of the unit are still in correct place on the lower unit or on the stacking adapter.

Fix the anti-tilt anchor to a solid part of the building.

Floor stand incubators



Heratherm floor stand incubators must always be attached to the wall using two (2) retaining brackets on the outer left and right side on the back of the unit.



Remove the screws.

Attach the end of the retaining bracket that is facing downward to the unit.

Align the device at roughly 90°, +/-20° to the retaining bracket.

Affix the retaining bracket to the wall.



Unsafe part of the building!

Install the anti-tilt anchor to a solid part of the building, which is able for shoring loads.

The installation has to be carried out by qualified personnel only.

The connection to the building must be carried out with appropriate screws and dowels according to the consistence of the building part.

Additionally, the following caution notes must be heeded at all times:



Risk of overheating with stacked devices

To avoid the risk of electrical components and the outer enclosure overheating or temperature control failing due to insufficient ventilation, do not exceed the specified stacking height!



Risk of tipping and dropping of stacked devices

You should be aware at all times that stacked devices do not form a stable unit, even when the stacking pads and frames are correctly used. The top device may tip over and drop down when being transported in a stack. To avoid injury to persons and damage to equipment, do not attempt to move stacked devices as a unit! Separate and move each device one by one, then restack them.

3-12 Heratherm Thermo Scientific

Floor stand unit spacers

The spacer on the electrical module must be pulled out and fixed in place before installing the floor stand unit at its installation location.





- 1 Loosen the 2 screws
- 2 Pull out the spacer and slide the screws into the recesses
- 3 Tighten the two screws

Product Description

This section describes Heratherm advanced protocol microbiological incubators for high-end laboratory applications, which are available in two different versions distinguished by safety level:

- Heratherm IMH Series advanced protocol microbiological fan convection incubators (see "Heratherm IMH Series Incubator Overview" on page 4-1);
- Heratherm IMH-S Series advanced protocol security microbiological fan convection incubator for more exacting safety level requirements (see "Heratherm IMH-S Incubator Overview" on page 4-4).

Heratherm IMH Series Incubator Overview

IMH Series advanced protocol microbiological incubators come equipped with the following features:

- high-precision work space temperature control, adjustable in steps of one-tenth of a degree up to 105°C (221°F)
- a variable-speed work space fan, adjustable on the control panel



At high ambient temperatures the maximum speed is reduced.

- countdown, fixed-time, and weekly timers for timed process control
- two perforated shelves
- an access port for tubing, sensor leads, etc.
- power outlet

The individual features of IMH Series incubators are shown in the figures below.

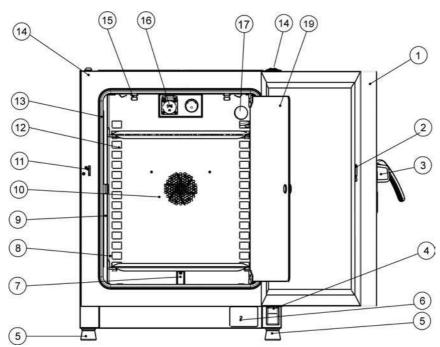


Figure 4-1 Heratherm IMH 60/ IMH 100/ IMH 180 Front View

- [1] Outer door
- [2] Door latch cutout
- [3] Door latch and handle
- [4] Door hinge, lower
- [5] Levelling foot
- [6] Nameplate
- [7] Temperature sensor
- [8] Support rail for perforated shelf
- [9] Shelf support
- [10] Fan opening, air baffle
- [11] Door hook catch
- [12] Air baffle
- [13] Door seal
- [14] Stacking pad
- [15] Spring
- [16] Power outlet
- [17] Access port
- [18] -
- [19] Glass door

4-2 Heratherm Thermo Scientific

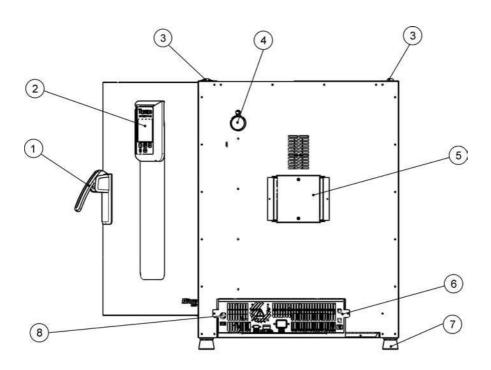


Figure 4-2 Heratherm IMH 60/ IMH 100/ IMH 180 Rear View

- [1] Door latch and handle
- [2] Control panel
- [3] Stacking pad
- [4] Access port
- [5] Fan
- [6] Electronics compartment
- [7] Levelling foot
- [8] Sample sensor connection (IMH-S only)

Heratherm IMH-S Incubator Overview

IMH-S Series advanced protocol security microbiological mechanical convection incubators come equipped with the features also found in IMH Series devices, including:

- high-precision work space temperature control, adjustable in steps of one-tenth of a degree up to 105 °C (221 °F)
- a variable-speed work space fan, adjustable on the control panel (at high ambient temperatures the maximum speed is reduced)



At high ambient temperatures the maximum speed is reduced.

- countdown, fixed-time, and weekly timers for timed process control
- two perforated shelves for sample containers
- an access port for tubing, sensor leads, etc.

Additionally, IMH-S Series incubators offer the following extra functionality:

- Monitoring of lower temperature
- Connection for Acceptable temperature sensor (accessory, must be ordered separately)
- a lockable door to secure a running process against unauthorized access
- a door switch and indicator on the control panel to indicate that the door is open
- a built-in, fully automatic decontamination routine

4-4 Heratherm Thermo Scientific

14 13 12 10 9 8 7

The individual features of IMH-S Series incubators are shown in the figures below.

Figure 4-3 Heratherm IMH 60-S/ IMH 100-S/ IMH 180-S Series Front View

- [1] Outer door
- [2] Door switch
- [3] Door latch and handle, with lock
- [4] Door hinge, lower
- [5] Levelling foot
- [6] Nameplate
- [7] Temperature sensor
- [8] Support rail for perforated shelf
- [9] Shelf support
- [10] Fan opening, air baffle
- [11] Door hook catch
- [12] Air baffle
- [13] Door seal
- [14] Stacking pad
- [15] Spring
- [16] -
- [17] Access port
- [18] -
- [19] Glass door

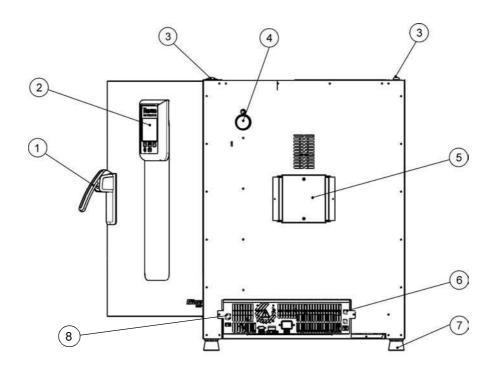


Figure 4-4 Heratherm IMH 60-S / IMH 100-S / IMH 180-S Series Rear View

- [1] Door latch and handle
- [2] Control panel
- [3] Stacking pad
- [4] Access port
- [5] Fan
- [6] Electronics compartment
- [7] Levelling foot
- [8] Sample Sensor connection

4-6 Heratherm Thermo Scientific

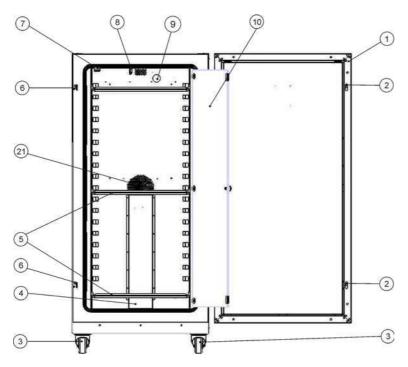


Figure 4-5 Heratherm IMH 400-S Series Front View

- [1] Outer door
- [2] Door latch
- [3] Unit caster
- [4] Air baffle
- [5] Perforated shelf
- [6] Door hook catch
- [7] Glass door latch
- [8] Temperature sensor
- [9] Access port
- [10] Glass door
- [11] -
- [12] -
- [13] -
- [14] -
- [15] -
- [16] -
- [17] -
- [18] -
- [19] *-* [20] *-*
- [21] Fan opening, air baffle

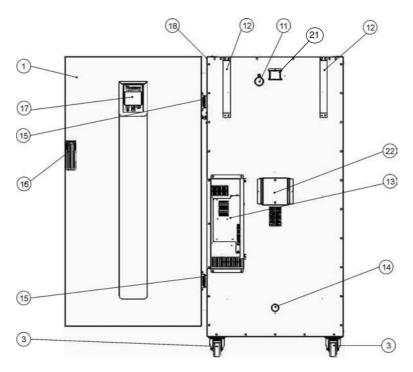


Figure 4-6 Heratherm IMH 400-S Rear View

- [1] Outer door
- [2] -
- [3] Unit caster
- [4] -
- [5] -
- [6] -
- [7] -
- [8] -
- [9] -
- [10] -
- [11] Access port
- [12] Anti-tilt anchor
- [13] Electronics compartment
- [14] Inlet air tube
- [15] Hinge
- [16] Door handle
- [17] Display
- [18] Nameplate on sidewall
- [19] -
- [20] -
- [21] Exhaust air tube
- [22] Fan

4-8 Heratherm Thermo Scientific

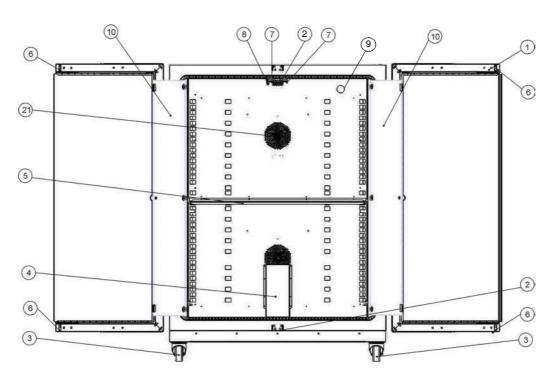


Figure 4-7 Heratherm IMH 750-S Front View

- [1] Outer door
- [2] Door hook catch
- [3] Unit caster
- [4] Air baffle
- [5] Perforated shelf
- [6] Door latch
- [7] Glass door latch
- [8] Temperature sensor
- [9] Access port
- [10] Glass door
- [11] -
- [12] -
- [13] -
- [14] -
- [15] -
- [16] -
- [17] *-* [18] *-*
- [19] -
- [20] -
- [21] Fan opening, air baffle

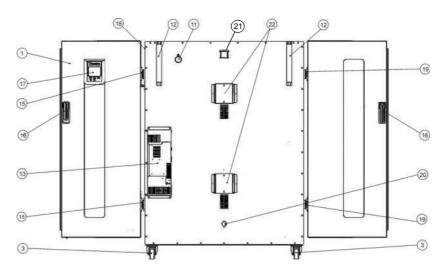


Figure 4-8 Heratherm IMH 750-S Rear View

- [1] Outer door
- [2] -
- [3] Unit caster
- [4] -
- [5] -
- [6] -
- [7] -
- [8] -
- [9] -
- [10] -
- [11] Access port
- [12] Anti-tilt anchor
- [13] Electronics compartment
- [14] -
- [15] Hinge, right
- [16] Door handle
- [17] Display
- [18] Nameplate on sidewall
- [19] Hinge, left
- [20] Inlet air tube
- [21] Exhaust air tube
- [22] Fan

4-10 Heratherm Thermo Scientific

Safety Devices

The incubators are equipped with the following safety features:

- a sample protection feature that safeguards the samples against destruction through overheating in case of contoller failure;
- an overheat protection cut-out feature that shuts down the incubators IMH-S completely when excessive temperatures occur in the workspace;
- an alarm relay that is energized, along with audible and visual alarms, to indicate errors encountered during operation;
- dual fuses rated at 16 amperes.

Work Space Atmosphere

To ensure undisturbed operation, the ambient temperature in the operating room must be at least 18 °C (64.4 °F).

The heating system uses this temperature threshold to control the ambient temperature plus 5 °C (41 °F) up to the maximum of 105 °C (221 °F).

Door Switch

Heratherm IMH 60/100/180/400/750 S incubators come with a door switch [1] integrated into the latch mechanism. If the door switch is activated by opening the door, heating operations in the work space are suspended and an icon is illuminated in the display window (see D4 in figure 7-1 on page 7-1).

If the door remains open for more than 30 seconds, a short audible alarm is sounded in addition to the icon in the display window. If the door is left open for more than 10 minutes, an audible alarm is sounded, a "door open" (E001) alarm message appears in the display and is issued through the RS-232 interface, and the alarm relay is energized.

Sensing and Control System

The PT 100-type sensor for the control of the work space temperature and for the thermal protection [1] is installed on the bottom of table-top incubators and in the top of floor stand incubators.

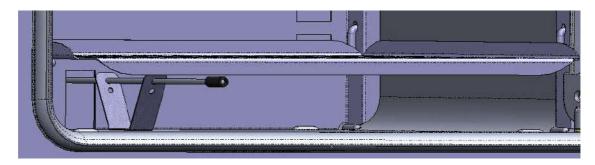


Figure 4-9 Sensor System (for table-top incubators)

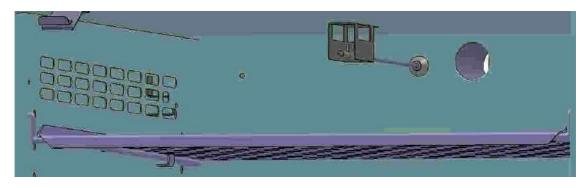


Figure 4-10 Sensor System (for floor stand incubators)

The work space temperature sensor provides the inputs to the incubator's built-in controller, which continuously compares the measured values to the user-specified set value and adjusts the heaters according to the result.



The unit features a thermal protection function that is factory-preprogrammed and not adjustable. It protects the cultures in the work space from overheating: Thermal protection kicks in on a brief violation of the upper limit, based on the defined setpoint temperature, at between 2 and 3 °C (35.6 °F and 37.4 °F) (37 °C (99 °F): 2 °C (35.6 °F), > 50 °C (122 °F): 3 °C (37.4 °F)), automatically reducing the work space temperature to the user-specified set value and allowing the incubation process to be continued even in case of a controller malfunction. If the thermal protection is activated, the error message (E111) "Temperature too high" appears in the display window and an audible alarm is sounded.

When the user acknowledges the error message, the red alarm icon (D4 in figure 7-1 on page 7-1) is illuminated and the Temperature Set Value icon (see table 7-3 on page 7-4) is highlighted by a red border to indicate that thermal protection has kicked in.


4-12 Heratherm Thermo Scientific

Data Communications & Alarm Interface

All signal connections are installed in the electrical interface panel at the rear of the incubator.

RS-232 Interface

The RS- 232 interface (item 2 in figure 4-11 below) may be used to connect Heratherm incubators to the serial interface port of a computer to allow for the computer-aided acquisition and documentation of major operating parameters (temperature, error codes, etc.).

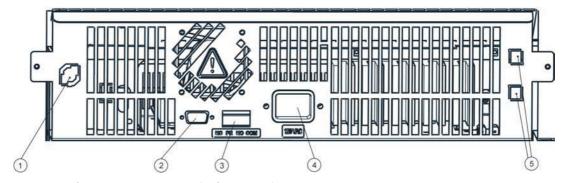

- [1] Sample Sensor connection (only IMH-S)
- [2] RS 232 interface
- [3] Alarm contact
- [4] Power socket
- [5] 2 fuses 2 A, for socket and lamp

Figure 4-11 Signal Interfaces and Power Socket

Alarm Contact

The incubator can be connected to an on-site, external alarm system (such as a private branch telephone exchange, a facility monitoring system, visual or audible alarm indicators). For this purpose, the devices come with a pre-wired potential-free alarm contact (item 3 in figure 4-11). The alarm contact is energized whenever an error occurs in an internal control loop or the incubator's electrical circuits or hardware.

AC Power Socket

The incubator is connected to the AC supply mains via the socket (item 4 in figure 4-11), which accepts a power cord with an IEC standard plug.

Fuses

Two 16 A slow-blow fuses mounted on the incubator's main electronic circuit board protect internal circuitry from the impact of excessive power consumption.



Replacement should only be carried out by skilled and authorized qualified personnel of electrotechnology/signal engineering!

NOTE

Fuse replacement

The device fuses are not user-serviceable. When the incubator exhibits the typical signs of a blown fuse (no response to pressing the On/Off button, control panel remains extinguished, no heating operation), call Thermo Fisher Scientific Customer Service to have the fuses replaced.

Work Space Components

Inner Chamber

All components of the work space are made of corrosion-resistant stainless steel and have an absolutely smooth and easy-to-clean surface. Any embossings have a large radius.

Connecting Nozzles for Fresh-air Filters (Accessory Items)

Observe the following note for IMH Series and IMH-S Series incubators with connecting nozzles for fresh-air filters (accessory item).



Connecting nozzle for fresh-air filters

The connecting nozzle for the fresh-air filter (accessory item) may neither be extended with nor be diverted through other tubing.

A fresh-air filter can be attached to the connecting nozzle. After puncturing the plastic cap the fresh-air filter can be screwed onto the nozzle.

The connecting nozzle can be closed off using the 1/8" closure plug delivered with the unit after opening the plastic cap.

The connecting nozzle for the Heratherm IMH 60/100/180 and IMH 60-S/100-S/180-S incubators has a G 1/8" threaded fitting.

4-14 Heratherm Thermo Scientific

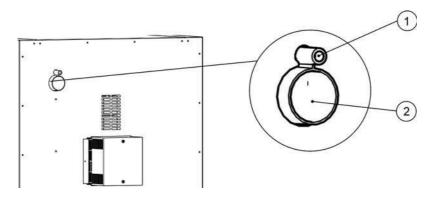


Figure 4-12 Connecting Nozzle for Fresh-air Filters (Accessory Item)

Access Port

A re-sealable, capped access port (can be closed off using the plugs delivered with the unit) (Heratherm IMH 60/100/180 and IMH 60-S/100-S/180-S/400-S/750-S only, see figure 4-7) allows cables, hoses or additional sensor leads to be routed into the work space of the incubator.


The access port [2] has a diameter of 42 mm.



When accessories are to be operated in the work space of the incubator, the ambient condition requirements must be observed (see table below). The energy introduced into the work space has an impact on the lower end of the temperature control range. When additional heating sources are introduced into the work space, temperature control may be adversely affected.

Shelf System

The incubator is supplied with two perforated shelves. The shelf support rails [1] have an alternating pattern of oblong and round perforations spaced evenly at 30 mm, allowing the shelf support brackets [8] to be inserted without any room for error, yet in a very flexible way to accommodate any required height of sample container. The shelves [2] have an integrated tilt protection and pull-out stop. For details on using the shelf system, see "Installing the Shelf System" on page 5-1.

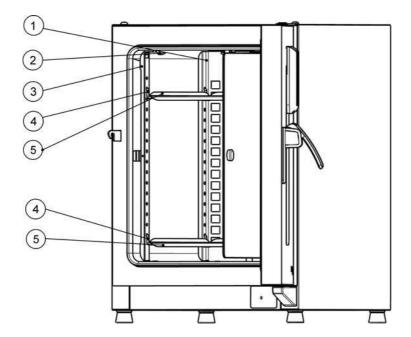

- [1] Air Baffles
- [2] Retaining Springs (only for table-top incubators)
- [3] Support Rails
- [4] Shelf Support
- [5] Shelves

Figure 4-13 Shelf System

4-16 Heratherm Thermo Scientific

Tube Access Ports

Heratherm incubators may be equipped with additional tube access ports in the side and top panels.

Available tube access port options are listed in table 4-1 below.

Tabelle 4-1 Tube Access Ports for Heratherm Incubators (Option)

Model	Side Panel Mounted Port, dia. in mm/inch	Top Panel Mounted Port, dia. in mm/inch
IMH and IMH-S	24/0.95 or 58/2.28	24/0.95 or 58/2.28

The tube access ports are mounted in fixed locations in the side and top panels (see figure 4-14).

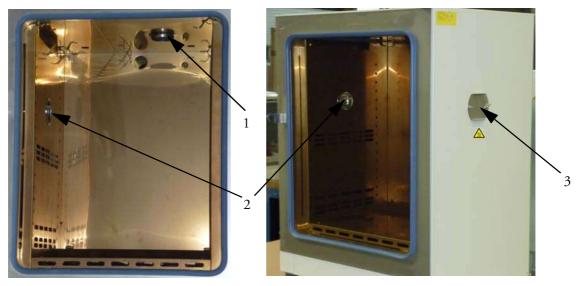


Figure 4-14 Tube Access Ports

- [1] Top panel mounted tube access port
- [2] Side panel mounted tube access port
- [3] Sealing cap for side panel mounted tube access port

Once the cables, tubes or other conduits have been inserted, the tube access ports must be padded with the heat-resistant fiber pads shipped with the device and the cap must be mounted to seal the port as far as possible.





Risk of Burning on Hot Surface

The surface around the tube access ports may become extremely hot.

Installation procedures

Table-top incubators

Installing the Shelf System

The installation of the shelf system does not require any tools. The support rails are secured in place by spring action. Once the shelf support brackets have been inserted into the rails, the perforated shelves can be simply pushed onto their support hooks to complete the installation.



Initial Installation

- 1. Peel off the protective foil from the support rails.
- 2. Push the retaining spring [1] into the guide on the support rail [2], making sure that the locking nub [3] on the retaining spring safely engages with the matching hole in the support rail.


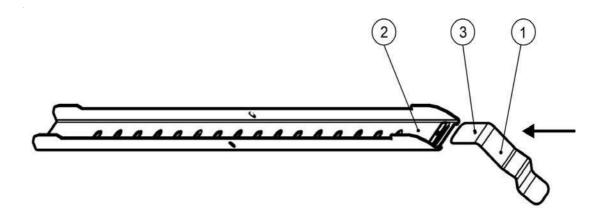


Figure 5-1 Sliding the Retaining Spring into the Support Rail

Installing the Perforated Shelves

The illustration below shows the placement of the shelf system elements.

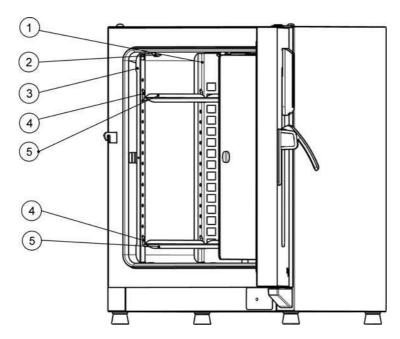

- [1] Air Baffles
- [2] Retaining Springs (only for table-top incubators)
- [3] Support Rails
- [4] Shelf Support
- [5] Shelves

Figure 5-2 Installing the Shelving

Preparing the useable space

Upon delivery, Heratherm incubators are not in a sterile state. Before the initial start-up, the incubator must be decontaminated.

The following work space components should be checked for cleanliness and disinfected prior to use:

- Support rails,
- Shelf support,
- Perforated shelves,
- Useable space surfaces,
- Work space seals and gaskets,
- Glass door

5-2 Heratherm Thermo Scientific



Cleaning and disinfection

For details about the cleaning and disinfection of the incubator, please refer to "Cleaning" on page 9-1.

Installation or Removal of the Support Rails

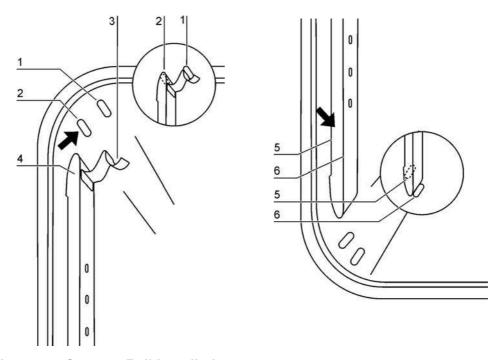


Figure 5-3 Support Rail Installation

The embossings at [2] and [5] act as lateral guides for the support rails, while the embossings at [1] and [6] secure the support rails in place. For the support rails to install correctly the retaining spring [3] must be facing upwards.

- 1. Place the support rail [4] on the lower embossing [6] and tilt it upwards against the work space side wall so that the rail is positioned over the two embossings at [5] and [2].
- 2. Clamp the retaining spring [3] behind the upper embossing [1].
- 3. To remove the support rails, pull the retaining spring tab down out of the embossing and remove the support rail assembly.

Installing and Uninstalling the Air Baffle

Heratherm IMH and IMH-S Series incubators are shipped from the factory with the air baffle readily pre-installed. Before the air baffle can be removed from the back wall, the support rails need to uninstalled.

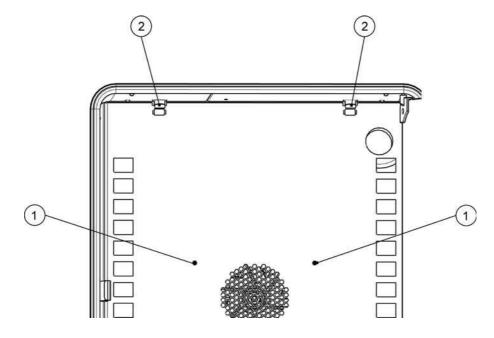


Figure 5-4 Removing the Air Baffle

- 1. Losen the two screws [1] that hold the air baffle to the interior container back wall.
- 2. Grab the two retaining springs [2] at their tabs and pull them downwards out of the embossings, then take off the air baffle from the back wall.
- 3. For the air baffle to install correctly, the retaining springs must be facing upwards. Place the air baffle on the lower embossings and tilt it upwards against the back wall of the work space.
- 4. Clamp the two retaining springs [2] into the upper embossings.
- 5. Secure the air baffle against the interior container back wall by fastening the two screws at [1].

Levelling the Incubator

- 1. Position a bubble level onto the center shelf.
- 2. Manually adjust the levelling feet until the shelf is horizontally aligned in all directions. Perform the adjustment of the levelling feet from left to right and from rear to front.

5-4 Heratherm Thermo Scientific

Commissioning of floor stand incubators

Installing/Removing air baffles (IMH-S Series)

The section below describes how to install/remove the bottom plate.

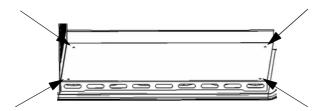


Figure 5-5 Removing the bottom plate

1. Loosen and remove the four (4) screws in the bottom plate and then remove the bottom plate completely.

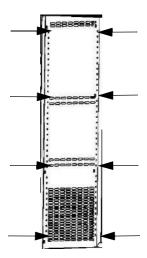


Figure 5-6 Removing the left and right support profiles

Loosen and remove the eight (8) screws for the left and right support profiles and then take out the lateral air baffles.

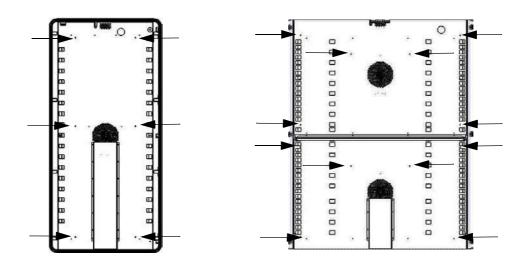


Figure 5-7 Removing the rear air baffle IMH 400-S / 750-S

On the IMH-S 400 model loosen and remove the six (6) screws for the rear air baffle and for model IMH-S 750 loosen and remove the six (6) screws for the top and bottom air baffles and then remove the air baffle(s).

Commissioning, general

Installing the Shelf Support Brackets

- 1. Insert the shelf support bracket [3] into the perforations [1] of the support rail and tilt it downwards.
- 2. Make sure that the two vertical elements [2] of the shelf support bracket butt against the support rail.


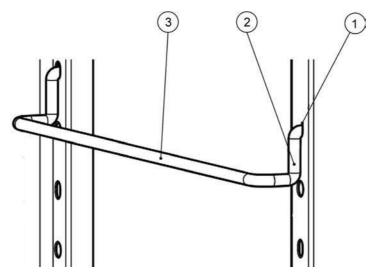


Figure 5-8 Shelf Support Bracket Installation

5-6 Heratherm Thermo Scientific

Installing the Perforated Shelves

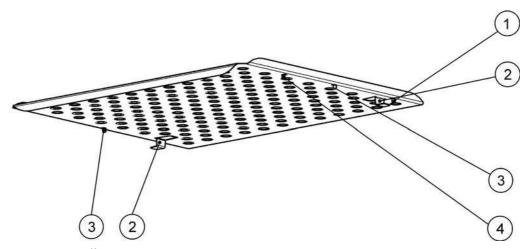

- [1] Rear Pull-out Stop
- [2] Tilt Protection Device
- [3] Front Pull-out Stop
- [4] Shelf

Figure 5-9 Perforated Shelf

- 1. Push the shelf [4] onto the shelf support brackets with the tilt protection devices [2] facing the rear panel of the incubator.
- 2. Slightly raise the perforated shelf so that the pull-out stops [1] and [3] can slide over the shelf support brackets.
- 3. Make sure that the shelves and both of their tilt protection devices are free to move over the shelf support brackets.

Mains connection





Electric shock

Contact with live electrical components may cause a lethal electric shock.

Before connecting the incubator to the power supply, check the power cord and the plug for damage. Do not use damaged cables for connecting the incubator to the power supply!

The incubator has a class I, protection-earthed enclosure. To minimize the risk of electrical shock, use the AC power cord supplied to connect the incubator to a correctly installed and protection-earthed power supply source, with the following features in place for each incubator:

- T 16 A slow-blow fusing
- B 16 circuit breaker

Connecting to the Power Supply Source

- 1. Before connecting the incubator to the power source, check to see if the power supply voltage corresponds with the specifications on the nameplate on the front of the incubator. If the voltage (V) and current (A) ratings given are not as required, do not connect the incubator to the power source!
- 2. Connect the IEC connector to the socket at the rear of the incubator.
- 3. Route the power cord along a path that does not cross exhaust air piping or passageways and aisles.
 - With stacked devices, keep the power cord away from hot spots on the other incubator in the stack.
- 4. Connect the protection-earthed plug of the power cord to a correctly protection-earthed and earth leakage circuit breaker fused power socket.
- 5. Make sure the power cord is not subjected to tensile or compressive force.



Keep the power outlet accessible!

To allow a rapid disconnection of power in case of an emergency, make sure that power outlets remain freely accessible at all times!



Figure 5-10 AC Power Supply Socket



Condensation

When taking the incubator into operation for the first time allow some time before switching on for stabilization to avoid condensation forming on live parts.

5-8 Heratherm Thermo Scientific

Connecting the RS-232 Interface



Replacement should only be carried out by skilled and authorized qualified personnel of electrotechnology/signal engineering!

The RS-232 data communication interface supports the querying of status information and temperature data from the incubator by entering basic commands in a standard terminal window provided by your computer's operating system. The interconnection requires a standard RS-232 cable with 9-pin connectors and a straight "1:1" pinout without any crossed wires, which is not supplied with the incubator.

Users may employ the RS-232 command inventory listed in table 5-1 below for automating process data logging - for example, by embedding these commands in scripts that run on a remote computer.



RS-232 interface compatibility

To avoid overloading and damaging the RS-232 interface check the interfacing parameters against the pin-out description given above and make sure that computer's interface port works with a signal level of +/-5V DC.

Interconnecting the Incubator with a Computer

- 1. Turn the computer off.
- 2. Route the serial interface cable along a path that does not cross hot exhaust air piping, tables, aisles or passageways.
 - With stacked devices, keep the serial interface cable away from hot spots on the other incubator in the stack.
- 3. Connect one connector of the serial interface cable (cable length, 5 to max. 10 m, not supplied as a standard item) to the socket labeled **RS 223** in the computer and alarm interface section at the rear of the incubator (see "Signal Interfaces and Power Socket" on page 4-13).
- 4. Connect the second connector to an unused COM 1 /COM 2 or other serial port on the computer.
- 5. Boot the computer.
- 6. Launch your standard terminal program and set up the connection with the following parameters:
 - 57600 bits per second

- 8 data bits
- 1 stop bit
- No Parity
- 7. Once your terminal indicates that serial communication has been established successfully, enter any of the commands listed in table 5-1 below, depending on what type of information you want to query.
- 8. Use the following generic command syntax:

?:aaaa:bb::cc<CR> , where:

- ?: identifies the command line as a query;
- **aaaa:** is the parameter address;
- **bb::** is a query, that must be left at "00" for technical reasons;
- cc is for a command specific checksum listed in the table below.
- **<CR>** is for carriage return.

You will receive a response of the following general format:

!:aaaa:bb:XXXXX:cc<CR> , where:

- —!: identifies the line as a response to a query;
- **aaaa:** is the parameter address entered with the query;
- **bb:** is the number of payload bytes in hexadecimal code— for example, **1F** for the decimal value **31**:s
- **XXXXXX:** is the significant status information queried;
- cc: is a check sum (technically an inverted XOR of all bytes returned, excluding the check sum bytes and the <CR> character);
- **<CR>** is for carriage return.

Table 5-1 Terminal Commands for Querying Data

Command Syntax	Response Example	
Combined Date and Time		
?:0010:00::c1	!:0010:11: 31.07.10;01:02:23 :e2 Date Time	
Date only		
?:0011:00::c0	!:0011:08: 31.07.10 :d2 Date	
Time only		
?:0012:00::c3	!:0012:08: 01:02:23 :dc Time	

5-10 Heratherm Thermo Scientific

Table 5-1 Terminal Commands for Querying Data

Command Syntax	Response Example	
Temperature Set Value (T1); Current Work Space Temperature (T2); Reference Temperature (T3); Sample Sensor Temperature (T4)		
?:3010:00::c2	!:3010:1f:+125.00;+124.96;+000.000;+000.00:b0 T1 T2 T3 T4	

Wiring the Alarm Contact



Skilled work

Thermo Scientific warrants the operational safety and the operativeness of the incubator only if installation and repairs are performed by skilled personnel.

The connection of the incubator to an external alarm system must only be carried out by adequately trained and authorized electrical engineering or telecommunications expert personnel!

Functional Description

When system errors and failures occur in the temperature control circuits, an alarm message is issued to the connected alarm monitoring system. The potential-free contact (single changeover-type contact) has been designed for the circuit configuration specified below.



Switching behavior

The alarm relay is energized by all error conditions reported by internal control loops.

Alarm Relay Specifications

Circuit	Voltage	External fusing	
Circuits with system voltage	max. 250 V ~	max. 2 A	
SELV circuits (cf.	25 V ~	max. 2 A	
VDE 0100, Part 410)	60 V =	max. 1 A	
SELV-E circuits (cf. VDE 0100,	50 V ~	max. 1 A	
Part 410)	120 V =	max. 0.5 A	



Alarm contact electrical compatibility considerations

To avoid overloading and damaging the alarm contact, check the electrical interfacing parameters of the alarm-receiving system for compatibility with the alarm relay specifications given above.

Connection Example

The connector [5] for the interface cable is supplied with the incubator as a standard item. Specifications for the operating voltage and the fusing of external alarm circuitry are given in the table above.

- 1. Wire the individual conductors [1] through [4] of the interface cable as shown in the wiring diagram.
- Route the alarm cable along a path that does not cross hot exhaust air piping, tables, aisles
 or passageways.
 With stacked devices, keep the alarm cable away from hot spots on the other incubator in
 the stack.
- 3. Plug the alarm system interface cable connector into the interface port [5] in the rear panel of the incubator.

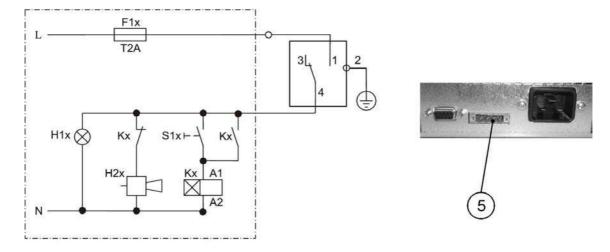


Figure 5-11 Alarm Contact Connection Example

The circuit diagram shown above represents the undisturbed condition of operation. In case of an error condition - including a power outage - contact closure occurs on the path between contacts 1-4.

5-12 Heratherm Thermo Scientific

Operation

Preparing the Incubator

The incubator must not be released for operation before all major start-up activities have been completed (see "Installation procedures" on page 5-1).

Device Check

Prior to starting operation, the following incubator components must be checked for their correct function:

- The door seal in the front frame must not be damaged.
- The glass door must not be damaged.
- The shelving components must be installed safely.
- Disinfecting the Incubator's Work Space

Run the decontamination routine (Heratherm IMH 60/100/180/400/750-S devices only; see section "Decontamination" on page 7-16) or disinfect the work space according to the operator-specified hygiene guidelines.

Starting Operation

- 1. Turn the incubator on using the control panel.
- 2. Adjust the temperature set value on the control panel.
- 3. Launch decontamination (optional with Heratherm IMH 60/100/180/400/750-S devices only; see "Decontamination" on page 7-16).
- 4. The temperature controller starts adjusting the work space to the user-specified temperature set value now.

5.





To avoid any risk of explosion or fire

- refrain from loading the incubator with any of the substances listed in the section "Incorrect Use" on page 1-6
- make sure that the ambient air is free of any solvents
- do not operate the incubator in areas with an explosion hazard
- 6. Load the work space with samples.





Hot surfaces

The screen of the glass door, the interior panel of the outer door as well as the surfaces of the shelving and the work space become hot while the incubator is running through its heating cycles and need some time to cool down.

When removing samples from a running or recently completed heating cycle, always wear safety gloves and other appropriate personal protection equipment to avoid burns on hot surfaces!



Risk of overloading

Overloading may damage the shelves or cause the shelves and/or the incubator to tilt when the shelves are being drawn out, ultimately destroying the samples. To avoid overloading the incubator or its shelving be sure to observe the sample weight limits specified in chapter 13, "Technical Data."



Proper loading

To ensure sufficient air circulation and uniform heating of the samples, do not use more than 70% of the maximum surface area of the work space. Bulky objects in the work space that dissipate heat may impair heat distribution.

6-2 Heratherm Thermo Scientific

Handling and Control

Heratherm IMH Series and IMH-S Series incubators come with a front panel mounted control unit consisting of a multifunctional display, four control buttons, and an on/off button. The four control buttons interact with the display window to let users access all of the user control functions and adjustments of the incubator, including - for example, the temperature set value, timer, energizing/de-energizing the AC outlet in the work space, as well as a variety of other functions.

Under normal operating conditions the display presents user with the work space temperature. The display returns to its default mode upon completion of the adjustments or whenever no entries have been made for a period of 30 seconds.

The graphic below shows the Heratherm IMH 60/100/180 and IMH 60/100/180/400/750-S control panel with all of its visualization elements and controls.



Figure 7-1 Control Panel for Heratherm IMH Series and IMH-S Series Incubators.



The table below contains brief descriptions of the buttons on the control panel (items K1 through K5 in figure 7-1).

Table 7-1 Control Buttons

Icon	Item	Function
MENU	K1	Menu/Enter button First key press: Activates the menu, highlighting the first menu item with a red border (see D3). Second key press: Selects the currently activated menu item (as highlighted by the red border). Depending on the currently selected function, pressing this button enables entries with item D2, D5 or D6. Third key press (once a setting has been changed): Confirms a previous entry or selection.
«	K2	After the first press of the Menu/Enter button: - Moves the selection in the menu (see item D3) to the next icon on the left. Once a menu item has been selected: - Decreases an adjustable parameter value - for example, the temperature set value in D5 or the fan speed level at D6. Holding this button depressed for a few seconds changes the selected value in quick run mode. - Moves the selection in the multifunctional display pane in D2 in the currently activated menu item to the next option on the left for example, from the Off state of the timer to On.
	K3	On/Off Button Holding this button depressed for 2 seconds switches the incubator off. The display window goes out, except for the readiness indicator icon in the status display area at item D4. The temperature display field D1 provides a dimmed readout of the work space temperature, provided that the temperature exceeds 50 °C (122 °F).
>	K4	Right button After the first press of Menu/Enter button: - Moves the selection in the menu (see item D3) to the next icon on the right. Once a menu item has been selected: - Increases an adjustable parameter value - for example, the temperature set value in D5 or the fan speed level at D6. Holding this button depressed for a few seconds changes the selected value in quick run mode Moves the selection in the multifunctional display pane in D2 in the currently activated menu item to the next option on the left for example, from the Off state of the timer to On.
ESC	K5	Escape button Returns to the previous level of the menu or standard display. Upon exiting from the current menu item the user may be prompted to save any previously made settings.

The table below contains brief descriptions of the display features of the control panel (items D1 through D6 in figure 7-1; the identifiers K1 through K4 refer to the buttons shown in that figure).

7-2 Heratherm Thermo Scientific

Table 7-2 Display Features

Feature	Item	Function
2 4.0 _°	D1	Display field showing a permanent readout of the actual temperature in the work space either in °C or °F (depending on the user's preferences, see "Temperature Display Unit" on page 7-27). At temperatures below 105 °C or 221 °F the temperature readout has one digit after the decimal point, while temperatures beyond are shown without any decimal places. Alternatively, a flashing time entry prompt of the general format hh:mm (hours:minutes, both with two digits) appears in this place while the user is setting the incubator's built-in clock.
2010-03-29 12:59 pm 4 On Off) 29.03.2010 12:59 heating relay error (E109)	D2	Four-line multifunctional display pane with fields for date and time, a display area for the specific options of the selected menu item, alarm codes, progress indicators for continuous processes (for example, program-controlled temperature ramping), etc.
	D3	Menu bar with iconized representations of adjustable parameters. A red border is used to highlight the current menu item, as selected using the Menu (K1) and arrow buttons Left (K2) and Right (K4). Brief descriptions of the individual menu items are given in table 7-3 below. Note If a menu item cannot be selected, then the function it represents is not part of the equipment configuration of your unit.
	D4	Status display area with three icons representing specific statuses of the incubator (from left to right): - The Door Open icon appears when the front door of the incubator is open or has not been closed correctly (see "Door Switch" on page 4-11). Note The Door Open icon is only functional with IMH-S Series devices. - Upon occurrence of an error condition, the red alarm icon will be illuminated. At the same time the current error code will flash in the display field at D2. The alarm may be acknowledged by pressing the button. - The readiness indicator icon appears when the incubator has been switched off using the On/Off button (item K3 in figure 7-1).
Set 889.818	D5	Settings pane labeled Set for temperature set value in either °C or °F (depending on the user's preferences; see "Temperature Display Unit" on page 7-27). At temperatures below 105 °C or 221 °F the temperature readout has one digit after the decimal point, while temperatures beyond are shown without any decimal places.

Table 7-2 Display Features

Feature	Item	Function
	D6	The left one of the two vertical bar graphs belongs to the Fan icon directly beneath it and displays the current fan speed level.
		Fan speed level for table-top incubators Bar graph for fan speed setting (in 6 steps: 0 – 5) - 0% (no chevron illuminated) - 20% (chevron 1 illuminated) - 40% (chevrons 1 and 2 illuminated) - 60% (chevrons 1 through 3 illuminated) - 80% (chevrons 1 through 4 illuminated) - 100% (chevrons 1 through 5 illuminated)
(((((((((((((((((((((Fan speed level for floor stand incubators Bar graph for fan speed setting (in 2 steps) Minimum fan speed level (chevrons 1 through 3 illuminated) Maximum fan speed level (chevrons 1 through 5 illuminated)

The table below contains brief descriptions of the menu bar icons (item D3 in figure 7-1).

Table 7-3 Menu Bar Icons

lcon	Function
	Temperature Set Value Allows for changing the temperature set value (factory-preset to 37 °C/99 °F) within the permissible temperature range. The set value can be changed by pressing the Left and Right (item K2 or K4) and you can, after confirming your changes with the Menu/Enter button (item K1), track the impact on the actual temperature in the temperature display field at D1. Instructions: "Temperature Set Value" on page 7-7.
	Timer Allows for having the incubator turn on or off upon expiry of a user-specified countdown period or at a fixed on or off time, or having it operate on a complete weekly schedule of daily on and off times. When the user enables an "on timer" the display goes out. A rotating hand in the Timer icon and the illuminated readiness indicator icon in the status display area indicates that the timer is running. Instructions: "Timer" on page 7-8.
#	Power Outlet (IMH Series only, table-top incubators only) Toggles the AC outlet in the work space of the incubator on and off. Instructions: "Power Outlet (table-top incubators only)" on page 7-15.

7-4 Heratherm Thermo Scientific

Table 7-3 Menu Bar Icons

Function Icon **Decontamination (IMH-S Series only)** Activates the built-in decontamination routine. **Instructions:** "Decontamination" on page 7-16. Fan Turns the fan on and allows for choosing the fan speed levels described under D6. The current setting is shown by the bar graph located directly above the icon and spelled out as a numeric value in the display pane at D2. Instructions: "Fan" on page 7-19. **Settings** Invokes a submenu with the following functions: - Read access to error log - Calibrating the incubator - Setting date and time - Toggling the temperature display unit between °C and °F - Selecting the timer mode of operation (countdown / fixed time of day / weekday timer) - Entering a configuration control code (Instructions: "Settings" on page 7-22)

Powering Up

1. Plug the power plug of the incubator into a suitable protection-earthed AC power outlet.



In the display window on the front panel the readiness indicator icon (rightmost icon in the status display area at D4 in figure 7-1 on page 7-1) is illuminated.



2. Keep the **On/Off** button depressed for two seconds.

An initialization routine will be run after the incubator has been powered up. On completion of the initialization, the display will light up and the current work space temperature will appear in the temperature display field (item D1 in figure 7-1 on page 7-1). The incubator is ready for use now.

Switching the Incubator Off / Powering Down



1. Keep the **On/Off** button depressed for two seconds.



The display window goes out, except for the readiness indicator icon (rightmost icon in the status display area at D4 in figure 7-1 on page 7-1) and a residual heat temperature readout in case the work space temperature is still higher than 50 °C (122 °F). The incubator is switched off now.

2. If required, unplug the AC power plug to power down the incubator completely.

7-6 Heratherm Thermo Scientific



Temperature Set Value

Heratherm incubators allow for setting the desired work space temperature directly using only a few button presses. After confirming the new temperature set value in the settings pane **Set** (item D5 in figure 7-1 on page 7-1), the user may trace the resulting temperature change in the temperature display field (item D1 in figure 7-1 on page 7-1).

Table 7-4 Adjusting the Temperature Set Value

	Press to activate the menu bar, then use to select the Temperature icon and press to confirm.
Set 888.818	In the flashing settings pane Set, press or to adjust a new temperature set value, then press to confirm your settings.
	The display returns to its default mode. The actual temperature measured in the work space and shown in the temperature display area starts to change until it reaches the newly adjusted set value.



Timer

The **Timer** feature from the menu bar enables the user to turn the incubator on and off at scheduled times. The timer supports three different modes of operation, depending on the user's preferences:

- Countdown-type on or off timer: Turns the incubator on or off after a user-specified period of time. Instructions on setting the preferences for this option are given in table 7-5, while its use as an off timer and on timer are described in table 7-6 on page 7-9 and table 7-7 on page 7-9, respectively.
- Fixed-time on or off timer: Turns the incubator on or off at a scheduled time. Instructions on setting the preferences for this option are given in table 7-8 on page 7-10, while its use as an off timer and on timer are described in table 7-9 on page 7-10 and table 7-10 on page 7-10, respectively.
- Weekly timer: Turns the incubator on or off at scheduled times on specific days of the week. The process for setting the preferences for this option is described in table 7-11 on page 7-11, while instructions for programming the daily turn-on and turn-off times are given in table 7-12 on page 7-11.

Programming a turn-on time causes the incubator to switch off until it is scheduled to restart, while a turn-off time keeps the device running before it shuts down at the user-specified time. The timer starts running immediately as soon as the user confirms his or her entries.

The decontamination routine (see "Decontamination" on page 7-16) overrides timer schedules, that is, a pre-programmed timer will not start before decontamination is finished.

Table 7-5 Presetting the Countdown Timer Mode of Operation

۶	Press to activate the menu bar, then use to select the Settings icon and press to confirm.
	Press to switch to the Timer menu item and confirm the selection with
2010-04-12 10:14 _{PM} ♦ Count Down	Press or to select the Count Down timer mode of operation, then use to confirm the selection.
	The display returns to its default mode.

7-8 Heratherm Thermo Scientific

Table 7-6 Setting a Countdown-type off Timer

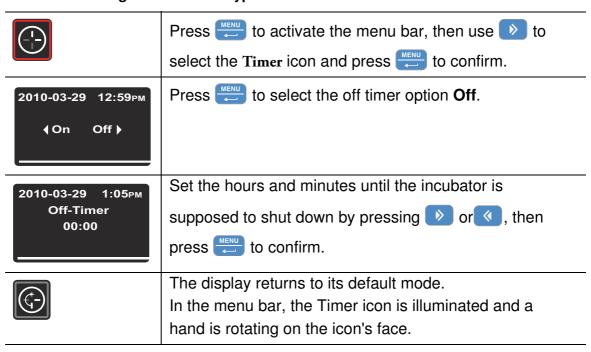


Table 7-7 Setting a Countdown-type On Timer

	Press to activate the menu bar, then use to select the Timer icon and press to confirm.
2010-03-29 12:59 _{РМ}	Press to select the on timer option On , then to confirm.
2010-03-29 1:05рм Оп-Timer 00:00	Set the hours and minutes until the incubator is supposed to turn on by pressing or , then press to confirm.
	The incubator switches off. The display goes out, the Timer icon is illuminated in the menu bar with a rotating hand on its face. Additionally, the readiness indicator icon is illuminated.

Setting a Fixed-time On or Off Timer

Table 7-8 Presetting the "Fixed-time" Timer Mode of Operation

۶	Press to activate the menu bar, then use to select the Settings icon and press to confirm.
	Press to switch to the Timer menu item and confirm the selection with .
2010-04-12 10:14 _{PM} Timer (Absolute)	Press or to select the Absolute fixed-time timer mode of operation, then to confirm the selection.
	The display returns to its default mode.

Table 7-9 Setting a Fixed-time Off Timer


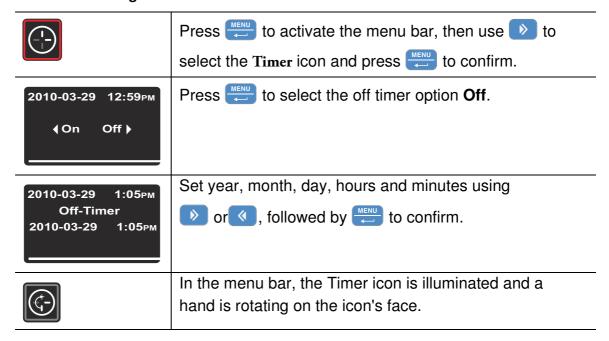


Table 7-10 Setting a Fixed-time On Timer

	Press to activate the menu bar, then use to select the Timer icon and press to confirm.
2010-03-29 12:59 _{РМ}	Press to select the on timer option On , then to confirm.

7-10 Heratherm Thermo Scientific

Table 7-10 Setting a Fixed-time On Timer

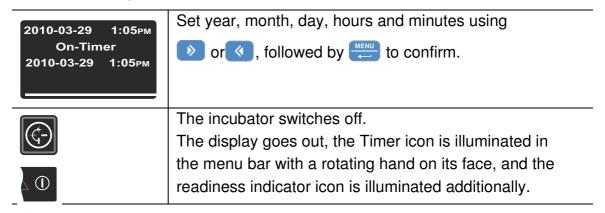

Setting a Weekly Timer


Table 7-11 Presetting the Weekly Timer Mode of Operation

۶	Press to activate the menu bar, then use to select the Settings icon and press to confirm.
	Press to switch to the Timer menu item and confirm the selection with .
2010-04-12 10:14 _{PM} Timer (Weekly)	Press or to select the Weekly timer mode of operation, then to confirm the selection.
	The display returns to its default mode.

Table 7-12 Setting a Weekly Timer

	Press to activate the menu bar, then use to select the Timer icon and press to confirm.
2010-03-29 12:59 _{PM} Timer (On)	In the selection screen shown at left, press to switch from the default On to the Edit option.
2010-03-29 12:59 _{РМ} Timer • Edit •	Select the Edit option by pressing

Table 7-12 Setting a Weekly Timer

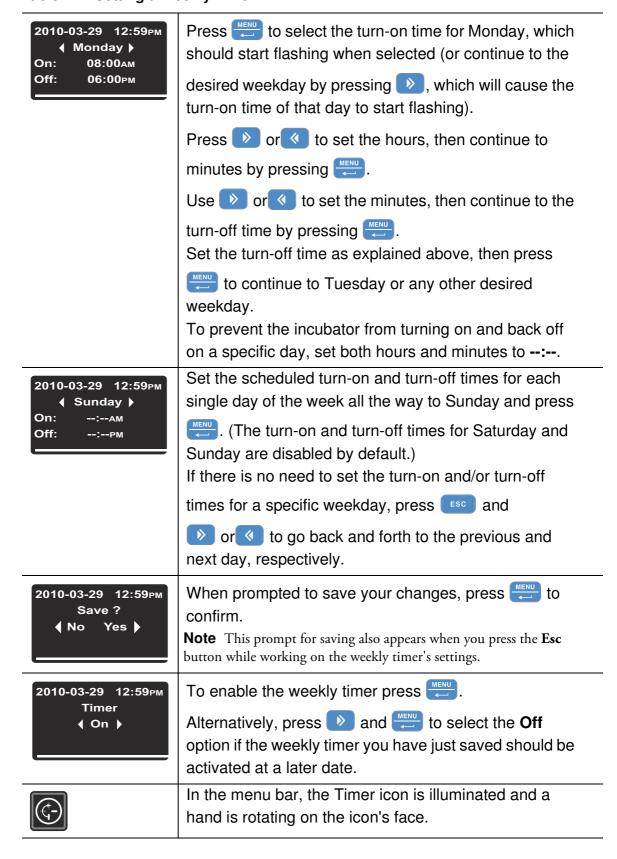

7-12 Heratherm Thermo Scientific

Stopping a Timer


Table 7-13 Stopping an Off Timer Before It Expires

	Press to activate the menu bar, then use to select the Timer icon and press to confirm.
2010-03-29 1:12 _{PM} Stop Timer ? 《 No Yes 》	Press to confirm the Yes default selection.
2010-03-29 4:05 _{РМ} Timer stopped!	The message Timer stopped! appears as a confirmation.
	The display will return to the main menu. In the menu bar, the Timer icon will go out.

Table 7-14 Stopping an On Timer Before it Expires

	To cancel a pre-programmed on timer while the incubator is switched off, hold the On/Off button depressed for a few seconds.
2010-03-29 1:12 _{PM} Stop Timer ? ♦ No Yes ▶	In the Stop Timer? prompt that appears, confirm the default selection Yes by pressing
2010-03-29 4:05 _{РМ} Timer stopped!	The message Timer stopped! appears as a confirmation.
<u>-</u>	The display will return to the main menu. In the menu bar, the Timer icon will go out.

Table 7-15 Stopping a Weekly Timer Before It Expires

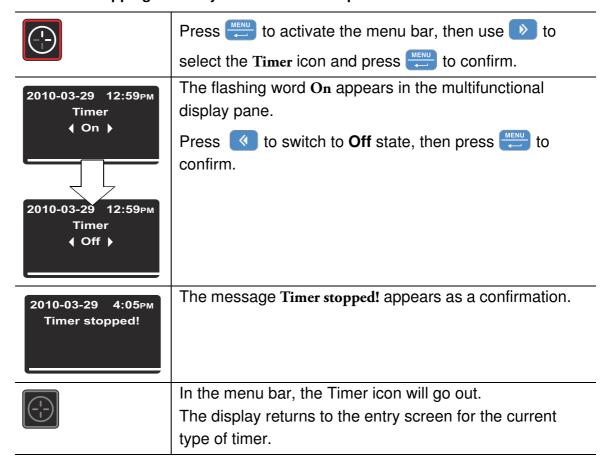

7-14 Heratherm Thermo Scientific



Power Outlet (table-top incubators only)

This menu item (IMH Series only) toggles the built-in AC outlet in the work space of the incubator on and off. While the outlet is in live state the **Power Outlet** icon is illuminated in the menu bar.

Table 7-16 Turning the AC Outlet On

#	Press to activate the menu bar, then use to select the Power Outlet icon and press to confirm.
2010-03-31 1:34 _{PM} 《 On Off 》	In the selection screen that appears, press or to switch to the On option and confirm the selected On option by pressing.
2010-03-31 1:34 _{РМ} Оп	The selected option On remains in the display pane for a few seconds to let you track your action.
\tau	The display returns to its default mode. The AC Outlet icon in the menu bar is illuminated to indicate that the outlet is live.

Table 7-17 Turning the AC outlet off

#	Press to activate the menu bar, then use to select the Power Outlet icon and press to confirm.
2010-03-31 1:34 _{РМ} (On Off)	In the selection screen that appears, press or to switch to the OFF option and confirm the selected Off option by pressing.
2010-03-31 1:35 _{РМ} Оff	The selected option Off remains in the display pane for a few seconds to let you track your action.
#	The display returns to its default mode. The AC Outlet icon in the menu bar is extinguished to indicate that the outlet is switched off.




Decontamination

This menu item (available on IMH-S Series devices only) launches the built-in decontamination routine of the incubator. The decontamination routine consists of a factory-preprogrammed, multi-phase workflow (heating, holding, cooling). It works with a temperature set value of 140 °C (284 °F), which is held over a period of six hours. The overall duration of the process depends on the set temperature that is in place before decontamination is started, on the initial temperature of the work space and miscellaneous other factors. The incubator will be unavailable during that period of time. Be sure to close off the access port with the provided closure plug before starting the decontamination routine.

Table 7-18 Launching Decontamination



Press to activate the menu bar, then use to select the **Decontamination** icon and press to confirm.

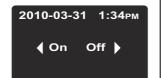

In the selection screen that appears, confirm the selected On option by pressing ...



If a Stop Timer prompt appears while decontamination is launching:

If a timer is running, an additional **Stop Timer** prompt appears at this point (see "Stopping an On Timer Before it Expires" on page 7-13). When you answer that prompt with **Yes**, the timer will be aborted and decontamination be launched instead. Choosing **No** allows the timer to continue. However, if decontamination and a timer overlap, decontamination has priority over the timer, that is, the timer will be put on hold and start later after the decontamination routine has finished.

2010-03-31 1:34_{PM} Decontamination Open door and check interior After pressing to launch decontamination, you will be prompted to open the door and perform a safety check to make sure that all cultures or other temperature-sensitive items have been removed from the work space.

Perform the check as requested, close the door and finally launch decontamination.

7-16 Heratherm Thermo Scientific

Table 7-18 Launching Decontamination



A progress bar and a residual time readout for the decontamination process appear in the multifunctional display pane.



Opening the door while decontamination is running

Opening the front door during the heating and holding phases causes the decontamination routine to stop and wait for the door to be closed. At the same time, the **Door Open** icon (middle item of the group "D4" on page 7-1) is lit in the display window. As soon as the door is closed, decontamination starts over, cycling through the entire process from the start. This may extend the overall duration of the decontamination effort substantially.

Power Outage during a Decontamination Run

A power outage relaunches the decontamination routine automatically. The Alarm icon (leftmost item of the group "D4" on page 7-1) is illuminated in the display window, and a E007 Power Down Error message

appears at the end of the process. Press the button to acknowledge that alarm.



Put up bio-hazard warning signs!

Under specific circumstances, there is a risk that bacteria, viruses, fungi, prions, and other biological substances survive when the decontamination routine is aborted due to a power outage. Normally, the incubator starts over with the decontamination cycle without any need for user intervention upon return of power. Users who fail to notice the power outage, may open the door and/or load the incubator with samples, thus exposing themselves or their samples to bio-hazards without being aware of that risk.



The **Decontamination** icon in the menu bar is illuminated to indicate that the decontamination process is running.

Table 7-18 Launching Decontamination

2010-03-31 4:13PM Decontamination finished!	When decontamination is completed, the message Decontamination finished! appears in the multifunctional display pane. Press the button to acknowledge that message.
	The Decontamination icon in the menu bar will go out. The settings that were in place before the decontamination routine was started - for example, fan speed, will be restored.

Table 7-19 Stopping Decontamination Prematurely

	Press to activate the menu bar, then use to select the Decontamination icon and press to confirm.
2010-03-31 1:34pm Quit Decon? ♦ No Yes ▶	In the selection screen that appears, press to switch to Yes and confirm the selection with
	The display returns to its default mode. The Decontamination icon in the menu bar will go out. The settings that were in place before the decontamination routine was started - for example, fan speed, will be restored.

7-18 Heratherm Thermo Scientific



Fan

This menu item allows for turning on the fan inside the work space and gradually adjusting its speed in one go. The current setting of the fan is indicated by the illuminated **Fan** icon in the menu bar and the five-level bar graph (see item D6 in figure 7-1 on page 7-1) for the fan speed setting located directly above the icon.

Table-top incubators

The fan can be adjusted in 6 levels:

- 0% (all chevrons extinguished)
- 20% (1 chevron illuminated)
- 40% (2 chevrons illuminated)
- 60% (3 chevrons illuminated)
- 80% (4 chevrons illuminated)
- 100% (5 chevrons illuminated)

Table 7-20 Turning on the Fan


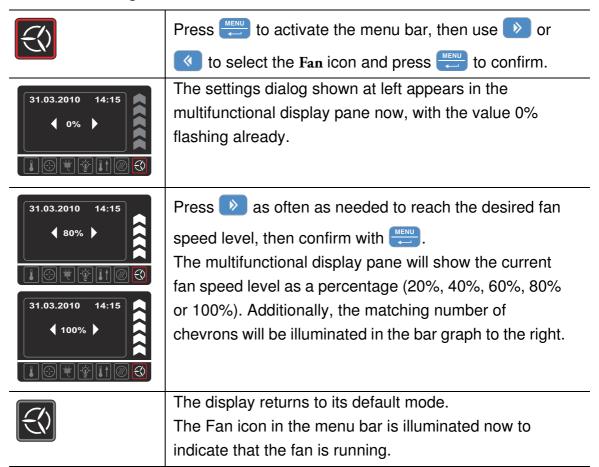


Table 7-21 Adjusting Fan Speed or Turning the Fan Off

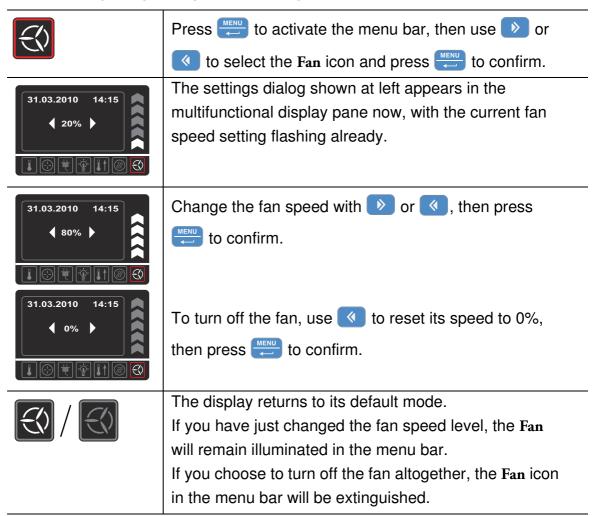


Floor stand incubators

The unit contains a regulated fan; fan speed can be lowered (adjusted) in 2 stages:

Minimum fan speed level (chevrons 1 through 3 illuminated)

Maximum fan speed level (chevrons 1 through 5 illuminated)

Table 7-22 Turning on the Fan

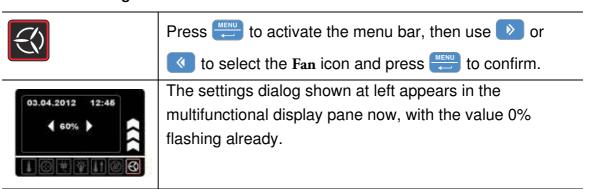

7-20 Heratherm Thermo Scientific

Table 7-22 Turning on the Fan


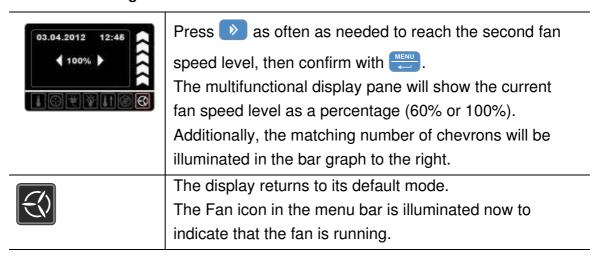


Table 7-23 Adjusting Fan Speed or Turning the Fan Off (table-top incubators only)



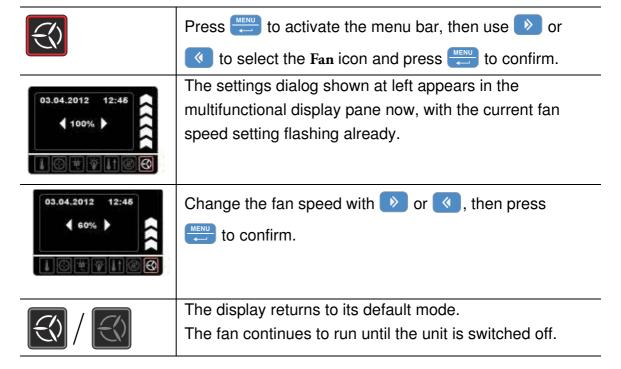

Settings

The **Settings** menu item opens a submenu populated with various commands for viewing general status information on the Heratherm unit and setting for the operation of the incubator or its display window:

- Read access to error log
- Calibrating the incubator
- Setting date and time
- Toggling the temperature display unit between °C and °F
- Entering a configuration control code

Instructions for using these features are given in the following. Also supported is an option for presetting the timer mode of operation, as explained previously in the section "Timer" on page 7-8.

Error Log

Users calling customer service for support may be asked by the Thermo Fisher Scientific agent to supply information from the error log of the incubator. This internal memory may be accessed by choosing the **Settings** -> **Error** menu item. It enables the user to browse through the most recent 22 alarm messages that were caused by hardware or control loop errors. Each error is displayed with the date and time of its occurrence, a brief clear text description and an internal error code.

Error codes and instructions for clearing alarm conditions appear in the section "Error Codes" on page 12-1.

Table 7-24 Reading the Error Log

F	Press to activate the menu bar, then use or to select the Settings icon and press to confirm.
2010-04-06 1:33рм (Error)	Press to select the Error item from the Settings submenu.
2010-04-06 1:36рм Error 0 2010-04-06 1:31рм Fan Error (E009)	The first entry of the error log is displayed, named "Error 0."

7-22 Heratherm Thermo Scientific

Table 7-24 Reading the Error Log

2010-04-06 1:37рм Error 1 2010-04-06 1:34рм Heat Relay (Е109)	Press to go to the next entry (or to go back to the previous one). After the entry numbered 21 the display wraps and returns to the beginning of the error log, displaying Error 0 again.
F	To exit from the error log and return to normal display mode press twice. The Settings icon in the menu bar will go out.

Calibration

The **Settings** -> **Calibration** menu item enables the user to initiate a temperature calibration process (see "Temperature Calibration Procedure" on page 10-4) for the built-in temperature sensors and choose whether calibration should be accomplished manually or automatically:

• The **Manual** option allows for entering an absolute temperature directly, as measured - for example, using an external reference sensor.

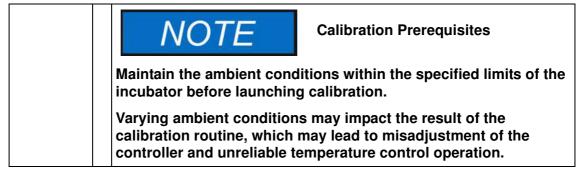


Table 7-25 Entering the Calibration Reference Temperature Manually

	Prepare for temperature calibration (see "Preparing Temperature Calibration" and "Comparison Measurement Procedure" on page 10-3).
F	Press to activate the menu bar, then use or to select the Settings icon and press to confirm.
2010-04-06 1:33рм 4 Error)	The Error menu item from the Settings submenu appears in the multifunctional display pane.

Thermo Scientific Heratherm 7-23

Table 7-25 Entering the Calibration Reference Temperature Manually


2010-04-06 1:33 _{РМ}	Press to switch to the Calibration menu item and confirm the selection with .
2010-04-06 1:33 _{PM} Calibration ◀ Manual ▶	In the Calibration selection screen, press to choose the preselected option Manual.
2010-04-06 1:33 _{РМ} Calibration 36.9°С	In the settings dialog that appears, set the temperature measured with the external reference sensor by using or and confirm the settings with.
	The newly entered value will be stored and used to calibrate the internal temperature sensors with the value measured by the reference sensor. The display returns to its default mode. The Settings icon in the menu bar will go out.

Date and Time

The **Settings -> Time / Date** option allows for switching between the international time and date display formats and for setting the time and date of the internal clock. There are two display formats to choose from:

- European date format *DD.MM.YYYY* and 24-hours time format. Example: 07.04.2010 and 15:05.
- US standard date format *YYYY-MM-DD* and 12-hour time format with *AM/PM* suffix. Example: 2010-04-07 and 3:05 PM.

Table 7-26 Setting the Date Format

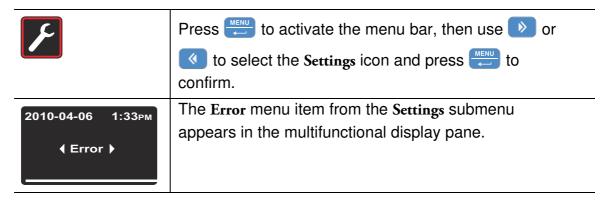

7-24 Heratherm Thermo Scientific

Table 7-26 Setting the Date Format


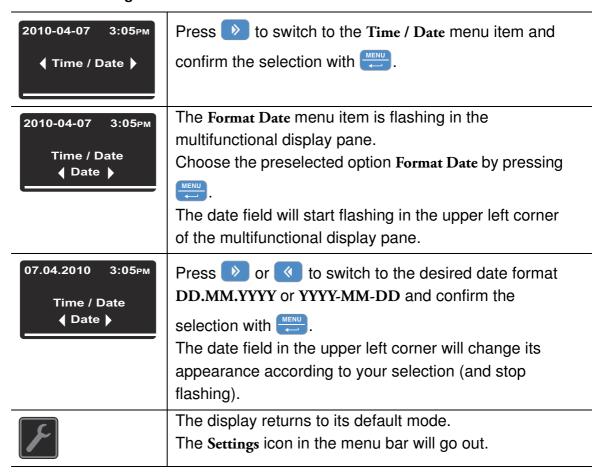


Table 7-27 Setting the Time Format

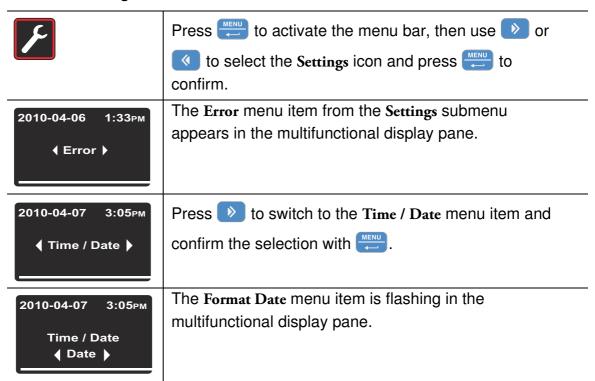

Thermo Scientific Heratherm 7-25

Table 7-27 Setting the Time Format



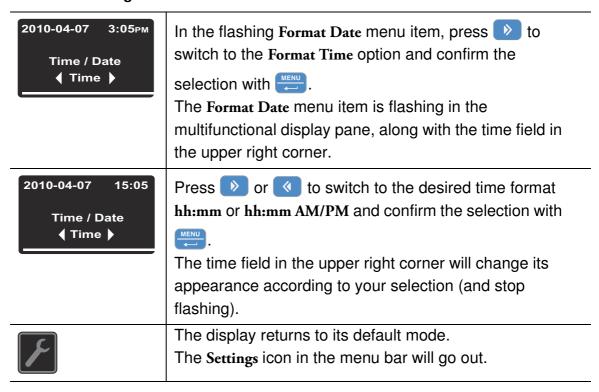


Table 7-28 Setting Date and Time

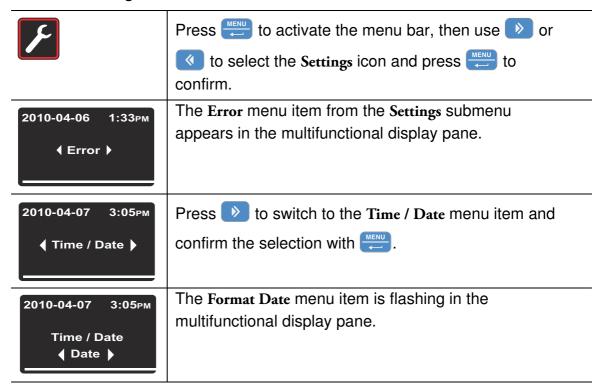

7-26 Heratherm Thermo Scientific


Table 7-28 Setting Date and Time

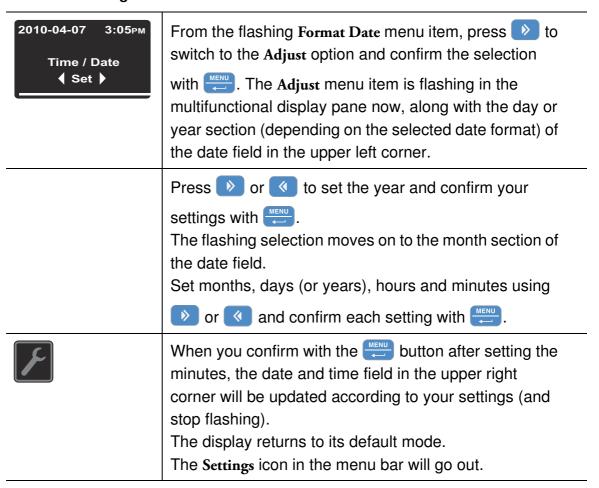

Temperature Display Unit

The **Settings** -> °C / °F menu item allows for toggling the incubator used for displaying temperatures between degrees Centigrade and Fahrenheit.

Table 7-29 Toggling the Temperature Display Unit

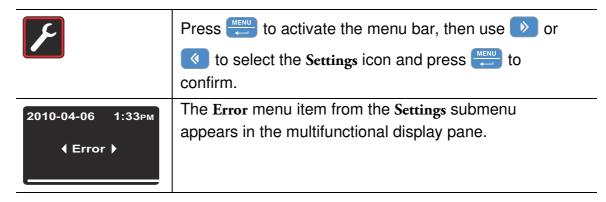


Thermo Scientific Heratherm 7-27

Table 7-29 Toggling the Temperature Display Unit



Press to switch to the °C / °F menu item and confirm the selection with ...

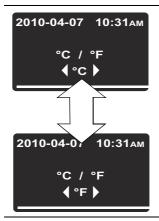

The multifunctional display pane changes to the following selection screen, with the currently unused temperature display unit (factory setting is $^{\circ}$ C) flashing.

Confirm the selection with



The temperatures in the display field (item D1 in figure 7-1 on page 7-1) and settings pane Set (item D5 in the same figure) will be displayed with the newly selected unit.

The display returns to its default mode.

The Settings icon in the menu bar will go out.

7-28 Heratherm Thermo Scientific

Shut-down

Shutting the Incubator Down

This chapter provides instructions for shutting the incubator down for prolonged periods of time, that is, at least for several days in a row.





Contamination hazard

If the work space surfaces are contaminated, harmful biological material may spread to the environment of the incubator.

To rule out any risk for subsequent users, perform a full cleaning, disinfection and decontamination cycle to the standards set out in the section "Cleaning and Disinfection" on page 9-1 if you suspect (or if you are positive) that harmful biological material has been processed with the incubator.

- 1. Remove the containers with the cultures, all accessories, and other objects from the work space.
- 2. Clean and disinfect the work space, as explained in the section "Cleaning and Disinfection" on page 9-1 or if the above warning note applies run the dry decontamination routine at 140 °C / 284 °F (see "Decontamination" on page 7-16).
- 3. When cleaning and disinfection and/or decontamination are done, turn the incubator off using the control panel.
- 4. Unplug the power cord and secure it against accidental reconnection.
- 5. Until the incubator is shut down, the work space must be continuously ventilated. Leave the glass door and the outer door open and secure them against accidental closure.

Thermo Scientific Heratherm 8-1

Cleaning and Disinfection

Cleaning





Incompatible cleaners

Some incubator components are made of plastics. Solvents may attack plastics. Strong acids or alkaline solutions may cause embrittlement of plastics.



Moisture-sensitive components

Do not spray cleaning agent onto the control panel and the interfaces at the rear of the incubator. When wiping the incubator clean, always make sure that no moisture enters into these components.

Wipe the display window down with a slightly dampened cloth, then wipe dry with a rag made of 100% microfiber.

Cleaning Exterior Surfaces

Remove dirt residues and depositions thoroughly using a solution of lukewarm water and commercial detergent.

Wipe the surfaces clean using a clean cloth and clear water.

Then, wipe the surfaces dry using a clean cloth.

Wipe / Spray Disinfection

The manual wipe and spray disinfection is a three-stage process:

- predisinfection,
- cleaning,
- final disinfection.

Thermo Scientific Heratherm 9-1





Alcoholic disinfectants!

Disinfectants having an alcohol content of more than 10% may form, in combination with air, easily combustible and explosive gas mixtures.

When using such disinfectants, avoid open flames or exposure to excessive heat during the entire disinfection process!

Use such disinfectants only in adequately ventilated rooms.

After the disinfectant has been allowed to react, wipe the cleaned incubator components thoroughly dry.

Observe safety regulations to avoid fire and/or explosion hazard caused by alcohol-containing disinfectants.





Chloride-containing disinfectants!

Chloride-containing disinfectants can corrode stainless steel.

Use only disinfectants that do not affect stainless steel!

Preparing the Manual Wipe/Spray Disinfection





Health hazard

The surfaces of the work space may be contaminated. Contact with contaminated cleaning liquids may cause infections. Disinfectants may contain harmful substances.



When cleaning and disinfecting, always observe the safety instructions and hygiene guidelines!

- · Wear safety gloves.
- · Wear safety goggles.
- Wear mouth and respiratory system protection gear to protect your mucous membranes.
- Observe the safety instructions of the disinfectant's manufacturer and the hygiene supervisor.

9-2 Heratherm Thermo Scientific

Predisinfection

- 1. Remove all samples from the work space and store them in a safe place.
- 2. Spray disinfectant onto the surfaces of the work space and of the accessories or wipe the surfaces clean using disinfectant.



Risk of injury caused by breaking of glass panel

The glass panel may only be removed by properly trained and authorized specialists.

The glass panel must be held by two persons.

3. Allow time for the disinfectant to act as specified by the manufacturer.



Disinfecting hard-to-reach components

Spray the sensor and other hard-to-reach components with disinfectant!

Thermo Scientific Heratherm 9-3

Cleaning

- 1. Remove all internals from the specimen chamber.
- 2. Wipe off the specimen chamber surfaces and the internals that have been removed from the chamber using lukewarm water mixed with standard rinsing agents. Completely remove any tenacious impurities using rinsing agent and warm water.
- 3. Re-rinse the cleansed surfaces 3 to 5 times with autoclaved water to completely remove any cleaning agent residues.
- 4. After this, wipe the cleansed surfaces and internals dry with a soft, sterile cloth.

Final Disinfection

- 1. Spray the specimen chamber surfaces and the internals removed from this chamber again with disinfectant, or wipe them down.
- 2. Let the disinfectant work on the surfaces/internals as detailed in the manufacturer's instructions.
- 3. Re-install the internals in the specimen chamber.

Starting the Decontamination Routine

For detailed instructions, please refer to "Decontamination" on page 7-16 (only applicable for IMH-S Series). Be sure to close off the access port with the provided closure plug before starting the decontamination routine.



Recommendations for decontamination

The user should be familiar with the internationally recognized "Laboratory Biosafety Manual" (from the World Health Organization (WHO)) and with national, pertinent recommendations. These manuals/recommendations provide information about decontaminants, their use, dilution, properties and possible applications.



Opening the door during the decontamination routine

If the door will be opened and closed again whilst decontamination is running, the routine returns to a phase which will ensure a continuation without faults.

9-4 Heratherm Thermo Scientific





Bio-hazard

Be sure to determine the current operating condition of the incubator before you open the door!

Under specific circumstances, there is a risk that bacteria, viruses, fungi, prions, and other biological substances survive when the decontamination routine is aborted due to a power outage. Normally, the incubator starts over with the decontamination cycle without any need for user intervention upon return of power. Users who fail to notice the power outage, may open the door and/or load the incubator with samples, thus exposing themselves or their samples to bio-hazards without being aware of that risk.

Decontamination must be performed immediately if any biohazardous material is spilled in or on the incubator.

Thermo Scientific Heratherm 9-5

Maintenance

Maintenance and inspection at regular intervals of the features and components listed below are mission-critical to maintain the product in a fully operative and safe condition and avoid malfunctions due to ageing and wear. Failure to perform maintenance on a regular basis may result in:

- deviations in heating performance
- damage to samples
- loss of control over temperature distribution throughout the work space

Inspections and Checks

To ensure the operational performance and safety of the incubator and its functions, the components listed below must be checked at regular intervals.

Regular Checks

- Check the incubator for overall cleanliness and remove any residues from previous processes.
- Check the locking screw for the glass panel
- To avoid incubator operation without an appropriate fresh air supply, check the air filter (optional) in the air inlet path for contamination.

Semi-annual Inspection

- Check integrity and proper seating of the seal.
- Swap the fresh-air filter (accessory) in air inlet.
- Perform functional check of the control panel and of the incubator's built-in controller.
- Perform electrical safety check in accordance with the relevant national regulations.

Thermo Scientific Heratherm 10-1



Functional Check

If safety devices were removed or disabled for inspections, the incubator must not be operated before the safety devices have been reinstalled and checked for their correct function.



Spare Parts and User Modifications

To avoid major malfunctions of the incubator and associated safety hazards that may result in death, serious injuries, or damage to the incubator and other equipment, use spare parts approved by Thermo Scientific only. Third-party spares without Thermo Scientific approval void the limited warranty.

Do not modify the incubator in any way without obtaining the prior written authorization from Thermo Scientific. Unauthorized modifications may compromise operational safety and give rise to hazards that may result in death, serious injuries, or damage to the incubator and other equipment.

Service Intervals

During running operation, the following service works must be performed:

3-monthly service

• Perform the comparative temperature measurement outlined in the following section.

Annual Service

• Have the incubator inspected and services by a Thermo Scientific authorized Technical Service agent.



Service contract

Thermo Electron LED GmbH offer a product-specific service contract comprising all necessary tests and service work.

10-2 Heratherm Thermo Scientific

Preparing Temperature Calibration

To determine the exact measured value of the incubator's integral temperature sensor, a temperature comparison measurement must be performed every three months. If a major temperature deviation is found during this check, temperature calibration is required. During this process, the temperature controller of the incubator is set to the value measured during the temperature comparison measurement.

Use a calibrated measuring instrument with an accuracy of < ± 0.1 °C (0.18 °F) for this test.

To minimize temperature variations during the measurement, put the measuring sensor in an isothermal container (such as a bowl filled with glycerol) before placing it in the work space. Use the center of the work space as the reference location for the comparison measurement.

NOTE

Isothermal container

Do not use a container filled with water as an isothermal container because the evaporation of water will result in a lower temperature reading.



Excessive work space temperatures

Excessive work space temperatures after the calibration may be reduced by leaving the doors open for approx. 30 seconds.


Comparison Measurement Procedure

- 1. Turn the incubator on using the power switch.
- 2. Set the temperature set value and allow the incubator to stabilize. This may take several hours.
- 3. Place the measuring device in the center area of the work space. Alternatively, a temperature sensor may be positioned in this location. Route the connecting cable between the glass panel and the interior tank.
- 4. Close the doors.
- 5. Wait until the temperature value displayed on the measuring instrument has stabilized.
- 6. Use the temperature reading from the measuring device to calibrate temperature control manually, as explained "Entering the Calibration Reference Temperature Manually" on page 7-23.

Thermo Scientific Heratherm 10-3

Temperature Calibration Procedure

For detailed instructions on how to perform a manual temperature calibration, please refer to the instructions in the section "Calibration" on page 7-23.

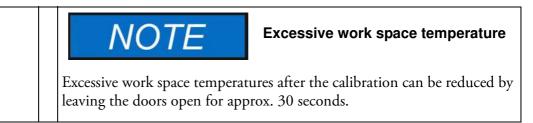

Replacing the Door Seal

The door seal of the outer door is located in the retaining slot.

The door seal should be inspected for any signs of embrittlement at half-yearly intervals.

No tools are required to replace the seal.

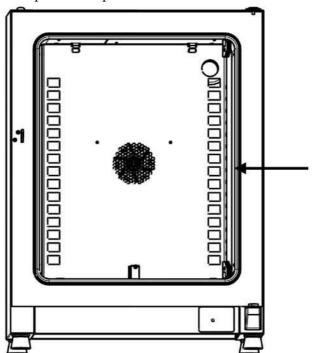


Figure 10-1 Door Seal Replacement

- 1. Pull the seal out of the guide slot.
- 2. Starting on the hinge side of the door, position the end of the new seal at the location indicated by the arrow in figure 10-1 above.
- 3. Gently press the seal into the slot, working around the circumference of the door. In corner areas in particular ensure that the seal lip is installed without any wrinkles and that the seal is not stretched or compressed.
- 4. Make sure that the seal is positioned correctly in the slot and flush with the door frame. Correct, if necessary.

10-4 Heratherm Thermo Scientific

Replacing the Power Cord

If the power cord is damaged, it must be replaced by an original spare part. Using a standard power cord with a lower thermal withstand capabilities is prohibited.

Returns for Repair

Prior to returning any materials, please contact our Customer Service Department for a "Return Materials Authorization" number (RMA).

Material returned without an RMA number will be refused.





Contamination hazard

The incubator may have been used for treating and processing infectious substances, which may have caused contamination of the incubator and its components.

Prior to return shipment, it is therefore mandatory that all incubator components be properly decontaminated..

- Clean the incubator components thoroughly, then disinfect or decontaminate them (depending on application).
- Fill in and attach a safety declaration with details on decontamination activities performed to the items that are to be repaired.

Thermo Scientific Heratherm 10-5

Disposal





Contamination hazard

The incubator may have been used for treating and processing infectious substances, which may have caused contamination of the incubator and its components.

Prior to disposal, it is therefore mandatory that all incubator components be properly decontaminated.

Clean the incubator components thoroughly, then disinfect or decontaminate them (depending on application).

Attach a declaration of decontamination with details on decontamination activities performed to the items that are to be disposed of.

Overview of Materials Used

Component	Material
Thermal insulation components	Glass wool
Printed circuit boards	Coated electrical components contain various plastics materials. Components mounted on circuit boards containing epoxy resin bonder.
Plastic components, general	see material labelling
Exterior housing	Galvanized steel sheet, painted
Device rear panel	Galvanized steel sheet
Outer door	Galvanized steel sheet, painted, + stain- less steel (optional)
Door inner panel	Stainless steel 1.4301
Control panel and display window protective foil	Polyethylene
Heater	Silicone-sheathed resistance heater wires, stainless-steel-jacketed resistance heater wires with the IMH-S model

Thermo Scientific Heratherm 11-1

Component	Material
Work space containers, installed components and shelves	Stainless steel 1.4301
Door frame seal	Silicone
Glass screen	Sodium silicate glass
Fan wheel	Stainless steel 1.4016
Cables	Plastic-sheathed stranded copper wire
Packaging	Corrugated board, polyethylene film, and styrofoam, chemically untreated wood

11-2 Heratherm Thermo Scientific

Error Codes

Table 12-1 below lists the error messages that may appear in the control panel display window (see "Error Log" on page 7-22) and provides instructions for clearing such alarms.

Table 12-1 Heratherm Incubator Error Codes

Error Message & Code	Root Cause	Alarm Response	Alarm Clearing Instructions*		
Door Open Error (E001)	The door switch (IMH-S Series only) has triggered an alarm because the door has been open for more than 10 minutes.	Audible alarm activated, alarm relay energized, message shown on display.	Close the door.		
Display Error (E002)	Display communication error. The built-in controller was unable to restore communication with the control panel.	error. The built-in control- er was unable to restore alarm relay energized, message shown on dis- communication with the play. Reset after 30 s.			
Mirrored Parameter Loaded (E003)			Check the latest settings, for example the set value.		
Factory Parameter Loaded (E004)	The controller was unable to read the mirrored parameter set and had to resort to factory-preset parameters.	Fallback to factory-preset parameters. Audible alarm activated, alarm relay energized, message shown on display. User-specific settings may be lost - for example, the temperature display unit preference, or user programs.	Acknowledge by pressing Esc . Re-enter cus- tomer-specific settings.		
Default Parameter Loaded (E005)	The controller was unable to read the factory-preset parameters.and had to resort to default settings	Fallback to default parameters. Audible alarm activated, alarm relay energized, message shown on display. The device is completely inoperative.	Call service.		
Disinfection Routine Error (E006)	Process error in decontamination routine. Controller, heater, or fan error, etc.	Abort decontamination. Audible alarm activated, alarm relay energized, message shown on dis- play.	Acknowledge by pressing Esc. Relaunch decontamination. If this doesn't solve the problem, call service.		

Thermo Scientific Heratherm 12-1

Table 12-1 Heratherm Incubator Error Codes

Error Message & Code	Root Cause	Alarm Response	Alarm Clearing Instructions*	
Power Down Error (E007)	Power has been cut off (power outage) while the device was running.	Audible alarm activated, alarm relay energized, message shown on display.	Check the power supply. Power up then device, then acknowledge the alarm by pressing	
Fan Error (E009)	Fan speed out of range.	Audible alarm activated, alarm relay energized, message shown on display. Heater disabled by controller.	Acknowledge by pressing ESC. If this doesn't solve the problem, call service.	
Config Error (E012)	General device configuration error.	Audible alarm activated, alarm relay energized, message shown on display. The device is completely inoperative.	Call service.	
OTP error (E013)	Klixon contact not closed.	Overtemperature Protection fault. Audible alarm activated, alarm relay energized, message shown on display. Bridging across Klixon contact has failed.	Restart the device. If this doesn't solve the problem, call service.	
Incorrect voltage (E014)	The applied voltage is too high or too low.	Audible alarm activated, error message shown on display.	Apply the correct voltage as indicated on the name-plate and acknowledge the error.	
Sensor Error (E100)	Process sensor damaged. The actual measured value is out of range.	Audible alarm activated, alarm relay energized, message shown on display. Control transferred to reference sensor. If both sensors are defective, disable all control circuits.	Call service.	
Temperature Too High (E101)	Actual measured value exceeds permissible range. The Triac is defective.	Process protection activated, control continues on set value. Audible alarm activated, alarm relay energized, E101 message shown on display.	If the error occurs repeatedly, call service.	
Temperature Too Low (E102)	Actual measured value falls short of permissible range . Only on devices equipped with a door switch.	Audible alarm activated, alarm relay energized, message shown on display. Temperature control continued.	Check the AC mains supply for an undervoltage condition and have the problem remedied, if necessary. Check whether the incubator has been overloaded with samples. If this is not the root cause of the problem, call service.	

12-2 Heratherm Thermo Scientific

Table 12-1 Heratherm Incubator Error Codes

Error Message & Code	Root Cause	Alarm Response	Alarm Clearing Instructions*
Temperature Not Plausible (E103)	The difference between the control and reference sensors exceeds the maximum permissible deviation, rendering the measurement implausible.	Device uses the sensor that indicates the higher temperature for servo control. Audible alarm activated, alarm relay energized, message shown on display. Error can be acknowledged and doesn't reset.	If the error occurs repeatedly, call service.
Calibration Value Too High (E104)	The calibration reference value calculated on the basis of the user input falls short of the upper limit for calibration references.	Fallback to previous calibration reference. Audible alarm activated, alarm relay energized, message shown on display.	Check the external reference sensor for proper function and replace, if necessary. If this does not solve the problem, call service.
Calibration Value Too low (E105)	The calibration reference value calculated on the basis of the user input exceeds the lower limit for calibration references.	Fall back to previous calibration reference. Audible alarm activated, alarm relay energized, message shown on display.	Check the external reference sensor for proper function and replace, if necessary. If this does not solve the problem, call service.
Constant Sensor Sig- nal (E106)	None of the decimal places of the A/D-converter output for the process sensor has changed over a specific time period.	ne A/D-converter out- for the process sensor changed over a spe- ence sensor, audible alarm activated, alarm relay ener- gized, message shown on	
Constant Reference Sensor Signal (E107)	None of the decimal places of the A/D-converter output for the reference sensor has changed over a specific time period.	he A/D-converter out- for the reference sen- has changed over a cess sensor, audible alarm activated, alarm relay ener- gized, message shown on	
Constant Sample Sensor Signal (E108)	None of the decimal places of the A/D-converter output for the sample sensor has changed over a specific time period.	Audible alarm activated, alarm relay energized, message shown on display.	Call service.
Heating Relay Error (E109)	The voltage measurement has indicated a defect in the heater circuit relay.	Device completely inoperative, audible alarm activated, alarm relay energized, message shown on display.	Be sure to pull the power plug and disconnect the device from the AC mains. Call service.

Thermo Scientific Heratherm 12-3

Table 12-1 Heratherm Incubator Error Codes

Error Message & Code	Root Cause	Alarm Response	Alarm Clearing Instructions*
Heating Triac Error (E110)	The voltage measurement has indicated a defect in the triac	Audible alarm activated, alarm relay energized, message shown on display. Overheat protection activated to prevent destruction of the samples. Audible alarm returns upon acknowledgement.	Call service and switch the device off.
Temperature Too High (E111)	The actual measured value exceeds the upper limit of the permissible error range.	e actual measured value Audible alarm activated, alarm relay energized,	
Sensor Error (E112)	The measured actual value is out of range.	Audible alarm activated, alarm relay energized, message shown on display. Control continues on process sensor. If both sensors are defective, all control circuits will be disabled.	If error does not go away, call service.
Sensor Error (E113)	Sample sensor damaged. The actual measured value is out of range.	Audible alarm activated, alarm relay energized, message shown on display.	Call service.
ADC Error (E114)	A/D converter does not supply a plausible output. Measurement across reference resistor has failed. ADC converter may be defective.	Audible alarm activated, alarm relay energized, message shown on display. All control circuits disabled.	Call Service.
Watchdog error (E115)	Watchdog test failed on power-up.	Audible alarm activated, alarm relay energized, message shown on display.	Call service.

^{*}Clearing should mute the audible alarm, de-energize the alarm relay, and clear the message from the control panel display.

12-4 Heratherm Thermo Scientific

Technical Data

The technical data are valid only for an empty device equipped with three shelves, a spray-painted outer enclosure and a power line voltage 120 V/60 Hz. Options may have an impact on the specified performance.

Table 13-1 Technical Data - Series IMH 60/100/180 and IMH 60-S/100-S/180-S

Parameter	Unit	IMH 60	IMH 100	IMH 180	IMH 60-S	IMH 100-S	IMH 180-S
Process (first value is with fan off, second with fan running.)							
Temperature deviation from set value at 37 °C (99 °F), spatial.	12	10.0/10.0	10.0/10.0	10.0/10.4	10.0/10.0	10.0/10.0	10.0/10.4
Typical value	K	±0.6/±0.2	±0.6/±0.3	±0.6/±0.4	±0.6/±0.2	±0.6/±0.3	±0.6/±0.4
Max. value	K	±0.6/±0.3	±0.6/±0.4	±0.6/±0.5	±0.6/±0.3	±0.6/±0.4	±0.6/±0.5
Temperature deviation from set value at 37° C (99 °F), over time.	K	±0.1	±0.1	±0.1	±0.1	±0.1	±0.1
Work space atmosphere, Min.	°C/°F	plus 5 ℃/41°F	plus 5 ℃/41℉	plus 5 ℃/41 ℉	Amb. temp. plus 5 °C/41 °F	Amb. temp. plus 5 °C/41 °F	Amb. temp. plus 5 °C/41 °F
Max.	°C/°F	105 ℃/221 ℉	105 ℃/221 °F	105 ℃/221 °F	105 ℃/221 F	105 ℃/221 °F	105 ℃/221 ℉
Heat-up time (work space unoccupied, from 25 °C (77 °F) to 98% of set temperature of 37 °C/98.6 °F). Max. value/typical value	min	25 / 22	35 / 42	35 / 40	25 / 22	35 / 42	35 / 42
Recovery time (work space unoccupied, door open for 30 s, to set temperature of 37 °C/98.6 °F). Max. value/typical value.	min	6/4	5/4	5/4	5 / 2.5	4 / 2.5	4 / 2.5
Heat dissipation to environment (at set temperature of 37 °C/98.6 °F and room temperature of 25 °C/77 °F)	W	23/65 ±10%	30/68 ±10%	36/78 ±10%	23/65 ±10%	30/68 ±10%	36/78 ±10%
Overall dimensions							
Height	mm/in	720/ 28.3	820/ 32.3	920/ 36.2	720/ 28.3	820/ 32.3	920/ 36.2
Width	mm/in	530/ 20.8	640/ 25.2	640/ 25.2	530/ 20.8	640/ 25.2	640/ 25.2

Thermo Scientific Heratherm 13-1

Table 13-1 Technical Data - Series IMH 60/100/180 and IMH 60-S/100-S/180-S

Parameter	Unit	IMH 60	IMH 100	IMH 180	IMH 60-S	IMH 100-S	IMH 180-S	
Depth	mm/in	565/ 25.2	565/ 25.2	738/ 29.1	565/ 25.2	565/ 25.2	738/ 29.1	
Overall weight	kg/lbs	45/99	56/123	70/154	45/99	56/123	70/154	
Loading capacity per shelf	kg/lbs		25/55			25/55		
Max. overall loading capacity per device	kg/lbs	50/110	50/110	75/165	50/110	50/110	75/165	
Electrical data								
Power consumption	W	600	840	1020	1390	1390	1390	
Maximum current	Α	5.0	7.0	8.5	11.6	11.6	11.6	
Earthing system (e. g. 1/N/PE)		1/N/PE	1/N/PE	1/N/PE	1/N/PE	1/N/PE	1/N/PE	
Power line frequency	Hz		60			60		
Power line voltage +/- 10 %	V		120			120		
IP protection system			IP 20			IP 20		
Protection class			I		I			
Overvoltage category			II		II			
Device fusing, building side	Α		16			16		
Device fusing, on PCB	Α	2 x 16			2 x 16			
Environmental conditions	S							
Min. ambient temperature	°C/°F	18/65			18/65			
Max. ambient temperature	°C/°F	32/90			32/90			
Max. humidity in service, non condensing	% r.F./ % r.H.	80, non condensing			80, non condensing			
Min. storage temperature	°C/°F		20/68 20/68					
Max. storage temperature	°C/°F		60/140			60/140		
Max. humidity in storage, non condensing	% r.F./ % r.H.	90, r	non conden	sing	90,	non conder	sing	
Post-transport acclimation time	h		2			2		
Noise level	dB(A)		45		45			
Pollution degree			2			2		
Site conditions								
Maximum altitude above sea level	m/y NN		2000/2187			2000/2187		
Minimum side clearance	mm/in	50/2				50/2		
Minimum front clearance	mm/in	590 / 23.2	590 / 23.2 690 / 27.2 814 / 32			690 / 27.2	814 / 32	
Minimum back wall clearance	mm/in		80/3.2			80/3.2		

13-2 Heratherm Thermo Scientific

Table 13-1 Technical Data - Series IMH 60/100/180 and IMH 60-S/100-S/180-S

Parameter	Unit	IMH 60	IMH 100	IMH 180	IMH 60-S	IMH 100-S	IMH 180-S
Minimum bottom clearance	mm/in		200/8			200/8	
Minimum top clearance	mm/in		300/12			300/12	

Table 13-2 Technical Data - Series IMH 400-S / IMH 750-S

Parameter	Unit	IMH 400-S	IMH 750-S
Process			
Work space atmosphere, Min. Max.	°C/°F	Amb. temp. plus 5 °C/41°F 105 °C/221 °F	Amb. temp. plus 5 °C/41°F 105 °C/221 °F
Temperature deviation from set value at 37 °C (99 °F), spatial. Max. value/typical value	K	±0.3/±0.2	±0.4/±0.3
Temperature deviation from set value at 37 °C (99 °F), over time	К	±0.2	±0.2
Heat-up time (work space unoccupied, from 25 °C (77 °F) to 98% of set temperature of 37 °C/99 °F))	min	35	30
Recovery time (work space unoccupied, door open for 30 s. Max. value/Typical value	min	<3/<2	<3/<2
Heat dissipation to environment (at set temperature of 37 °C (99 °F) and room temperature of 25 °C/77 °F)).	W	87 ±10%	149 ±10%
Overall dimensions			
Height (with casters)	mm/in	1655/ 65.2	1655/ 65.2
Width	mm/in	755/ 29.7	1215/ 47.8
Depth	mm/in	770/ 30.3	770/ 30.3
Overall weight	kg/lbs	145/320	205/452
Loading capacity	kg/lbs	30/66	30/88
Max. overall loading capacity per device	kg/lbs	75/165	150/330
Electrical data			
Power rating	W	1380	1560
Max. Current	Α	11.5	13.0
Earthing system (e. g. 1/N/PE)		1/N/PE	1/N/PE
Power line frequency	ower line frequency Hz 60		0

Thermo Scientific Heratherm 13-3

Table 13-2 Technical Data - Series IMH 400-S / IMH 750-S

Parameter	Unit	IMH 400-S	IMH 750-S
Power line voltage +/- 10 %	V	120	
IP protection system		IP 20	
Protection class		I	
Overvoltage category as per IEC 60364-4-443		II	
Device fusing, building side	А	16	i
Device fusing, on PCB	А	2 x	16
Environmental conditions			
Min. ambient temperature	°C/°F	18/6	65
Max. ambient temperature	°C/°F	32/9	90
Max. humidity in service, non condensing	% r.F./ % r.H.	80, nicht kondensierend	
Min. storage temperature	°C/°F	20/68	
Max. storage temperature	°C/°F	60/140	
Max. humidity in storage, non condensing	% r.F./ % r.H.	90, nicht kondensierend	
Post-transport acclimation time	h	2	
Noise level	dB(A)	36 dB	
Degree of pollution as per IEC EN 61010-1		2	
Site conditions			
Maximum altitude above sea level	m/y ASL	2000/2187	
Minimum side clearance	mm/in	50/2	
Minimum front clearance	mm/in	810 / 31.9 670 / 26.4	
Minimum back wall clearance	mm/in	120/4.7	
Minimum top clearance	mm/in	200/7.9	

13-4 Heratherm Thermo Scientific

Spare Parts and Accessories

Material No.	Description
50126665	Stacking adapter Heratherm 60L
50126666	Stacking adapter Heratherm 100L
50126667	Stacking adapter Heratherm 180L
50127146	Fresh air filter IMH / IMH-S
50127431	Outer door for Heratherm IGS 60, IMH 60, IMH 60-S, OMS 60, OMH 60, OMH 60-S, OGS 60, OGH 60 and OGH 60-S with a door stop on the left side
50127432	Outer door for Heratherm IGS 100, IMH 100, IMH 100-S, OMS 100, OMH 100, OMH 100-S, OGS 100, OGH 100 and OGH 100-S with a door stop on the left side
50127433	Outer door for Heratherm IGS 180, IMH 180, IMH 180-S, OMS 180, OMH 180, OMH 180-S, OGS 180, OGH 180 and OGH 180-S with a door stop on the left side
50127434	Outer door for Heratherm IGS 60, IMH 60, IMH 60-S, OMS 60, OMH 60, OMH 60-S, OGS 60, OGH 60 and OGH 60-S with a door stop on the right side
50127435	Staple feet Heratherm incubators and heating & drying ovens
50127436	Kit door seal for Heratherm 60L
50127437	Kit door seal for Heratherm 100L
50127438	Kit door seal for Heratherm 180L
50127439	Kit door clips for Heratherm 60L / 100L / 180 L
50127440	Door inner panel seal for Heratherm IMH 60 and IMH 60-S
50127441	Door inner panel seal for Heratherm IMH 100 and IMH 100-S
50127442	Door inner panel seal for Heratherm IMH 180 and IMH 180-S
50127443	Levelling feet for Heratherm incubators and heating & drying ovens
50127444	Kit door handle right Heratherm incubators and heating & drying ovens
50127445	Kit door handle left Heratherm incubators and heating & drying ovens
50127446	Kit door handle with lock right Heratherm incubators and heating & drying ovens
50127447	Kit door handle with lock left Heratherm incubators and heating & drying ovens
50127449	Temperature sensor for Heratherm incubators

Thermo Scientific Heratherm 14-1

Material No.	Description
50127450	Upper door hinge for Heratherm incubators and heating & drying ovens
50127451	Lower door hinge for Heratherm incubators and heating & drying ovens
50127455	Outer door for Heratherm IGS 100, IMH 100, IMH 100-S, OMS 100, OMH 100, OMH 100-S, OGS 100, OGH 100 and OGH 100-S with a door stop on the right side
50127456	Outer door for Heratherm IGS 180, IMH 180, IMH 180-S, OMS 180, OMH 180, OMH 180-S, OGS 180, OGH 180 and OGH 180-S with a door stop on the right side
50127457	Kit operating panel Heratherm General Protocol incubators and heating & drying ovens
50127458	Kit operating panel Heratherm Advanced Protocol and Advanced Protocol Security incubators and heating & drying ovens
50127461	Kit electronic insert Heratherm General Protocol incubators and heating & drying ovens
50127462	Kit electronic insert Heratherm Advanced Protocol and Advanced Protocol Security incubators and heating & drying ovens without main board fan.
50127463	Mainboard cable for Heratherm incubators and heating & drying ovens
50127468	Glass door hinges for Heratherm incubators
50127469	Door switch for the right side of Heratherm incubators and heating & drying ovens
50127470	Door switch for the left side of Heratherm incubators and heating & drying oven
50127472	Kit glass door IGS 60, IMH 60, IMH 60-S
50127473	Kit glass door IGS 100, IMH 100, IMH 100-S
50127474	Kit glass door IGS 180, IMH 180, IMH 180-S
50127478	Temperatur limiter for IMH-S
50127480	Door hook catch for Heratherm incubators and heating & drying ovens with a door stop on the right side
50127481	Door hook catch for Heratherm incubators and heating & drying ovens with a door stop on the left side
50127482	Magnetic door hook catch for Heratherm incubators and heating & drying ovens with a door stop on the right side
50127483	Magnetic door hook catch for Heratherm incubators and heating & drying ovens with a door stop on the left sid
50127497	Heating coils for Heratherm IMH-S at a voltage of 120 V
50127498	Heating coils for Heratherm IMH-S at a voltage of 230 V
50127504	Kit Heating coils IMH 120 V
50127511	Kit fan system Heratherm IMH 60 IMH 60-S, OMH 60 and OMH 60-S, IMH 100, IMH 100-S, IMH 180 and OMH 180-S, 120 V.
50127515	Kit Heating coils IMH 230 V
50127531	Circulation fan IMH / IMH-S D = 160 mm (6.3 inch), H = 28 mm (1.1 inch)

14-2 Heratherm Thermo Scientific

Material No.	Description
50127555	Kit fan system Heratherm IMH 60 IMH 60-S, OMH 60 and OMH 60-S, IMH 100, IMH 100-S, IMH 180 and OMH 180-S, 230 V
50127567	Fresh air filter for IMH, IMH-S
50127741	Support stand with castors for Heratherm 60L
50127742	Support stand with castors for Heratherm 100L
50127743	Support stand with castors for Heratherm 180L
50127764	Wire mesh shelf IGS 60/100/180, IMH 60/100/180, IMH 60/100/180-S, including 2 shelf supports
50127768	Sample sensor for IMH 60-S / IMH 100-S / IMH 180-S
50127770	Stainless steel perforated shelf IGS 60, including 2 shelf supports
50127771	Stainless steel perforated shelf IGS 100, including 2 shelf supports
50127772	Stainless steel perforated shelf IGS 180, including 2 shelf supports
50127773	Stainless steel perforated shelf IMH 60/IMH 60-S/OMH 60/OMH 60-S/OMS 60/100/180, including 2 shelf supports
50127774	Stainless steel perforated shelf IMH 100 / IMH 100-S / OMH 100 / OMH 100-S/OMS 60/100/180, including 2 shelf supports
50127777	Stainless steel perforated shelf IMH 180 / IMH 180-S / OMH 180 / OMH 180-S/OMS 60/100/180, including 2 shelf supports
50127861	Retaining springs for Heratherm incubators and heating & drying ovens
50127862	Support rail for Heratherm IGS 60, IMH 60, IMH 60-S, OMS 60, OMH 60, OMH 60-S
50127863	Support rail for Heratherm IGS 100, IMH 100, IMH 100-S, OMS 100, OMH 100, OMH 100-S
50127864	Support rail for Heratherm IGS 180, IMH 180, IMH 180-S, OMS 180, OMH 180, OMH 180-S
50128179	Kit Fuses T2A Incubators 120 V
50128184	Sample sensor connection for Heratherm incubators and heating & drying ovens
50128197	Power socket for Heratherm IMH
50128203	Kit Fuses T2A Incubators 230 V
50128212	Fuse holder for Heratherm IMH
50128237	Kit Key for door handle with lock Heratherm
50128265	Lowenstein holder IGS 100 / IMH 10 / IMH 100-S / IGS 180 / IMH 180 / IMH 180-S
50128683	Drip tray IGS / IMH / IMH-S 60L
50128704	Kit Anti-tilt anchor
50128791	Drip tray IGS / IMH / IMH-S 100L
50128792	Drip tray IGS / IMH / IMH-S 180L

Thermo Scientific Heratherm 14-3

Glass door lock for Heratherm IGS, IMH und IMH-S 50128793 Petri dish holder 50 mm (2 inch) IGS 60 / IMH 60 / IMH 60-S 50128794 Petri dish holder 50 mm (2 inch) IGS 100 / IMH 100 / IMH 100-S 50128815 Petri dish holder 50 mm (2 inch) IGS 180 / IMH 180 / IMH 180-S 50128816 Petri dish holder 90 mm (3.54 inch) IGS 60 / IMH 60 / IMH 60-S 50128818 Petri dish holder 90 mm (3.54 inch) IGS 100 / IMH 100 / IMH 100-S 50128819 Petri dish holder 90 mm (3.54 inch) IGS 100 / IMH 100 / IMH 100-S 50128819 Petri dish holder 90 mm (3.54 inch) IGS 180 / IMH 180 / IMH 180-S 50128960 Kit Shelving system with glass door lock IGS 60, IMH 60, IMH 60-S 50128961 Kit Shelving system with glass door lock IGS 100, IMH 100, IMH 100-S 50128962 Kit Shelving system with glass door lock IGS 180, IMH 180, IMH 180-S 50130657 Kit Viton door seal 60 L Heratherm 50130658 Kit Viton door seal 100 L Heratherm 50130659 Kit Viton door seal 180 L Heratherm 50134116 Kit Heating coil IGS 400, 750 120 V, centre 50134117 Kit Heating coil IGS 400, 750 230 V, centre 50134120 Kit Heating coil IGS 750 120 V, outside 50134121 Kit Heating coil IMH 750 120 V, ring heater 50134122 Kit Heating coil IMH 750 120 V, ring heater 50134123 Kit pass door cpl 400 / 750 50134322 Kit glass door cpl 400 / 750 50134322 Kit glass door cpl 400 HTM 50134323 Kit glass door cpl 400 HTM 50134324 Kit glass door cpl 400 HTM 50134325 Kit glass door cpl 400 HTM 50134326 Door gasket 750 L HTM 50134327 Door gasket 750 L HTM 50134338 Kit castors 400 / 750 HTM 50134339 Kit castors 400 / 750 HTM 50134330 Kit castors 400 / 750 HTM 50134330 Kit electric insert cpl HTM 400 / 750 50135044 Kit electric insert cpl HTM 9 PH floor 50135056 Kit door lock 750 left cpl HTM	Material No.	Description
50128794 Petri dish holder 50 mm (2 inch) IGS 100 / IMH 100 / IMH 100-S 50128815 Petri dish holder 50 mm (2 inch) IGS 180 / IMH 180 / IMH 180-S 50128816 Petri dish holder 90 mm (3.54 inch) IGS 60 / IMH 60 / IMH 60-S 50128818 Petri dish holder 90 mm (3.54 inch) IGS 100 / IMH 100 / IMH 100-S 50128819 Petri dish holder 90 mm (3.54 inch) IGS 180 / IMH 180 / IMH 180-S 50128960 Kit Shelving system with glass door lock IGS 60, IMH 60, IMH 100-S 50128961 Kit Shelving system with glass door lock IGS 100, IMH 100, IMH 100-S 50128962 Kit Shelving system with glass door lock IGS 180, IMH 180, IMH 180-S 50130657 Kit Viton door seal 60 L Heratherm 50130658 Kit Viton door seal 180 L Heratherm 50134116 Kit Heating coil IGS 400, 750 120 V, centre 50134117 Kit Heating coil IGS 400, 750 120 V, centre 50134120 Kit Heating coil IGS 750 120 V, outside 50134121 Kit Heating coil IMH 750 120 V, ring heater 50134322 Kit plass door cpl 400 HTM 50134323 Kit glass door ri / le cpl 750 HTM 50134324 Door gasket 750 L HTM 50134325 Kit door conn. clips 400 / 750 HTM 50134328		Glass door lock for Heratherm IGS, IMH und IMH-S
50128815 Petri dish holder 50 mm (2 inch) IGS 180 / IMH 180 / IMH 180-S 50128816 Petri dish holder 90 mm (3.54 inch) IGS 60 / IMH 60 / IMH 60-S 50128818 Petri dish holder 90 mm (3.54 inch) IGS 100 / IMH 100 / IMH 100-S 50128819 Petri dish holder 90 mm (3.54 inch) IGS 180 / IMH 180 / IMH 180-S 50128960 Kit Shelving system with glass door lock IGS 60, IMH 60, IMH 60-S 50128961 Kit Shelving system with glass door lock IGS 100, IMH 100, IMH 100-S 50128962 Kit Shelving system with glass door lock IGS 180, IMH 180, IMH 180-S 50130657 Kit Viton door seal 60 L Heratherm 50130658 Kit Viton door seal 180 L Heratherm 501340659 Kit Viton door seal 180 L Heratherm 50134116 Kit Heating coil IGS 400, 750 120 V, centre 50134121 Kit Heating coil IGS 400, 750 230 V, centre 50134122 Kit Heating coil IGS 750 230 V, outside 50134123 Kit heating coil IMH 750 120 V, ring heater 50134322 Kit glass door rpl 400 HTM 50134323 Kit glass door ri / le cpl 750 HTM 50134326 Door gasket 750 L HTM 50134327 Door gasket 750 L HTM 50134328 Kit door conn. clips 400 / 750 HT	50128793	Petri dish holder 50 mm (2 inch) IGS 60 / IMH 60 / IMH 60-S
50128816 Petri dish holder 90 mm (3.54 inch) IGS 60 / IMH 60 / IMH 60-S 50128818 Petri dish holder 90 mm (3.54 inch) IGS 100 / IMH 100 / IMH 100-S 50128819 Petri dish holder 90 mm (3.54 inch) IGS 180 / IMH 180 / IMH 180-S 50128960 Kit Shelving system with glass door lock IGS 60, IMH 60, IMH 60-S 50128961 Kit Shelving system with glass door lock IGS 100, IMH 100, IMH 100-S 50128962 Kit Shelving system with glass door lock IGS 180, IMH 180, IMH 180-S 50130657 Kit Viton door seal 60 L Heratherm 50130658 Kit Viton door seal 180 L Heratherm 50130659 Kit Viton door seal 180 L Heratherm 50134116 Kit Heating coil IGS 400, 750 120 V, centre 50134120 Kit Heating coil IGS 400, 750 230 V, centre 50134121 Kit Heating coil IGS 750 230 V, outside 50134122 Kit Heating coil IMH 750 120 V, ring heater 50134323 Kit plass door cpl 400 / 750 50134324 Kit glass door ri / le cpl 750 HTM 50134325 Kit glass door ri / le cpl 750 HTM 50134326 Door gasket 750 L HTM 50134327 Door gasket 750 L HTM 50134328 Kit glass door ri / le cpl 750 HTM	50128794	Petri dish holder 50 mm (2 inch) IGS 100 / IMH 100 / IMH 100-S
50128818 Petri dish holder 90 mm (3.54 inch) IGS 100 / IMH 100 / IMH 100-S 50128819 Petri dish holder 90 mm (3.54 inch) IGS 180 / IMH 180 / IMH 180-S 50128960 Kir Shelving system with glass door lock IGS 60, IMH 60, IMH 60-S 50128961 Kir Shelving system with glass door lock IGS 100, IMH 100, IMH 100-S 50128962 Kir Shelving system with glass door lock IGS 180, IMH 180, IMH 180-S 50130657 Kir Viton door seal 60 L Heratherm 50130658 Kir Viton door seal 180 L Heratherm 50134059 Kir Viton door seal 180 L Heratherm 50134116 Kir Heating coil IGS 400, 750 120 V, centre 50134117 Kir Heating coil IGS 750 120 V, outside 50134120 Kir Heating coil IGS 750 120 V, outside 50134121 Kir Heating coil IMH 750 120 V, ring heater 50134122 Kir Heating coil IMH 750 230 V, ring heater 50134321 Kir glass door ri / le cpl 750 HTM 50134322 Kir glass door ri / le cpl 750 HTM 50134323 Kir glass door ri / le cpl 750 HTM 50134329 Kir door conn. clips 400 / 750 HTM 50134333 Kir castors 400 / 750 HTM 50134334 Kir shelf rack set cpl HTM 400 / 750	50128815	Petri dish holder 50 mm (2 inch) IGS 180 / IMH 180 / IMH 180-S
50128819 Petri dish holder 90 mm (3.54 inch) IGS 180 / IMH 180 / IMH 180-S 50128960 Kit Shelving system with glass door lock IGS 60, IMH 60, IMH 60-S 50128961 Kit Shelving system with glass door lock IGS 100, IMH 100, IMH 100-S 50128962 Kit Shelving system with glass door lock IGS 180, IMH 180, IMH 180-S 50130657 Kit Viton door seal 60 L Heratherm 50130658 Kit Viton door seal 180 L Heratherm 50130659 Kit Viton door seal 180 L Heratherm 50134116 Kit Heating coil IGS 400, 750 120 V, centre 50134117 Kit Heating coil IGS 400, 750 230 V, centre 50134120 Kit Heating coil IGS 750 120 V, outside 50134121 Kit Heating coil IMH 750 120 V, ring heater 50134122 Kit Heating coil IMH 750 120 V, ring heater 50134323 Kit glass door cpl 400 / 750 50134322 Kit glass door cpl 400 HTM 50134323 Kit glass door ri / le cpl 750 HTM 50134324 Door gasket 400 L HTM 50134325 Kit door conn. clips 400 / 750 HTM 50134333 Kit castors 400 / 750 HTM 50134334 Kit shelf rack set cpl HTM 400 / 750 50135044 Kit electric in	50128816	Petri dish holder 90 mm (3.54 inch) IGS 60 / IMH 60 / IMH 60-S
Kit Shelving system with glass door lock IGS 60, IMH 60, IMH 60-S 50128961 Kit Shelving system with glass door lock IGS 100, IMH 100, IMH 100-S 50128962 Kit Shelving system with glass door lock IGS 180, IMH 180, IMH 180-S 50130657 Kit Viton door seal 60 L Heratherm 50130658 Kit Viton door seal 100 L Heratherm 50130659 Kit Viton door seal 180 L Heratherm 50134116 Kit Heating coil IGS 400, 750 120 V, centre 50134117 Kit Heating coil IGS 400, 750 230 V, centre 50134120 Kit Heating coil IGS 750 120 V, outside 50134121 Kit Heating coil IGS 750 120 V, ring heater 50134122 Kit Heating coil IMH 750 120 V, ring heater 50134123 Kit heating coil IMH 750 120 V, ring heater 50134315 Kit DS bus cable cpl 400 / 750 50134322 Kit glass door cpl 400 HTM 50134323 Kit glass door ri / le cpl 750 HTM 50134326 Door gasket 400 L HTM 50134327 Door gasket 750 L HTM 50134328 Kit profile gasket 750 L HTM 50134330 Kit castors 400 / 750 HTM 50134331 Kit castors 400 / 750 HTM 50134333 Kit castors 400 / 750 HTM 50134334 Kit shelf rack set cpl HTM 400 / 750 50135043 Kit electric insert cpl HTM H floorstand Kit glass door hinges cpl 400/750 L HTM	50128818	Petri dish holder 90 mm (3.54 inch) IGS 100 / IMH 100 / IMH 100-S
Kit Shelving system with glass door lock IGS 100, IMH 100, IMH 100-S 50128962 Kit Shelving system with glass door lock IGS 180, IMH 180, IMH 180-S Kit Viton door seal 60 L Heratherm 50130658 Kit Viton door seal 100 L Heratherm 50130659 Kit Viton door seal 180 L Heratherm 50134116 Kit Heating coil IGS 400, 750 120 V, centre 50134117 Kit Heating coil IGS 400, 750 230 V, centre 50134120 Kit Heating coil IGS 750 120 V, outside 50134121 Kit Heating coil IGS 750 230 V, outside 50134122 Kit Heating coil IMH 750 120 V, ring heater 50134123 Kit heating coil IMH 750 230 V, ring heater 50134124 Kit glass door cpl 400 / 750 50134325 Kit glass door ri / le cpl 750 HTM 50134326 Door gasket 400 L HTM 50134327 Door gasket 750 L HTM 50134328 Kit profile gasket 750 L HTM 50134330 Kit castors 400 / 750 HTM 50134331 Kit castors 400 / 750 HTM 50134332 Kit door conn. clips 400 / 750 HTM 50134333 Kit castors 400 / 750 HTM 50134334 Kit electric insert cpl HTM H floorstand Kit electric insert cpl HTM S floorstand Kit electric insert cpl HTM 3 PH floor 50135056 Kit glass door hinges cpl 400/750 L HTM	50128819	Petri dish holder 90 mm (3.54 inch) IGS 180 / IMH 180 / IMH 180-S
50128962 Kit Shelving system with glass door lock IGS 180, IMH 180, IMH 180-S 50130657 Kit Viton door seal 60 L Heratherm 50130658 Kit Viton door seal 100 L Heratherm 50130659 Kit Viton door seal 180 L Heratherm 50134116 Kit Heating coil IGS 400, 750 120 V, centre 50134117 Kit Heating coil IGS 400, 750 230 V, centre 50134120 Kit Heating coil IGS 750 120 V, outside 50134121 Kit Heating coil IMH 750 120 V, ring heater 50134122 Kit Heating coil IMH 750 230 V, ring heater 50134123 Kit pass door cpl 400 / 750 50134324 Kit glass door cpl 400 HTM 50134325 Kit glass door rj le cpl 750 HTM 50134326 Door gasket 400 L HTM 50134327 Door gasket 750 L HTM 50134328 Kit profile gasket 750 L HTM 50134333 Kit castors 400 / 750 HTM 50134334 Kit shelf rack set cpl HTM 400 / 750 50135044 Kit electric insert cpl HTM 5 floorstand 50135056 Kit glass door hinges cpl 400/750 L HTM	50128960	Kit Shelving system with glass door lock IGS 60, IMH 60, IMH 60-S
50128962 Kit Shelving system with glass door lock IGS 180, IMH 180, IMH 180-S 50130657 Kit Viton door seal 60 L Heratherm 50130658 Kit Viton door seal 100 L Heratherm 50130659 Kit Viton door seal 180 L Heratherm 50134116 Kit Heating coil IGS 400, 750 120 V, centre 50134117 Kit Heating coil IGS 400, 750 230 V, centre 50134120 Kit Heating coil IGS 750 120 V, outside 50134121 Kit Heating coil IMH 750 120 V, ring heater 50134122 Kit Heating coil IMH 750 230 V, ring heater 50134123 Kit pass door cpl 400 / 750 50134324 Kit glass door cpl 400 HTM 50134325 Kit glass door rj le cpl 750 HTM 50134326 Door gasket 400 L HTM 50134327 Door gasket 750 L HTM 50134328 Kit profile gasket 750 L HTM 50134333 Kit castors 400 / 750 HTM 50134334 Kit shelf rack set cpl HTM 400 / 750 50135044 Kit electric insert cpl HTM 5 floorstand 50135056 Kit glass door hinges cpl 400/750 L HTM	50128961	
50130657 Kit Viton door seal 60 L Heratherm 50130658 Kit Viton door seal 100 L Heratherm 50130659 Kit Viton door seal 180 L Heratherm 50134116 Kit Heating coil IGS 400, 750 120 V, centre 50134117 Kit Heating coil IGS 400, 750 230 V, centre 50134120 Kit Heating coil IGS 750 120 V, outside 50134121 Kit Heating coil IGS 750 230 V, outside 50134122 Kit Heating coil IMH 750 120 V, ring heater 50134123 Kit beating coil IMH 750 230 V, ring heater 50134315 Kit DS bus cable cpl 400 / 750 50134322 Kit glass door cpl 400 HTM 50134323 Kit glass door ri / le cpl 750 HTM 50134326 Door gasket 400 L HTM 50134327 Door gasket 750 L HTM 50134328 Kit profile gasket 750 L HTM 50134333 Kit castors 400 / 750 HTM 50135043 Kit electric insert cpl HTM 400 / 750 50135044 Kit electric insert cpl HTM 5 floorstand 50135056 Kit glass door hinges cpl 400/750 L HTM		
50130658 Kit Viton door seal 100 L Heratherm 50130659 Kit Viton door seal 180 L Heratherm 50134116 Kit Heating coil IGS 400, 750 120 V, centre 50134117 Kit Heating coil IGS 400, 750 230 V, centre 50134120 Kit Heating coil IGS 750 120 V, outside 50134121 Kit Heating coil IMF 750 120 V, outside 50134122 Kit Heating coil IMH 750 120 V, ring heater 50134123 Kit heating coil IMH 750 230 V, ring heater 50134321 Kit glass door cpl 400 / 750 50134322 Kit glass door cpl 400 HTM 50134323 Kit glass door ri / le cpl 750 HTM 50134326 Door gasket 400 L HTM 50134327 Door gasket 750 L HTM 50134328 Kit profile gasket 750 L HTM 50134333 Kit castors 400 / 750 HTM 50134333 Kit castors 400 / 750 HTM 50135043 Kit electric insert cpl HTM H floorstand 50135055 Kit electric insert cpl HTM 3 PH floor 50135056 Kit glass door hinges cpl 400/750 L HTM		
50130659 Kit Viton door seal 180 L Heratherm 50134116 Kit Heating coil IGS 400, 750 120 V, centre 50134117 Kit Heating coil IGS 400, 750 230 V, centre 50134120 Kit Heating coil IGS 750 120 V, outside 50134121 Kit Heating coil IMF 750 230 V, outside 50134122 Kit Heating coil IMH 750 120 V, ring heater 50134123 Kit heating coil IMH 750 230 V, ring heater 50134315 Kit DS bus cable cpl 400 / 750 50134322 Kit glass door cpl 400 HTM 50134323 Kit glass door ri / le cpl 750 HTM 50134326 Door gasket 400 L HTM 50134327 Door gasket 750 L HTM 50134328 Kit profile gasket 750 L HTM 50134329 Kit door conn. clips 400 / 750 HTM 50134333 Kit castors 400 / 750 HTM 50135043 Kit electric insert cpl HTM 4 floorstand 50135044 Kit electric insert cpl HTM S floorstand 50135056 Kit glass door hinges cpl 400/750 L HTM		Kit Viton door seal 100 L Heratherm
50134116 Kit Heating coil IGS 400, 750 120 V, centre 50134117 Kit Heating coil IGS 400, 750 230 V, centre 50134120 Kit Heating coil IGS 750 120 V, outside 50134121 Kit Heating coil IGS 750 230 V, outside 50134122 Kit Heating coil IMH 750 120 V, ring heater 50134123 Kit heating coil IMH 750 230 V, ring heater 50134315 Kit DS bus cable cpl 400 / 750 50134322 Kit glass door rpl 400 HTM 50134323 Kit glass door ri / le cpl 750 HTM 50134326 Door gasket 400 L HTM 50134327 Door gasket 750 L HTM 50134328 Kit pofile gasket 750 L HTM 501343329 Kit door conn. clips 400 / 750 HTM 50134334 Kit shelf rack set cpl HTM 400 / 750 50135043 Kit electric insert cpl HTM H floorstand 50135044 Kit electric insert cpl HTM S floorstand 50135056 Kit glass door hinges cpl 400/750 L HTM		Kit Viton door seal 180 L Heratherm
50134117 Kit Heating coil IGS 400, 750 230 V, centre 50134120 Kit Heating coil IGS 750 120 V, outside 50134121 Kit Heating coil IGS 750 230 V, outside 50134122 Kit Heating coil IMH 750 120 V, ring heater 50134123 Kit heating coil IMH 750 230 V, ring heater 50134315 Kit DS bus cable cpl 400 / 750 50134322 Kit glass door cpl 400 HTM 50134323 Kit glass door ri / le cpl 750 HTM 50134326 Door gasket 400 L HTM 50134327 Door gasket 750 L HTM 50134328 Kit profile gasket 750 L HTM 50134332 Kit door conn. clips 400 / 750 HTM 50134333 Kit castors 400 / 750 HTM 50134334 Kit shelf rack set cpl HTM 400 / 750 50135043 Kit electric insert cpl HTM H floorstand 50135055 Kit electric insert cpl HTM 3 PH floor 50135056 Kit glass door hinges cpl 400/750 L HTM		
50134120 Kit Heating coil IGS 750 120 V, outside 50134121 Kit Heating coil IGS 750 230 V, outside 50134122 Kit Heating coil IMH 750 120 V, ring heater 50134123 Kit heating coil IMH 750 230 V, ring heater 50134315 Kit DS bus cable cpl 400 / 750 50134322 Kit glass door cpl 400 HTM 50134323 Kit glass door ri / le cpl 750 HTM 50134326 Door gasket 400 L HTM 50134327 Door gasket 750 L HTM 50134328 Kit profile gasket 750 L HTM 50134329 Kit door conn. clips 400 / 750 HTM 50134333 Kit castors 400 / 750 HTM 50135043 Kit electric insert cpl HTM 400 / 750 50135044 Kit electric insert cpl HTM S floorstand 50135055 Kit electric insert cpl HTM 3 PH floor 50135056 Kit glass door hinges cpl 400/750 L HTM		
Kit Heating coil IMH 750 120 V, ring heater 50134123 Kit heating coil IMH 750 230 V, ring heater 50134315 Kit DS bus cable cpl 400 / 750 50134322 Kit glass door cpl 400 HTM 50134323 Kit glass door ri / le cpl 750 HTM 50134326 Door gasket 400 L HTM 50134327 Door gasket 750 L HTM 50134328 Kit profile gasket 750 L HTM 50134329 Kit door conn. clips 400 / 750 HTM 50134331 Kit castors 400 / 750 HTM 50134334 Kit shelf rack set cpl HTM 400 / 750 50135043 Kit electric insert cpl HTM H floorstand 50135044 Kit electric insert cpl HTM S floorstand 50135055 Kit electric insert cpl HTM 3 PH floor 50135056 Kit glass door hinges cpl 400/750 L HTM		
50134123 Kit heating coil IMH 750 230 V, ring heater 50134315 Kit DS bus cable cpl 400 / 750 50134322 Kit glass door cpl 400 HTM 50134323 Kit glass door ri / le cpl 750 HTM 50134326 Door gasket 400 L HTM 50134327 Door gasket 750 L HTM 50134328 Kit profile gasket 750 L HTM 50134329 Kit door conn. clips 400 / 750 HTM 50134333 Kit castors 400 / 750 HTM 50134334 Kit shelf rack set cpl HTM 400 / 750 50135043 Kit electric insert cpl HTM H floorstand 50135055 Kit electric insert cpl HTM 3 PH floor 50135056 Kit glass door hinges cpl 400/750 L HTM	50134121	Kit Heating coil IGS 750 230 V, outside
50134315 Kit DS bus cable cpl 400 / 750 50134322 Kit glass door cpl 400 HTM 50134323 Kit glass door ri / le cpl 750 HTM 50134326 Door gasket 400 L HTM 50134327 Door gasket 750 L HTM 50134328 Kit profile gasket 750 L HTM 50134329 Kit door conn. clips 400 / 750 HTM 50134333 Kit castors 400 / 750 HTM 50134334 Kit shelf rack set cpl HTM 400 / 750 50135043 Kit electric insert cpl HTM B floorstand 50135055 Kit electric insert cpl HTM 3 PH floor 50135056 Kit glass door hinges cpl 400/750 L HTM	50134122	Kit Heating coil IMH 750 120 V, ring heater
50134322 Kit glass door cpl 400 HTM 50134323 Kit glass door ri / le cpl 750 HTM 50134326 Door gasket 400 L HTM 50134327 Door gasket 750 L HTM 50134328 Kit profile gasket 750 L HTM 50134329 Kit door conn. clips 400 / 750 HTM 50134330 Kit castors 400 / 750 HTM 50134334 Kit shelf rack set cpl HTM 400 / 750 50135043 Kit electric insert cpl HTM H floorstand 50135044 Kit electric insert cpl HTM S floorstand 50135055 Kit electric insert cpl HTM 3 PH floor 50135056 Kit glass door hinges cpl 400/750 L HTM	50134123	Kit heating coil IMH 750 230 V, ring heater
50134323 Kit glass door ri / le cpl 750 HTM 50134326 Door gasket 400 L HTM 50134327 Door gasket 750 L HTM 50134328 Kit profile gasket 750 L HTM 50134329 Kit door conn. clips 400 / 750 HTM 50134333 Kit castors 400 / 750 HTM 50134334 Kit shelf rack set cpl HTM 400 / 750 50135043 Kit electric insert cpl HTM H floorstand 50135044 Kit electric insert cpl HTM S floorstand 50135055 Kit electric insert cpl HTM 3 PH floor 50135056 Kit glass door hinges cpl 400/750 L HTM	50134315	Kit DS bus cable cpl 400 / 750
50134326 Door gasket 400 L HTM 50134327 Door gasket 750 L HTM 50134328 Kit profile gasket 750 L HTM 50134329 Kit door conn. clips 400 / 750 HTM 50134333 Kit castors 400 / 750 HTM 50134334 Kit shelf rack set cpl HTM 400 / 750 50135043 Kit electric insert cpl HTM H floorstand 50135044 Kit electric insert cpl HTM S floorstand 50135055 Kit electric insert cpl HTM 3 PH floor 50135056 Kit glass door hinges cpl 400/750 L HTM	50134322	Kit glass door cpl 400 HTM
50134327 Door gasket 750 L HTM 50134328 Kit profile gasket 750 L HTM 50134329 Kit door conn. clips 400 / 750 HTM 50134333 Kit castors 400 / 750 HTM 50134334 Kit shelf rack set cpl HTM 400 / 750 50135043 Kit electric insert cpl HTM H floorstand 50135044 Kit electric insert cpl HTM S floorstand 50135055 Kit electric insert cpl HTM 3 PH floor 50135056 Kit glass door hinges cpl 400/750 L HTM	50134323	Kit glass door ri / le cpl 750 HTM
50134328 Kit profile gasket 750 L HTM 50134329 Kit door conn. clips 400 / 750 HTM 50134333 Kit castors 400 / 750 HTM 50134334 Kit shelf rack set cpl HTM 400 / 750 50135043 Kit electric insert cpl HTM H floorstand 50135044 Kit electric insert cpl HTM S floorstand 50135055 Kit electric insert cpl HTM 3 PH floor 50135056 Kit glass door hinges cpl 400/750 L HTM	50134326	Door gasket 400 L HTM
50134329 Kit door conn. clips 400 / 750 HTM 50134333 Kit castors 400 / 750 HTM 50134334 Kit shelf rack set cpl HTM 400 / 750 50135043 Kit electric insert cpl HTM H floorstand 50135044 Kit electric insert cpl HTM S floorstand 50135055 Kit electric insert cpl HTM 3 PH floor 50135056 Kit glass door hinges cpl 400/750 L HTM	50134327	Door gasket 750 L HTM
50134333 Kit castors 400 / 750 HTM 50134334 Kit shelf rack set cpl HTM 400 / 750 50135043 Kit electric insert cpl HTM H floorstand 50135044 Kit electric insert cpl HTM S floorstand 50135055 Kit electric insert cpl HTM 3 PH floor 50135056 Kit glass door hinges cpl 400/750 L HTM	50134328	Kit profile gasket 750 L HTM
50134334 Kit shelf rack set cpl HTM 400 / 750 50135043 Kit electric insert cpl HTM H floorstand 50135044 Kit electric insert cpl HTM S floorstand 50135055 Kit electric insert cpl HTM 3 PH floor 50135056 Kit glass door hinges cpl 400/750 L HTM	50134329	Kit door conn. clips 400 / 750 HTM
50135043 Kit electric insert cpl HTM H floorstand 50135044 Kit electric insert cpl HTM S floorstand 50135055 Kit electric insert cpl HTM 3 PH floor 50135056 Kit glass door hinges cpl 400/750 L HTM	50134333	Kit castors 400 / 750 HTM
50135044 Kit electric insert cpl HTM S floorstand 50135055 Kit electric insert cpl HTM 3 PH floor 50135056 Kit glass door hinges cpl 400/750 L HTM	50134334	Kit shelf rack set cpl HTM 400 / 750
50135055 Kit electric insert cpl HTM 3 PH floor 50135056 Kit glass door hinges cpl 400/750 L HTM	50135043	Kit electric insert cpl HTM H floorstand
50135056 Kit glass door hinges cpl 400/750 L HTM	50135044	Kit electric insert cpl HTM S floorstand
	50135055	Kit electric insert cpl HTM 3 PH floor
50135058 Kit door lock 750 left cpl HTM	50135056	Kit glass door hinges cpl 400/750 L HTM
	50135058	Kit door lock 750 left cpl HTM

14-4 Heratherm Thermo Scientific

Material No.	Description
50135059	Kit door lock 750 right cpl HTM
50135060	Door outer casing left HTM 400
50135061	Door outer casing right HTM 400
50135062	Door outer casing left HTM 750
50135063	Door outer casing right HTM 750
50135150	Kit temp sensor cpl HTM incubator floor
50135153	Kit door lock 400 right cpl HTM
50135154	Kit door lock 400 left cpl HTM

Thermo Scientific Heratherm 14-5